

Tutorial: How to install the ink cartridge & the battery into the puck

- 1/ thanks to a flat screwdriver, remove the two screws at the top of the puck
- 2/ gently remove the hood of the puck (to avoid any destruction of the band of connectors)
- 3/ With the flat screwdriver, remove the 4 screws on the green base
- 4/ gently raise the green hood (to avoid any destruction of the band of connectors)
- 5/ Remove the ink cartridge of its pack
- 6/ Remove the protective film on the ink jet head
- 7/ place and clip the ink cartridge into the place provided on the green base
- 8/ Reposition the hood back on the base, then fix it with the 4 screws
- 9/ Take the battery and connect it to the connector, then place it between the 3 pins
- 10/ place the hood back then fix it with the 2 screws
- 11/ your inkjet air cushion puck is now ready to use !