Ex.A.O

Capteur Foxy / Tooxy / AirNeXT

Data logging

Sensor Foxy / Tooxy / AirNeXT

Réf :
482 093

Français – p 1

English – p 8

Version : 5204

Capteur Electrophy

Electrophy sensor

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09 69 32 02 10*

** appel non surtaxé*

You have just purchased an electrophy sensor and we thank you for your confidence. This document aims at guiding you in using your sensor.

1. Description

This electrophy sensor allows you to perform :

- heart rate measurements. These can be carried out both when the sensor is used in measurement mode, i.e. not connected to a PC, and in ExAO mode.
- ECG
- raw signal measurement (animal electrophysiology, human electrophysiology....)

2. Utilisation

This sensor can be used with any console in the Foxy range.

- Present the sensor facing the console connector.
- Slide the connector into place, following the console guide.
- Push until it clicks into place.

3. Measure

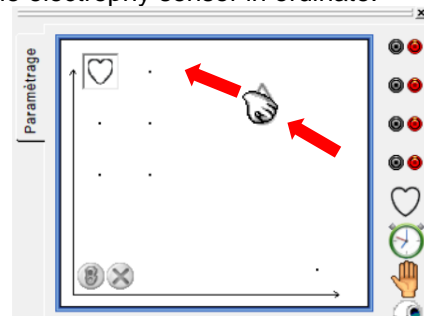
3.1 Using the sensor in standalone mode

Select the function using the orange button.


Unit is indicated by a triangle in the green band opposite the caliber.

3.2 Using the sensor with the Sientific Workshop software :

- Run the Standard Scientific Workshop software
- Drag and drop the electrophy sensor in ordinate:

The tabs specific to the sensor appear:

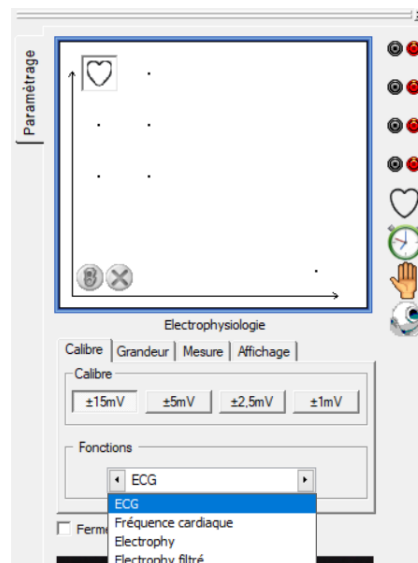3.3 Setting

The sensor is available in several functions.

Each function is adapted to a specific type of experiment.

3.3.1 Myotatic reflex

3.3.1.1 Connection :

Triangle electrodes

> 2 measurement electrodes placed on the chiefs of the gastrocnemius muscle

> 1 reference electrode on the gastrocnemius tendon.

Electrodes aligned

> 1 measuring electrode placed on one head of the gastrocnemius muscle, the other measuring electrode on the tendon part of this muscle.

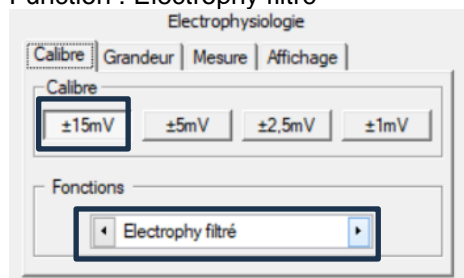
> The reference electrode is placed on the foot.


Connect the « Martoreflex » to the grounding terminal (black) and the synchro terminal (red) on the front of the Foxy console.

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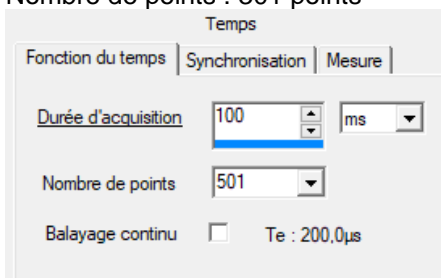
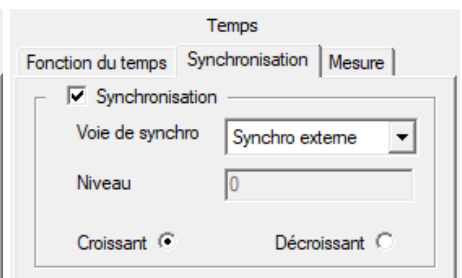
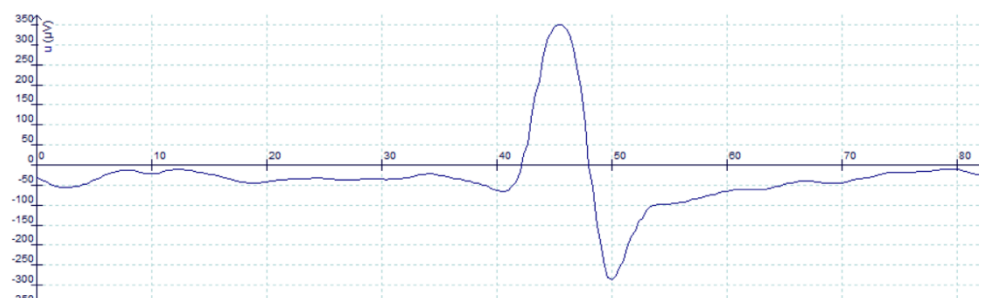
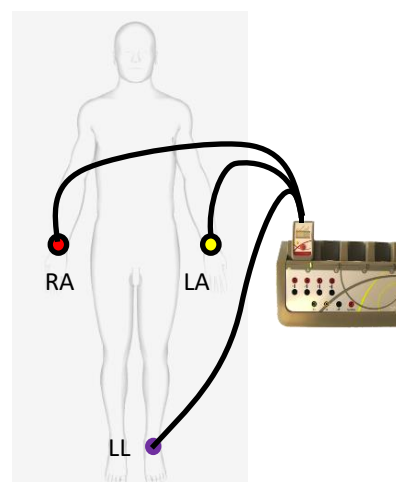
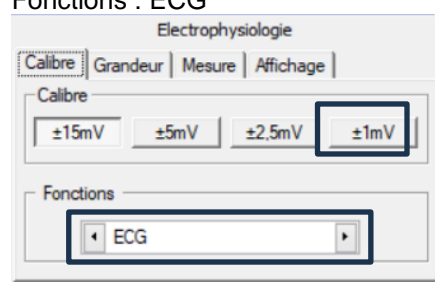
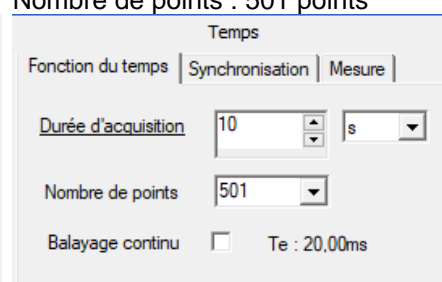
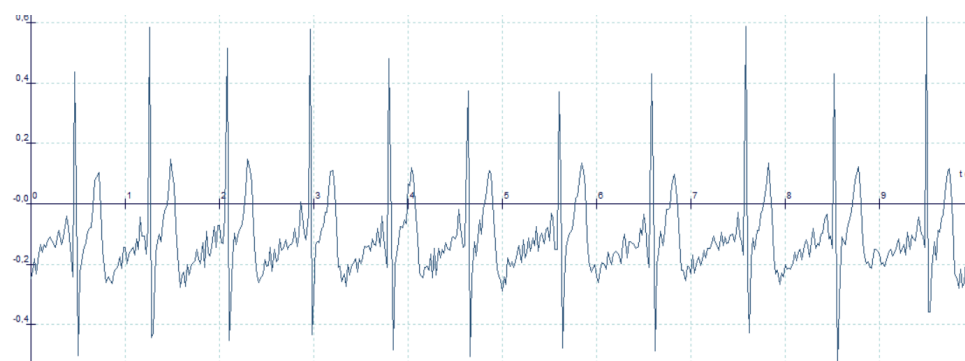
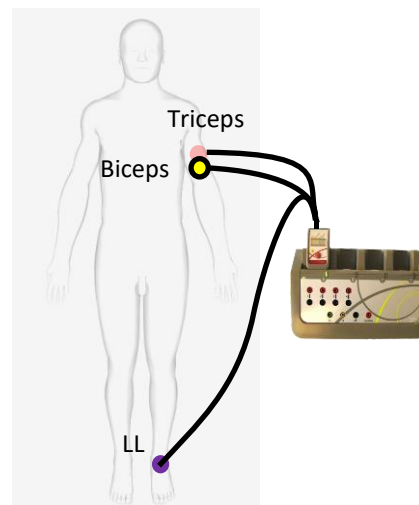
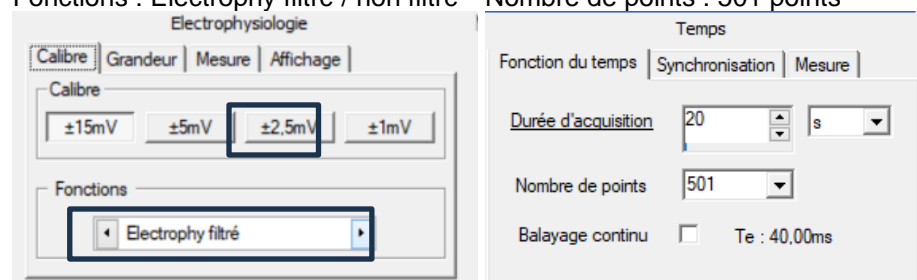
The filtered/unfiltered Electrophy function will produce similar results.

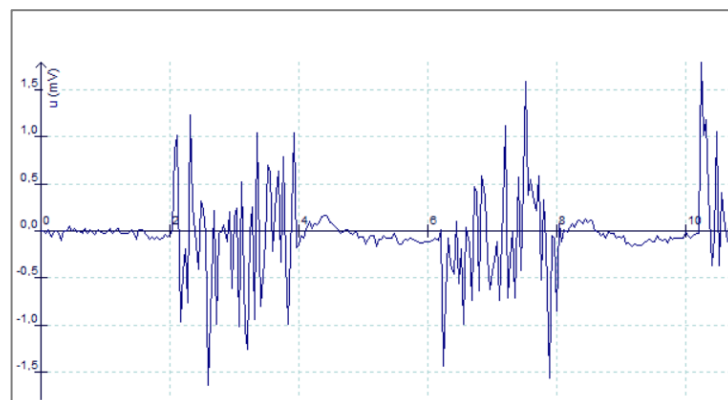3.3.3.3 Sample result :

4. Technical characteristics

Measurement range: 0 to 220 cpm

Max. sampling frequency: 135 μ s

Insulated to current standards.

Supplied with 50 adhesive ECG-EMG electrodes

5. After-Sale Services

The warranty is of 2 years.

For all settings, contact our **Technical Support** on **+33232294023**.

The material must be returned to our workshops and for all repairs or spare parts, kindly contact:

JEULIN – S.A.V.

468 Jacques Monod Street
CS 21900
27019 EVREUX CEDEX France

+33 2 32 29 40 23