Ultrasons

Ultrasounds

**Ref :
223 010**

Français – p 1

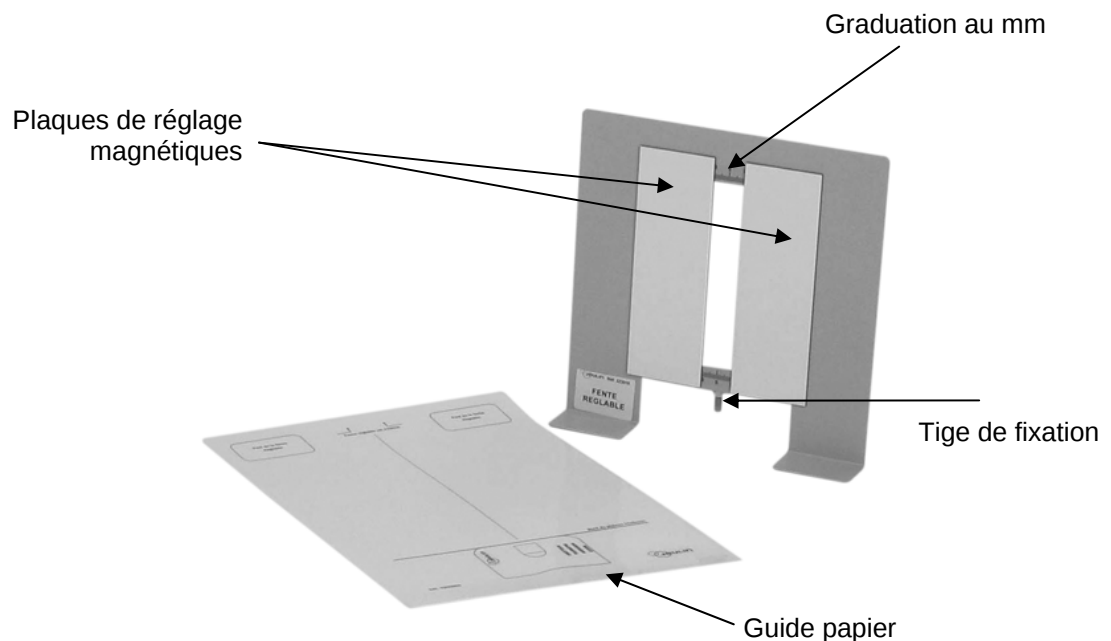
English – p 3

Version : 9003

Fente de diffraction réglable

Adjustable diffraction slit

1 Description

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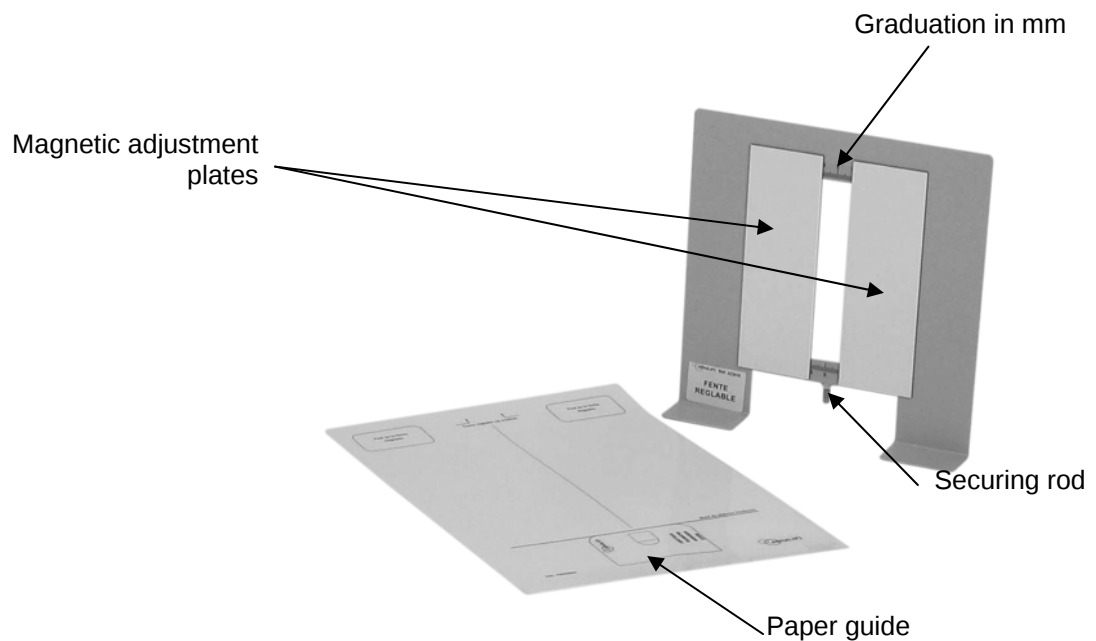
1 Description

Device used to view the minima and maxima of the diffraction pattern of ultrasonic waves: the width of the 40 x 120 mm slit can be adjusted from 0 to 40 mm by 2 magnetic plates.

The transmitter and the slit/plate assembly are placed on a plastic coated paper guide (to be fixed on the table) in perfect alignment.

Accessories required for the experiments (not provided with the device):

- | | |
|-----------------------|-------------------------|
| • Moduson Transmitter | Part no. 222 036 |
| • Moduson Receiver | Part no. 222 028 |
| • Moduson Plate | Part no. 223 008 |
| • Power supply R15+/- | |

2 Technical characteristics

- | | |
|----------------------|----------------------------|
| • Slit width: | adjustable from 0 to 40 mm |
| • Dimensions: | 111 x 222 x 333 mm |
| • Guarantee: | 2 years |

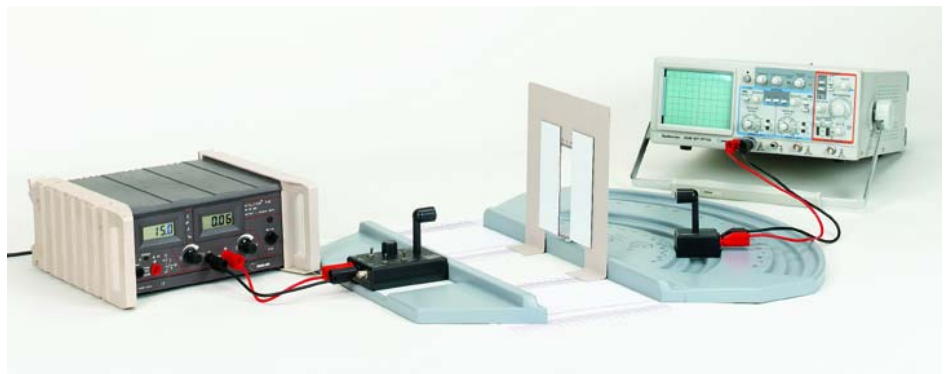
3 Experiment

❖ Demonstrating the diffraction of ultrasound

Fix the paper guide on a table then place the components as shown in the photograph below. The transmitter is powered by a voltage of 15V and constantly emits ultrasonic waves.

On the other side of the slit, the receiver moves on the graduated Moduson plate and sends the intensity of the signal received in function of the angle to an oscilloscope. We can thus construct the diffraction pattern for a determined slit width.

The 2 self-adhesive plates are used to vary the width of the slit and show its influence on the diffraction pattern.

Example of results

Width a (cm) of the slit	Angle θ (°) of the first minimum on the left	Angle θ (°) of the first minimum on the right	Theoretical angle θ (°)
2	- 25	24	25
3	- 18	18	16.5
4	-12	14	12.5

4 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
0825 563 563

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries