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## *Material properties*

Réf :  
251 009

Français – p 1

English – p 3

Version : 9101

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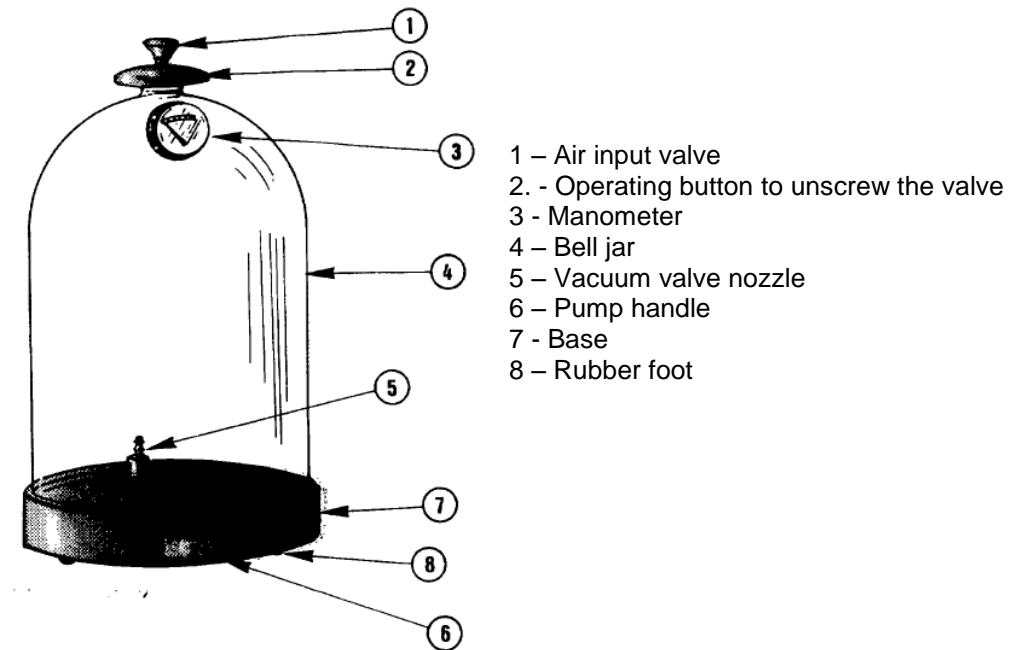
**0 825 563 563\***

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The device consists of a practically unbreakable transparent plastic vacuum chamber bell jar firmly fitted with a vacuum gauge and a dual shut-off valve. The bell jar is placed on a white plastic mounting plate consisting of a built-in pump and a retention valve.

The airtight joint of the bell jar to the base is done with a rubber ring inserted in a groove on the base. The bell jar can be used in complete safety as soon as it is placed on the rubber ring (do not use seal grease).

At the top of the bell jar is a double action valve. The flush-mounted red button can be pulled to let in small amounts of air into the bell jar. By unscrewing the valve completely, the large opening is freed for connecting other devices and instruments using its sleeve.

The vacuum valve located on the mounting plate is covered with a brass nozzle which, once the bell jar is removed, can receive a rubber tube that can be connected either to an independent mounting plate (Magdebourg hemispheres experiment) or a device that needs to be vacuumized (Dutch tears).

In this case the air input is obtained by unscrewing the nozzle slightly. To remove the tube it is recommended that the nozzle be completely unscrewed and the rubber be removed to avoid applying extra stress on the plastic threading.

The vacuum gauge consists of a relative pressure graduated scale of 0 to - 600 mm Hg. For example, if the air from the device is evacuated till the red line marking - 250, the pressure in the bell jar will be about  $760^* - 250 = 510$  mm Hg i.e. 0.67 atmosphere.

Caution: There is a risk of implosion if the vacuum level is exceeded or if the bell jar is dropped after creating a vacuum.

**\*average atmospheric pressure.**

**Additional accessories:**

- **Vacuum tube Part no.: 253 009**

## **After-Sales Service**

The device is under a 2-year guarantee, it must be sent back to our workshops.  
For any repairs, adjustments or spare parts please contact:

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## Direct connection for technical support

A team of experts  
at your disposal  
from Monday to Friday  
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request  
immediatly to provide you  
with the right answers regarding  
your activity field : Biology, Physics,  
Chemistry, Technology.

**Free service\***

**+33 2 32 29 40 50\*\***

\* Call cost not included.

\*\* Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux