Mesure

Initio® 2

Measure

Initio® 2

Réf :
251 180

Français – p 1

English – p 7

Version : 7105

Thermomètre Initio 2

Initio 2 Thermometer

1. Description

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0 825 563 563*

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
1. Description

The Initio 2 thermometer is an instrument that measures the temperature recorded using a Pt1000 probe. It can also be used to record data, store it and then restore it later on a computer.

A WIFI module (ref. 451 470) is also available for visualizing and saving measurements on a tablet or a PC equipped with WIFI.

Equipped with an analogue output, the Initio 2 thermometer can be used with a datalogging interface with direct inputs via banana jacks or via a Groomy® interface and jack.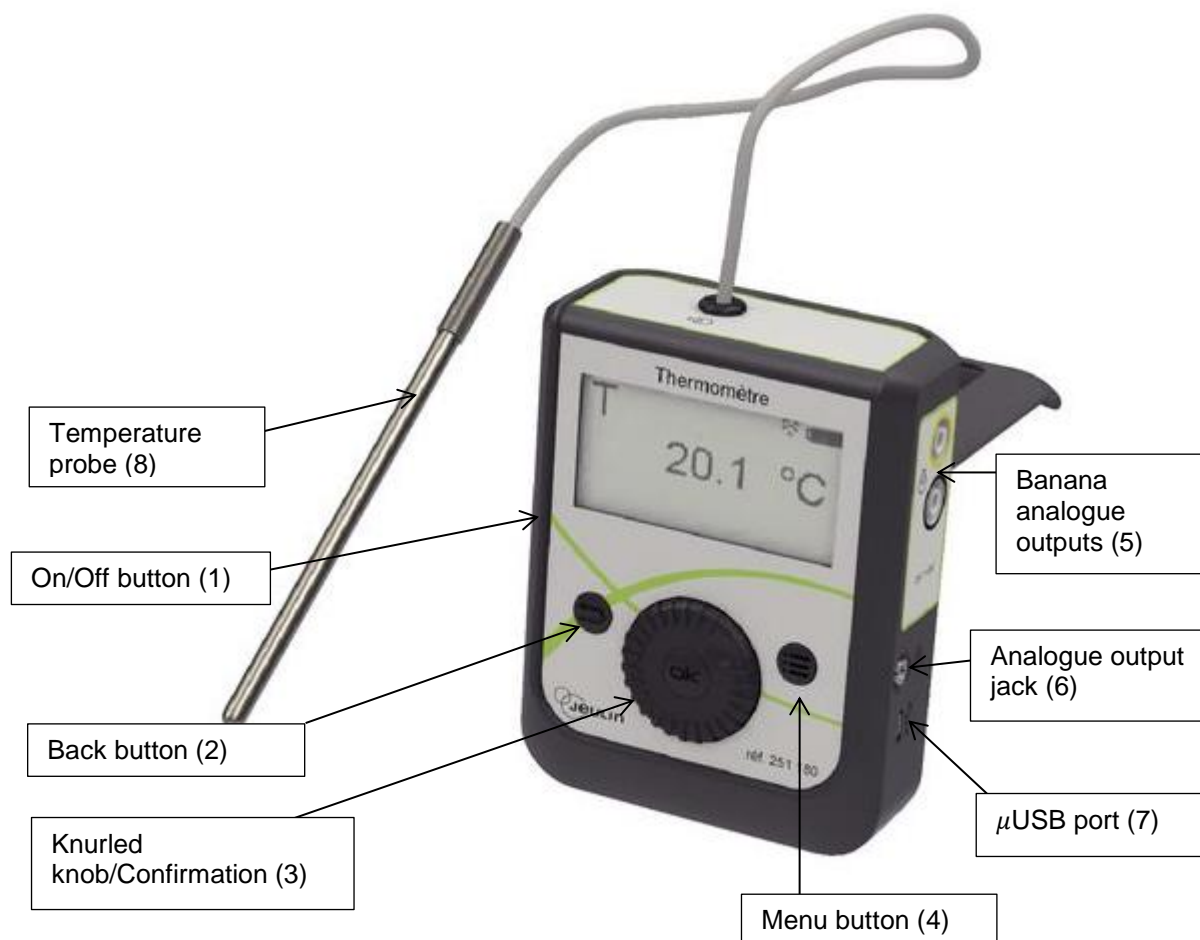

2. Utilisation

Upon first use, we recommend that the battery of the Initio 2 to be fully charged. For charging use the μ USB (7) charger supplied in separate packaging. In general, we recommend using the instrument in complete charge-discharge cycles (wait until it is completely discharged before fully recharging)

2.1 Putting into use

- Press the start button (1) situated on the left side of the instrument

The measurement displays directly on the instrument screen.

- To switch off the Initio 2, briefly press the start button (1) and confirm by pressing the knurled knob (3).

2.2 Navigation

Initio 2 has three navigation buttons:

Selectionconfirmation (3):

The central knurled knob (3) is for moving between the menus of the Initio 2. Briefly pressing the OK button in the centre of the knurled knob confirms the selection.

Back (2):

The "Back" button (2) makes it possible to change from one type of display to another from the main screen in the following order:

- Measurement
- Table
- Chart

Within a menu it is used to go back a step.

Menu (4):

This button displays the main menu of the Initio 2.

2.3 Data acquisition

The Initio 2 has an internal memory which allows it to save measurements. Two modes are available:

- Manual acquisition

Acquisition takes place on a point by point basis. Each measurement is manually saved by the user (see below)

- Time acquisition

Data acquisition takes place versus (duration entered by the user)

2.3.1 Manual acquisition

- Press the Menu button (4)
- Using the knurled knob (3), select **Acquisition** and confirm by pressing on the knurled knob (3).
- Select **Manual**
- To save a measurement, enframe the **Save** icon (white circle) and confirm by briefly pressing the knurled knob (3).
- To stop the acquisition enframe the **Stop** icon (black square) and confirm by pressing the knurled knob (3). Confirm the choice in the dialogue box.

2.3.2 Time acquisition

- Press the button Menu (4).
- Using the knurled knob (3), select **Acquisition** and confirm by pressing on the knurled knob (3)
- Select **Time**
- Using the knurled knob (3), enframe the number defining the duration and press on the knurled knob (3).
- Modify this duration by turning the knurled knob (3) then confirm by pressing on the centre of the knurled knob.
- If the data acquisition is to loop (return of the acquisition to zero at the end of a defined period) select **Loop**
- Confirm the final choice by selecting **OK**

Time data acquisitions last for a time between 60 seconds and 50 hours. The interval between two measurements adjusts automatically.

2.3.3 Recording the data

Once the acquisition is completed (time or manual) it must be saved to be able to retrieve the points on a computer.

- Press the Menu button (4)
- Select **File** and confirm
- Select **Save**

Only one data acquisition can be saved. By selecting **Open**, you open the last saved data acquisition.

2.3.4 Transfer of the data to the computer

Once the acquisition is completed, it is possible to transfer the data to a computer to process it:

- Install the Initio 2 software available under the address [www.jeulin.com/Initio 2 for an English version](http://www.jeulin.com/Initio_2_for_an_English_version)
- Connect the Initio 2 to the computer via the USB- μ USB cable
- Launch the Initio 2 software
- Load the saved acquisition by clicking on the icon:

The data then appears on the screen. Right-clicking the chart provides access to various chart tools.

3. Data visualisation

It is possible to view the data at any moment in tabular or chart format. To do so, simply press the Back button (2). The display then toggles between measurement, table and chart. Whatever the current display, the data acquisition can be monitored.

After 5 seconds without activity, table or chart is output as a full screen display. To again obtain the associated functions, turn the knurled knob.

In the table, the arrows situated at the bottom right are used to move between the rows of the table.

4. Analogue outputs

The Initio 2 has analogue outputs in the form of banana jacks and a 3.5 mm jack.


4.1 Banana jacks

- Connect the Initio 2 to an Ex.A.O. interface equipped with direct inputs using banana leads
- Configure the direct input in the following manner:
 - o -20 °C corresponds to -5 V
 - o 120 °C corresponds to +5 V


Example with the Foxy console

- Start the generalist

- Place the icon for a direct input  on the desired axis
- Select the range $\pm 5\text{ V}$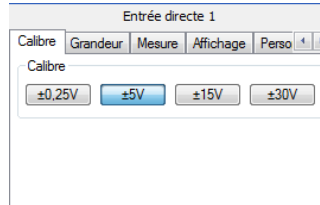

- Fill out the customization tab as below: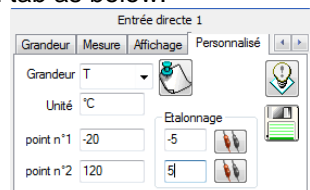

- Click on the diskette to save the configuration 

Configuration is complete, the value displayed is henceforth the temperature with symbol T and the units are °C.

4.2 Jack

The 3.5 mm jack is for connection of the Initio 2 to a Groomy interface (ref. 181 000) for example. The voltage supplied varies between 0 and 5 V. So that:

- -20 °C corresponds to 0 V
- 120 °C corresponds to 5 V

5. WIFI communication

Using the WIFI (ref. 451 470) module, it is possible to visualize the Initio 2 data on a tablet or similar.

5.1 Installation of the WIFI module

The WIFI module is fitted with the Initio 2 switched off.

- Raise the handle of the Initio 2
- Remove the back by sliding it down
- Position the WIFI module instead of the back
- Push the module firmly into place until it clicks. It must be possible to retract the handle without forcing it.

5.2 Visualisation on a tablet

Once the WIFI module is installed, the tablet must be configured:

- Press the Menu button
- Select Parameters
- Select Network

Information concerning the module is output to the screen.

- On the tablet, chose the name of the WIFI network displayed on the Initio 2

- Open a web browser
- Enter the address displayed by the Initio 2 in the address bar of the browser

6. Parameters

Various information concerning the Initio can be obtained by pressing on Menu and selecting Parameter → Information

This page displays the serial number of the instrument and the software version, the battery charge level and the operating language. The latter is changed by turning the knurled knob.

7. Update

On our website at the address <http://en.jeulin.fr/en-resources/english-resources/software/free-software.html>, the last update for your Initio® 2 is available. Its version is indicated in the designation.

In order to know the version of your Initio® 2, press the “Menu” button and then “Setup” “Informations”.

8. Technical characteristics

Temperature	-20 to +120 °C
Resolution	0.1 °C
Precision	±1 %
Probe	fixed aluminum Ø 6 x 120 mm with 1.6 m cable (no naked flames)
Analogue output	-5 V to +5 V via 4 mm safe banana jacks 0 V to +5 V via 3.5 mm jack
Communication	USB, WIFI (optional extra)
Power supply	Li-ion battery
Dimensions	95 x 123 x 44 mm

9. After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.
For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
468 rue Jacques Monod
CS 21900
27019 EVREUX CEDEX FRANCE

+33 (0)2 32 29 40 50