Mesure

Initio 2

Measure

Initio 2

Réf :
251 182

Français – p 1

English – p 7

Version : 7105

Oxymètre Initio 2
Initio 2 oxymeter

1 Description

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0 825 563 563*

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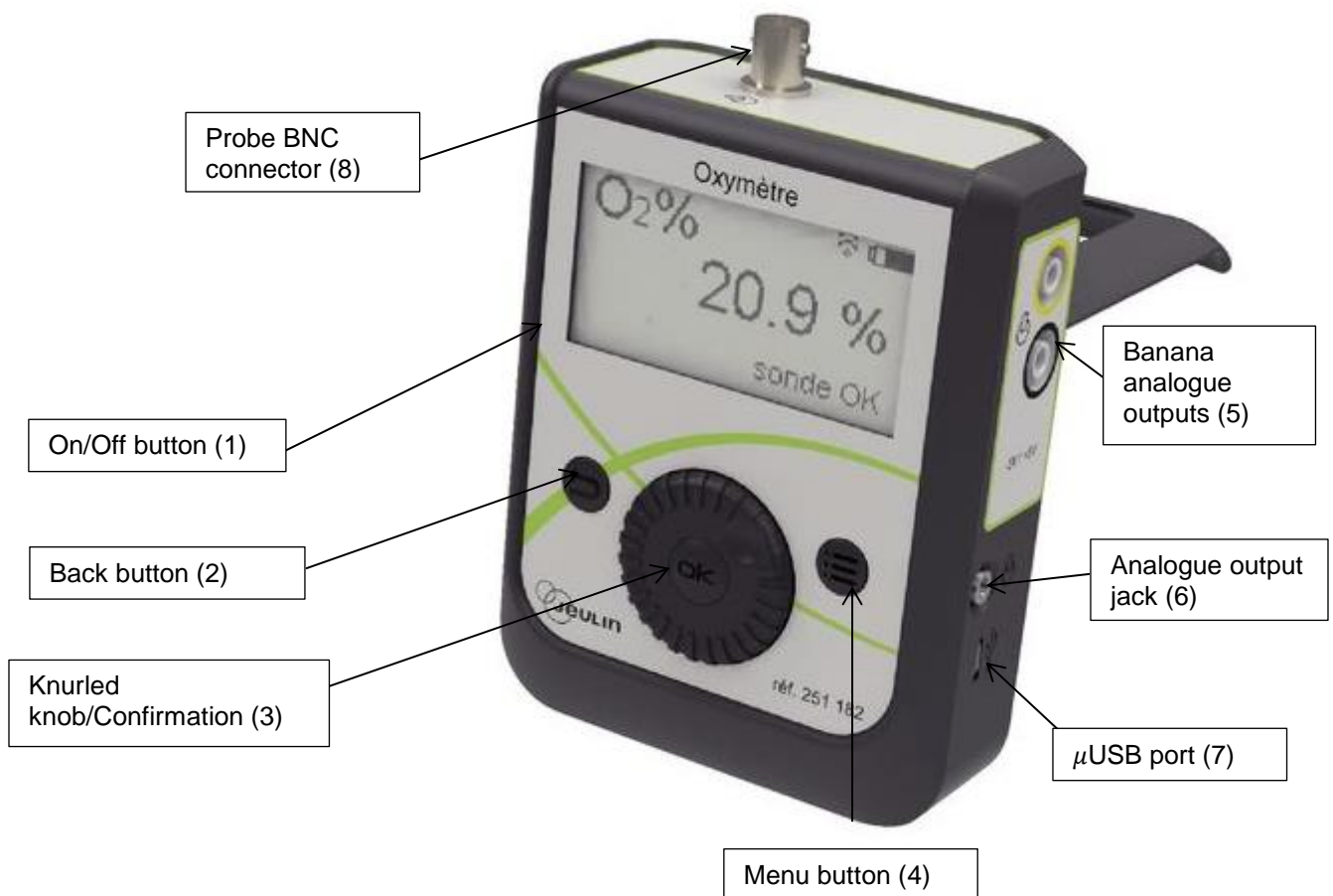
1 Description

The Initio 2 (a mixture of Initio 2 and Initio 2 is used in the source) oxygen meter is an instrument that measures the concentration of oxygen in air and water using an oxymetric probe, reference 453 052. It can also be used to record data, store it and then restore it later on a computer.

A WIFI module (ref 451 470) is also available for visualising and saving measurements on a tablet or a PC equipped with WIFI.

Equipped with an analogue output, the Initio 2 oxygen meter can be used with a datalogging interface with direct inputs via banana jacks or via a Groomy® interface and jack.

2 Utilisation

Upon first use, we recommend the battery of the Initio 2 to be fully charged. For charging use the µUSB (7) charger supplied in separate packaging. In general, we recommend using the instrument in complete charge-discharge cycles (wait until it is completely discharged before fully recharging).

2.1 Navigation

Initio 2 has three navigation buttons :

Selection-confirmation (3):

The central knurled knob (3) is for moving between the menus of the Initio 2. Briefly pressing the OK button in the centre of the knurled knob confirms the selection.

Back (2):

The "Back" button (2) makes it possible to change from one type of display to another from the main screen in the following order:

- Measurement
- Table
- Chart

Within a menu it is used to go back a step.

Menu (4):

This button displays the main menu of the Initio 2.

2.2 Putting into use

- Connect the probe (ref 453 052) to the Initio 2 via the BNC connector (8)
- Press the start button (1) situated on the left side of the instrument

The measurement displays directly on the instrument screen.

- To switch off the Initio 2, briefly press the start button (1) and confirm by pressing the knurled knob (3).

2.3 Calibration

- Press the menu button (4)
- Select Calibration
- Using the knurled knob, indicate the level of molecular oxygen in the surrounding air (for example 20.9%)
- Press on the knurled knob to validate.

2.4 Type of measurement

Using the probe ref 453 052, it is possible to measure the concentration of molecular oxygen in air or water. To switch from one function to the next

- From the measuring screen, press on the knurled knob
- Turn the knob to change from "Air" to "Water"
- Press on the knurled knob to confirm

2.5 Data acquisition

The Initio 2 has an internal memory which allows it to save measurements. Two modes are available:

- Manual acquisition

Acquisition takes place on a point by point basis. Each measurement is manually saved by the user (see below)

- Time acquisition

Data acquisition takes place versus time (duration entered by the user)

2.5.1 Manual acquisition

- Press the Menu button (4)
- Using the knurled knob (3), select **Acquisition** and confirm by pressing on the knurled knob (3)
- Select **Manual**
- To save a measurement, enframe the **Save** icon (white circle) and confirm by briefly pressing the knurled knob (3)
- To stop the acquisition enframe the **Stop** icon (black square) and confirm by pressing the knurled knob (3). Confirm the choice in the dialogue box

2.5.2 Time acquisition

- Press the button Menu (4).
- Using the knurled knob (3), select **Acquisition** and confirm by pressing on the knurled knob (3)
- Select **Time**
- Using the knurled knob (3), enframe the number defining the duration and press on the knurled knob (3)
- Modify this duration by turning the knurled knob (3) then confirm by pressing on the centre of the knurled knob
- If the data acquisition is to loop (return of the acquisition to zero at the end of a defined period) select **Loop**
- Confirm the final choice by selecting **OK**

Time data acquisitions last for a time between 60 seconds and 50 hours. The interval between two measurements adjusts automatically.

2.5.3 Recording the data

Once the acquisition is completed (time or manual) it must be saved to be able to retrieve the points on a computer.

- Press the Menu button (4)
- Select **File** and confirm
- Select **Save**

Only one data acquisition can be saved. By selecting **Open**, you open the last saved data acquisition.

2.5.4 Transfer of the data to the computer

Once the acquisition is completed, it is possible to transfer the data to a computer to process it:

- Install the Initio 2 software available under the address [www.jeulin.com/Initio 2](http://www.jeulin.com/Initio_2)
- Connect the Initio 2 to the computer via the USB- μ USB cable
- Launch the Initio 2 software
- Load the saved acquisition by clicking on the icon:

The data then appears on the screen. Right-clicking the chart provides access to various chart tools.

3 Data visualisation

It is possible to view the data at any moment in tabular or chart format. To do so, simply press the Back button (2). The display then toggles between measurement, table and chart. Whatever the current display, the data acquisition can be monitored.

After 5 seconds without activity, table or chart is output as a full screen display. To again obtain the associated functions, turn the knurled knob.

In the table, the arrows situated at the bottom right are used to move between the rows of the table.

4 Analogue outputs

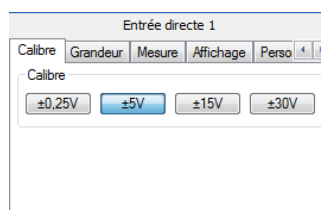
The Initio 2 has analogue outputs in the form of banana jacks and a 3.5 mm jack.

4.1 Banana jacks

- Connect the Initio 2 to a datalogging interface equipped with direct inputs using banana leads
- Configure the direct input in the following manner:
 - o In air
 - 0 % corresponds to -5V
 - 25 % corresponds to +5V
 - o In water
 - 0 mg/L corresponds to -5V
 - 20 mg/L corresponds to +5V

Example using the Foxy® console for an air measurement

- Start the generalist
- Place the icon for a direct input  on the desired axis
- Select the range $\pm 5V$

- Fill out the customization tab as below:

- Click on the diskette to save the configuration 

Configuration is complete, the value displayed is henceforth the concentration of oxygen with symbol [O2] and the unit is mg/L.

4.2 Jack

The 3.5mm jack is for connection of the Initio 2 to a Groomy interface (ref 181 000) for example. The voltage supplied varies between 0 and 5V. So that:

In air

- 0 % corresponds to 0V
- 25% corresponds to 5V

In water

- 0 mg/L corresponds to 0V
- 20 mg/L corresponds to 5V

5 WIFI communication

Using the WIFI (451 470) module, it is possible to visualize the Initio 2 data on a tablet or similar.

5.1 Installation of the WIFI module

The WIFI module is fitted with the Initio 2 switched off.

- Raise the handle of the Initio 2
- Remove the back by sliding it down
- Position the WIFI module instead of the back
- Push the module firmly into place until it clicks. It must be possible to retract the handle without forcing it.

5.2 Visualization on a tablet

Once the WIFI module is installed, the tablet must be configured:

- Press the Menu button
- Select Parameters
- Select Network

Information concerning the module is output to the screen.

- On the tablet, chose the name of the WIFI network displayed on the Initio 2.
- Open a web browser
- Enter the address displayed by the Initio 2 in the address bar of the browser.

6 Parameters

Various information concerning the Initio can be obtained by pressing on Menu and selecting Parameter → Information

This page displays the serial number of the instrument and the software version, the battery charge level and the operating language. The latter is modified by turning the knurled knob.

7 Update

On our website at the address <http://en.jeulin.fr/en-resources/english-resources/software/free-software.html>, the last update for your Initio® 2 is available. Its version is indicated in the designation.

In order to know the version of your Initio® 2, press the "Menu" button and then "Setup" "Informations".

8 Technical characteristics

Measurement in air

Measurement range 0 to 20 %

Resolution 0.1 %

Precision ± 1 %

Measurement in water

Measurement range 0 à 20 mg/L

Resolution 0.1 mg/L

Precision ± 1 %

Probe Ref 453 052 (Air/Water Probe)
Ref 453051 (Air Probe)

Analogue output -5 V to +5 V via 4 mm safe banana jacks
0 to +5 V via 3.5 mm jack

Communication USB, WIFI (optional extra)

Power supply Li-ion battery

Dimensions 95 x 123 x 44 mm

9 After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.
For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
468 rue Jacques Monod
CS 21900
27019 EVREUX CEDEX FRANCE

+33 (0)2 32 29 40 50