Electricité

Magnétisme

Electricity

Magnetism

Ref :
252 044

Français – p 1

English – p 3

Version : 9002

Spectre magnétique

Magnetic spectrum

1 Généralités

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0 825 563 563 *

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1 General information

Aesthetic and easy to use: no more iron filings to be handled. The perfect assistant to demonstrate magnetic field lines in 1, 2 or 3 dimensions. Is also suitable for demonstrations and experiments.

2 Use and composition

Several unique experiments may be performed with this new magnetic field demonstration kit.

The kit consists of 4 plastic plates containing a total of 376 magnetic slugs of 8 mm length. Each plate measures 153 x 77 x 6 mm.

The plates are transparent, which enables the experiments to be viewed with an overhead projector. The kit can be set up at your discretion by using the plates separately or in groups of 2, 3 or 4 in order to cover a larger area.

For demonstrations in 2 or 3 dimensions, you only need to put the necessary number end to end.

3 Possible experiments

- The plates are piled up and all 376 magnetic slugs represent elementary magnets in a demagnetised piece of steel. A magnetised bar is placed on top of the plates and the magnetic slugs position themselves along a unidirectional field with the North and South poles.
- Test with a magnetic needle. The "magnet" gets demagnetised if it is shaken or tapped and the magnetic slugs (elementary magnets) position themselves randomly.

4 Additional equipment

- | | |
|-----------------------------------|------------------|
| • Ticonal magnets | Part no. 263 007 |
| • Ferrite magnets | Part no. 263 016 |
| • Magnetised needle on agate yoke | Part no. 263 009 |

5 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries