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***Measurement and instrumentation***

**Réf :  
281 001**

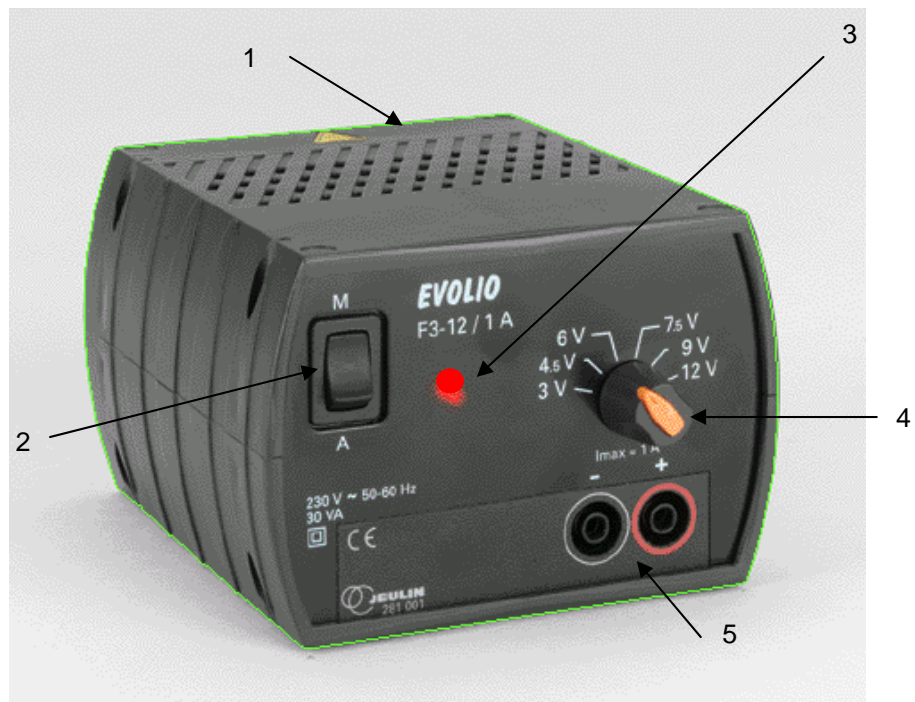
**Français – p 1**

**English – p 4**

**Version : 8010**

**Alimentation EVOLIO F3-12 / 1A**  
***EVOLIO F3-12 / 1A power supply***

## 1 Description

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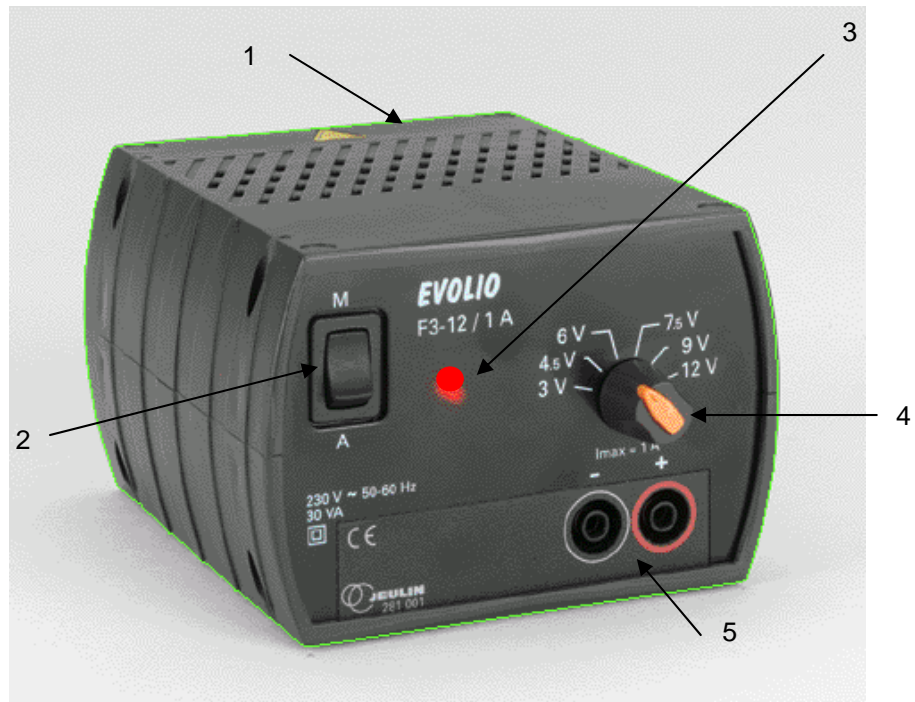
## 1 Description

- 1 – Plastic box with double insulation
- 2 – On/off switch
- 3 – Operating indicator
- 4 – 3 / 4.5 / 6 / 7.5 / 9 / 12 voltages switch
- 5 – CD outputs on safety sockets

## 2 Setting up

- Connect the power supply to the 230 V mains supply.
- Select the desired voltage by toggling the switch 3 / 4.5 / 6 / 7.5 / 9 / 12V to the corresponding position.
- Connect your circuit (lamp, motor, voltage divider, etc.) to the output sockets.
- Toggle the master switch so that the power-on indicator lamp comes on.

### Notes

If the current flow on the output used is greater than the maximum permissible value (1 A), the current remains at about 1 A and the voltage gradually reduces. Remove the source of the overload or the short circuit to return to normal operation.

## 3 Precautions for use

- Avoid using the power supply in direct sunlight or close to a heat source.

- Never block the ventilation vents when the power supply is operating.
- If this device is used in a manner not mentioned in these instructions, the safety ensured by the device may be compromised.

## 4 Technical characteristics

- **Voltage:** 3 / 4.5 / 6 / 7.5 / 9 / 12V fixed, DC
- **Outputs:** Safety sockets  $\varnothing$  4 mm
- **Maximum current:** 1 A
- **Stabilisation:** 1 %
- **Residual ripple:** Less than 10 mV
- **Power supply:** 230 V  $\pm$  10 %, 50-60 Hz
- **Power consumed**  
**from the mains:** 30 VA
- **Protection:** Primary: Auto-reset thermal protection in the transformer.  
Secondary: electronic
- **Dimensions:** 115 x 140 x 80 mm
- **Mass:** 950 g
- **Guarantee:** 2 years

*This device complies with:*

- *The low Voltage directive 73/23/CEE; applied standard: EN 61010-1*
- *The CEM 89/336/CEE directive; applied standards: EN 50081-1 and EN50082-1*

### 4.1 Environmental conditions for use of the device

- For indoor use
- Altitude less than 2000 meters
- Operating temperature from 15°C to 40°C
- Maximum relative humidity of 80% up to 31°C and 50% up to 40°C
- Mains supply variations of +/-10%
- Installation category II, pollution level 2

### 4.2 Sectioning

It is the power plug that is used as a sectioning device.

### 4.3 Symbols used

: Caution, refer to accompanying documents.

: Caution, hot surface.

## 5 Care and maintenance

In case of a malfunction:

- Check the mains voltage (230 V)
- Check the open circuit voltages, without connecting any load

**Note:** In case of use in high ambient temperatures, the thermal protection located in the primary of the transformer may get activated in exceptional cases. In this case, disconnect the power supply and wait for 1 hour so that the device cools down.

If the device still does not work, contact the After Sales Service.

In all cases, **do not open the device**.

Store the original packaging; it can be used to bring the device to our workshops for maintenance, if necessary.

## 6 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

**JEULIN - TECHNICAL SUPPORT**  
**Rue Jacques Monod**  
**BP 1900**  
**27 019 EVREUX CEDEX FRANCE**  
**0825 563 563**

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

## Direct connection for technical support

A team of experts at your  
disposal from Monday  
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

### Free service \*

**+ 33 (0)2 32 29 40 50\*\***

\* Call cost not included

\*\* Only for call from foreign countries