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## *Measurement and instrumentation*

Ref :  
281 002

Français – p 1

English – p 5

**Alimentation EVOLIO F15 ± / 0,4 A**  
***EVOLIO F15 ± / 0.4 A power supply***

## 1 Description

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## NOTES

## 1 Description

- 1 – Plastic box with double insulation
- 2 – On/off switch
- 3 – Operating indicator
- 4 – -15, 0, +15 V outputs on safety sockets

## 2 Setting up

- Connect the power supply to the 230 V mains supply.
- Toggle the master switch so that the power-on indicator lamp comes on.
- Connect your circuit (lamp, motor, voltage divider, etc.) to the selected output sockets.

### Notes

If the current flow on the output used is greater than the maximum permissible value (0.4 A), the current remains at about 0.4 A and the voltage gradually reduces. Remove the source of the overload or the short circuit to return to normal operation.

### 3 Precautions for use

- Avoid using the power supply in direct sunlight or close to a heat source.
- Never block the ventilation vents when the power supply is operating.
- If this device is used in a manner not mentioned in these instructions, the safety ensured by the device may be compromised.

### 4 Technical characteristics

• <b>Voltage:</b>	- 15, 0, + 15 V fixed, DC
• <b>Outputs:</b>	Safety sockets ø 4 mm
• <b>Maximum current:</b>	0.4 A at each output
• <b>Stabilisation:</b>	1 %
• <b>Residual ripple:</b>	Less than 10 mV
• <b>Power supply:</b>	230 V ± 10 %, 50-60 Hz
• <b>Power consumed from the mains:</b>	30 VA
• <b>Protection:</b>	Primary: fuse calibrated on I.C. Secondary: electronic
• <b>Dimensions:</b>	115 x 140 x 80 mm
• <b>Mass:</b>	950 g
• <b>Guarantee:</b>	2 years

This device complies with:

- The low Voltage directive 73/23/CEE; applied standard: EN 61010-1
- The CEM 89/336/CEE directive; applied standards: EN 50081-1 and EN50082-1

#### 4.1 Environmental conditions for use of the device

- For indoor use
- Altitude less than 2000 meters
- Operating temperature from 15°C to 40°C
- Maximum relative humidity of 80% up to 31°C and 50% up to 40°C
- Mains supply variations of +/-10%
- Installation category II, pollution level 2

## 4.2 Sectioning

It is the power plug that is used as a sectioning device.

## 4.3 Symbols used

: Caution, refer to accompanying documents.

: Caution, hot surface.

## 5 Care and maintenance

In case of a malfunction:

- Check the mains voltage (230 V)
- Check the open circuit voltages, without connecting any load

If the device still does not work, contact the After Sales Service.

In all cases, **do not open the device**. Replacing the mains fuse must be done at our workshops.

Store the original packaging; it can be used to bring the device to our workshops for maintenance, if necessary.

## 6 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

**JEULIN - TECHNICAL SUPPORT**  
**Rue Jacques Monod**  
**BP 1900**  
**27 019 EVREUX CEDEX FRANCE**  
**0825 563 563**

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# Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediately to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

## Free service \*

**+ 33 (0)2 32 29 40 50\*\***

\* Call cost not included

\*\* Only for call from foreign countries