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## *Test and Measurement*

Réf :  
**281 404**

Français – p 1

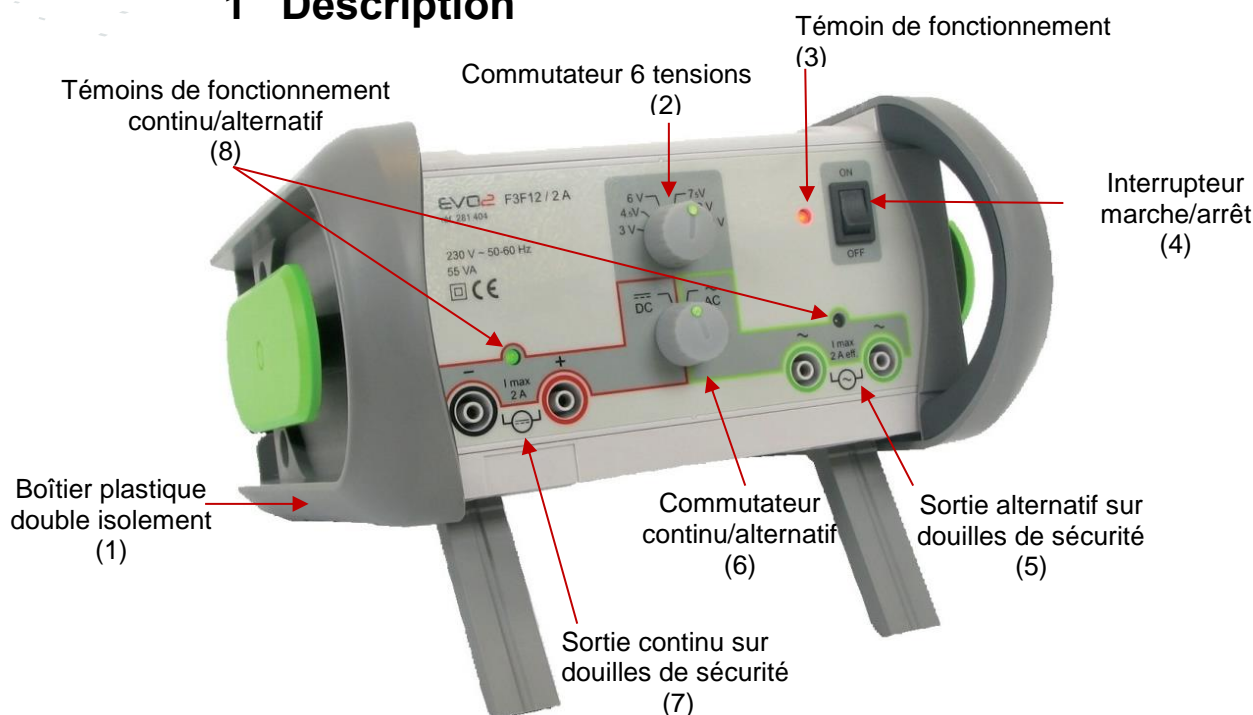
English – p 4

Version : 5106

**Alimentation EVO2<sup>®</sup> F3F12 / 2 A**

***EVO2<sup>®</sup> F3F12 / 2 A power supply***

## 1 Description

## 2 Mise en œuvre

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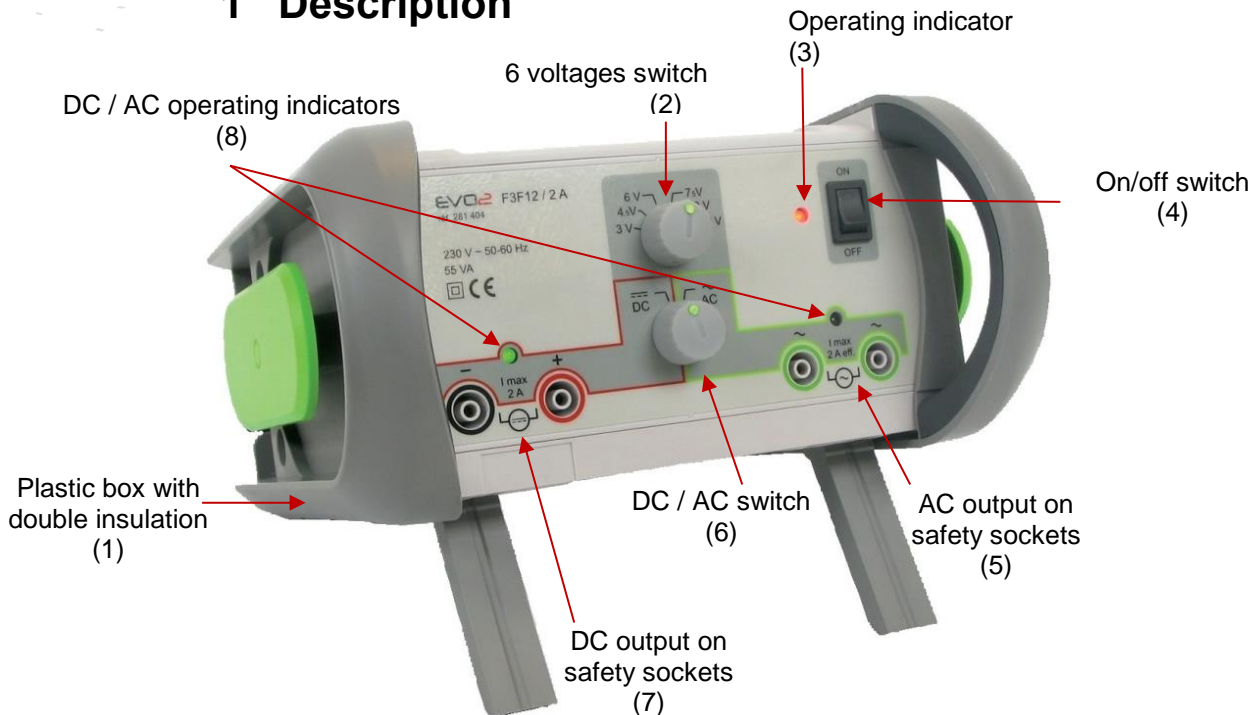
## 1 Description

## 2 Setting up

- Connect the power supply to the 230 V mains supply.
- Select the desired voltage by turning the switch 3 / 4.5 / 6 / 7.5 / 9 / 12V to the corresponding position (2).
- Select the type of current by turning the AC / DC switch (6). The lamp (8) corresponding to the activated output comes on.
- Connect your circuit (lamp, motor, voltage divider, etc.) to the selected output sockets.
- Toggle the master switch (4) to "On". The power-on indicator lamp (3) lights on.
- 

### Notes

In **DC**, if the current flow is greater than 2 A, the power supply trips. The resetting must be done manually by switching off the power supply for about fifteen seconds. Switch on the power supply after ensuring that the cause of the over-current is eliminated.

In **AC**, if the current flow is greater than 2 A RMS, the power supply trips. This protection is ensured by a thermal protector, which resets automatically after a few minutes once the source of the overload, or the short circuit is removed.

In **AC**, the power supply delivers a maximum voltage that fluctuates close to 12 V in function of the load of the electrical circuit and the mains voltage (between 11.5 V and 14 V).

## 3 Precautions for use

- Avoid using the power supply in direct sunlight or close to a heat source.
- Never block the ventilation vents when the power supply is operating.
- If this device is used in a manner not mentioned in these instructions, the safety ensured by the device may be compromised.

## 4 Technical characteristics

- **Voltage:** 3 / 4.5 / 6 / 7.5 / 9 / 12 V fixed, DC or AC
- **Outputs:** Safety sockets Ø 4 mm
- **Maximum current:** 2 A at each output
- **Stabilisation:** 3 %
- **Residual ripple:** less than 10 mV
- **Protection:** Primary: T250 mA / 250 V fuse, CEI127 approved, accessible from the back of the device,  
Secondary: Electronic switch-off in DC, Thermal with auto-reset in AC
- **Power supply:** 230 V  $\pm$  10 %, 50-60 Hz
- **Power consumed from the mains:** 55 VA
- **Dimensions:** 235 x 235 x 125 mm
- **Mass:** 2 kg
- **Warranty:** 2 years

This device complies with:

- The low Voltage directive 2006/95/CE ; applied standard : EN 61010-1:2010
- The CEM 2004/108/CEE directive ; applied standards : EN 60326-1

### 4.1 Environmental conditions for use of the device

- For indoor use,
- Altitude less than 2000 meters,
- Operating temperature from 15 °C to 40 °C,
- Maximum relative humidity of 80 % up to 31 °C and 50 % up to 40 °C,
- Mains supply variations of  $\pm$ 10 %,
- Installation category II, pollution level 2.

## 5 Care and maintenance

In case of a malfunction:

- Check the mains voltage (230 V),
- Check the open circuit voltages, without connecting any load,
- Wait 2 to 3 minutes for the protection against over-currents to be turned on again.

### 5.1 Replacing the fuse

Remove the fuse holder located at the rear face of the power supply. Install only a T250 mA / 250 V slow-blow fuse Ø 5 x 20 mm.

If the device still does not work, contact the After Sales Service.

In all cases, **do not open the device.**

Store the original packaging, it can be used to bring the device to our workshops for maintenance, if necessary.

## **6 After-Sales Service**

The device is under a 2-year guarantee, it must be sent back to our workshops.  
For any repairs, adjustments or spare parts please contact:

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## Direct connection for technical support

A team of experts  
at your disposal  
from Monday to Friday  
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request  
immediatly to provide you  
with the right answers regarding  
your activity field : Biology, Physics,  
Chemistry, Technology.

### Free service\*

**+33 2 32 29 40 50\*\***

\* Call cost not included.  
\*\* Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux