Electricité

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Electricity

Three-phase sinusoidal mode

**Ref :
302 189**

Français – p 1

English – p 4

Version : 8009

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1 Asynchronous three-phase motor

1.1 Objective

- To measure the power consumed by an asynchronous three-phase motor in a star-delta connection.
- To measure a frequency of rotation.
- To calculate the sliding $g = \frac{n_s - n}{n_s}$.

1.2 Equipment

- An extra-low voltage three-phase generator;
- An extra-low voltage three-phase asynchronous motor;
- A double-pole change-over switch;
- A set of digital multimeters;
- An extra-low voltage wattmeter WJ2000;
- Connecting wires;
- A stroboscope.

1.3 Experiment

- ✓ Make the star-delta connection (for this refer to the simulation performed on the extra- low voltage three phase terminal board).
- ✓ Adjust the extra-low voltage three-phase generator: $3 \times 9 \text{ V} - 50 \text{ Hz}$.
- ✓ Observe the rotational movement of the rotor. Interchange the two phase wires; observe the change in direction of the rotation of the rotor.
- ✓ Measure for each connection:

- Line current I_Y or I_Δ
- The voltage
- The total power consumed P_Y or P_Δ
- The rotational speed of the rotor n_Y or n_Δ
- The sliding g_Y or g_Δ .

For measuring the power, refer to the simulation performed on the extra-low voltage three-phase terminal board.

To measure the rotational speed n of the rotor using the stroboscope: increase the frequency of flashes; the rotational speed corresponds to the largest frequency for which the reference frame is immobile.

Check that the star circuit is balanced: $I_N \approx 0$ A.

Check that the motor rotates slower in star connection.

Verify the relations: $\frac{I_\Delta}{I_Y} = 3$ and $\frac{P_\Delta}{P_Y} = 3$.

Compare P measured with the value of P given by the relation:

$$P = U \times I \times \sqrt{3} \times \cos \varphi.$$

Examples of results:

$$n_s = \frac{f}{p} = \frac{50}{2} = 25 \text{ rps}$$

$$g = \frac{n_s - n}{n_s}$$

n_s : rotational frequency of the rotating field

f : Frequency of the current

p : number of pole pairs

g : sliding

n_s : rotational frequency of the rotating field

n : rotational frequency of the rotor

Connection	I measured	U measured	P measured	n measured	g
Star	78 mA	15.6 V	1.06 W	1378 rpm 23 rps	13%
Delta	171 mA	15.6 V	*	1554 rpm 25.9 rps	2%

* The extra-low voltage wattmeter WJ2000 enables measurements only between the neutral and phase. (star connection).

$$\cos \varphi = \frac{P}{S} \quad \cos \varphi = \frac{1,06}{15,6 \times 0,078 \times \sqrt{3}} \approx 0.50$$

2 Single phase asynchronous motor

2.1 Objective

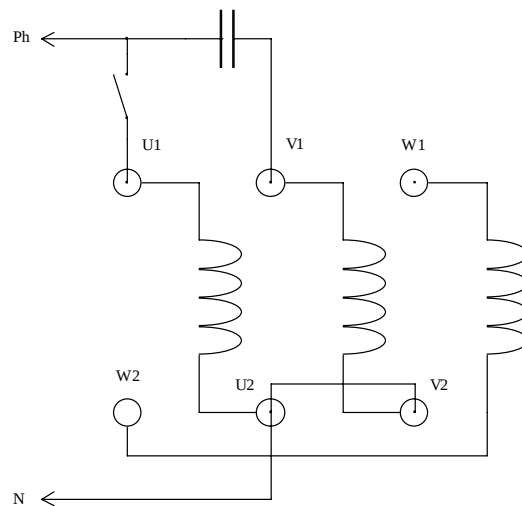
To measure the power consumed by a single-phase asynchronous motor.

2.2 Equipent

- An extra-low voltage A.C. supply
- An extra-low voltage three-phase asynchronous motor
- A 33 μF un-biased capacitor
- Two digital multimeters
- An extra-low voltage wattmeter WJ2000
- Connecting wires.

2.3 Experiment

Carry out the set-up as shown below

Two coils are set out of phase using a capacitor.

(Put on the switch; observe the direction of rotation of the rotor. Once the motor has started, remove the capacitor.

Measure the supply voltage, the line current and the power consumed; compare the power measured using the digital wattmeter to the value of the power given by the relation:

$$P = U \times I \times \cos \varphi .$$

3 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
0825 563 563

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries