# Mécanique

Statique

# *Mechanics*

*Statics*

Ref :  
322 008

Français – p 1

English – p 3

Version : 9007

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## 1 Contents

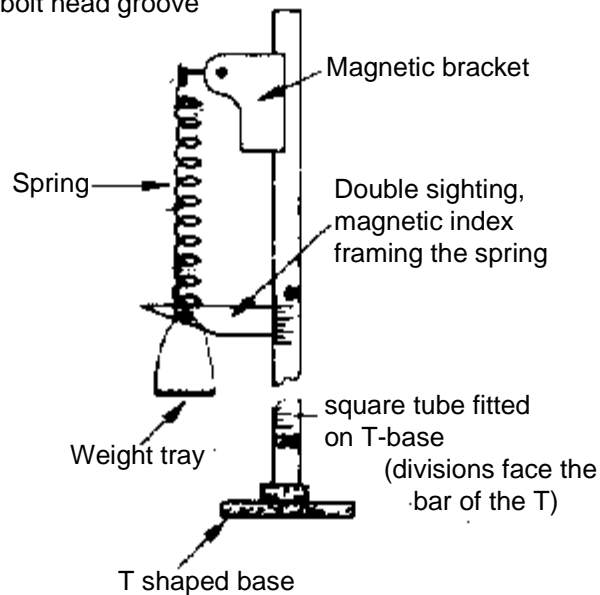
The apparatus consists of:

- 1 frame consisting of a T shaped base, highly stable, and a square, nickel steel tube having a printed scale on one side. For reasons of size, the two components are not pre-assembled.
- 1 double sighting index enabling balance points to be marked.
- 1 movable magnetic support, used as a base for attaching springs.
- 1 set of 3 springs, Ref. 323 006.
- 1 tray to take weights, Ref. 323 005.

The two last components can be purchased separately.

## 2 Assembly


Spring hooked to  
 bolt head groove

## 3 Practical

Principle of the sight.

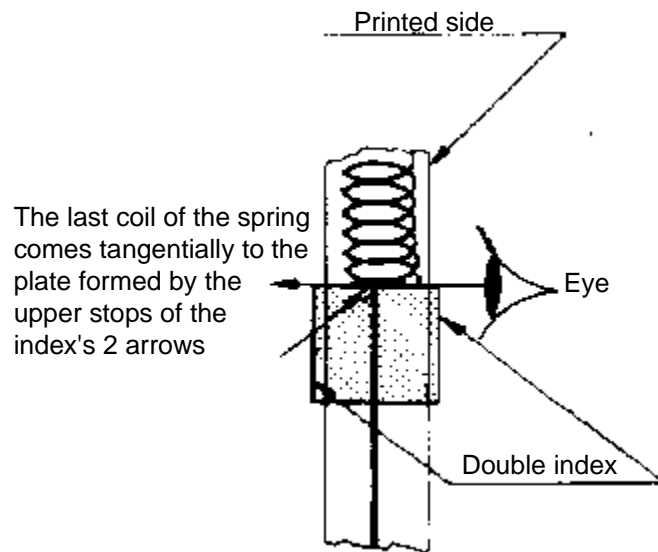The double sighting index eliminates errors due to parallax.

### 3.1 Zeroing

With the tray attached to the spring, slide the bracket until it aligns, as shown above, the index coinciding with the zero on the scale.

### 3.2 Comment

To increase or decrease the distance between spring and frame, adjust the lock bolt.

## 4 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts please contact:

**JEULIN - TECHNICAL SUPPORT**  
**Rue Jacques Monod**  
**BP 1900**  
**27 019 EVREUX CEDEX FRANCE**  
**+33 (0)2 32 29 40 50**

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

## Direct connection for technical support

A team of experts at your  
disposal from Monday  
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

### Free service \*

**+ 33 (0)2 32 29 40 50\*\***

\* Call cost not included

\*\* Only for call from foreign countries