Mécanique

Statique

Mechanics

Statics

Ref :
322 046


Français – p 1

English – p 5

Version : 6010

Plan incliné statique
Static incline

1 Principe et description

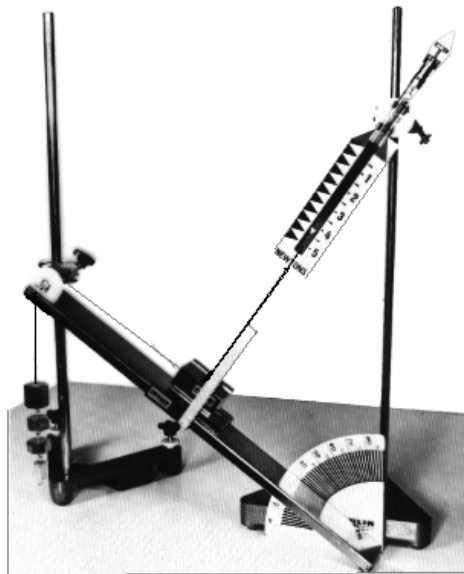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
1 Principle and description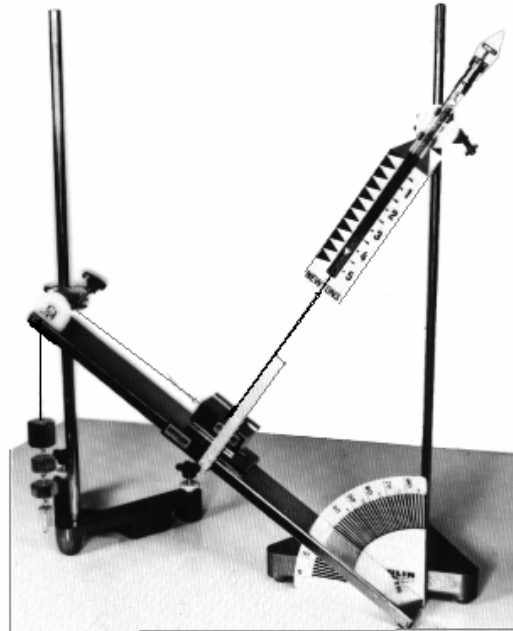

Simple set-up comprising 1 incline, 1 teaching dynamometer, 2 stands with rods, 2 fastening fixtures, 1 box of masses with hooks.


This apparatus is used to measure the forces applied to a moving object at equilibrium on an incline: useful component (parallel to the plane) and normal plane reaction.

This apparatus has been specifically designed for use both with hooked masses and with “teaching” or “lab work” dynamometers.

A further advantage is the possibility to demonstrate the plane reaction when measuring R . The plane can then be subtracted without suppressing the equilibrium.

The apparatus permits to :

- verify the relation :
 $F = P \sin \alpha$
- measure the normal plane reaction :
 $R = P \cos \alpha$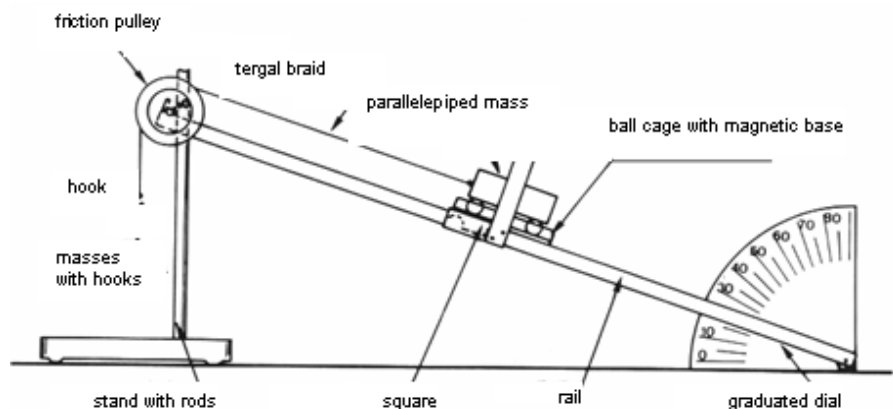

2 Composition

The apparatus comprises:

- a steel rail, 600 x 60 mm with a fixture for mounting a pulley;
- a (captive) ball cage on a magnetic base;
- a graduated sector dial for fitting on the rail;
- a sliding square for checking the wire perpendicularity during R measurements;
- a parallelepipedic carriage fitted with 2 hooks;
- two pulleys Ø. 50 mm mounted on a rod;
- two hooks and tergal braid for experiments with hooked masses.

3 Additional accessories required

The following additional accessories are necessary to perform the two above-mentioned experiment types:

<i>Designation</i>	<i>Quantity</i>	<i>Part no</i>
Universal rod stand	2	701031 or 701030
Fastening fixture	2	703 099
Hooked masses	1	703 016
Lab dynamometer	1	322 050
Teaching dynamometer	1	322 013

4 Set up

First mount the sliding square on the rail by inserting it from one end of the rail (notch side); the spring engages the lower side of the rail. Mount a pulley on the device provided for that purpose.

The spindle of this pulley is also used for maintaining the incline in the desired position.

Prepare a universal rod stand with fastening fixture. The pulley spindle engages the free groove in the fastening fixture.

Finish by mounting the graduated dial (insert the banana plug into the rivet located at the rear on the right of side of the rail) and install the ball cage on the rail.


5 Experiments

5.1 Measurement of the useful component

In this special case the square is not needed.

The experiment consists in establishing the equilibrium of the parallelepipedic mass under the action of the hooked masses by adjusting the value thereof.

For the 3 balls to be centered in their housing, it is best to proceed in two steps: first allow the balls to roll to the housing bottom then place the parallelepiped and raise it (see drawing below). Then use the braid to raise the parallelepiped until the balls are centered. If the equilibrium is correct, the mass should remain stationary.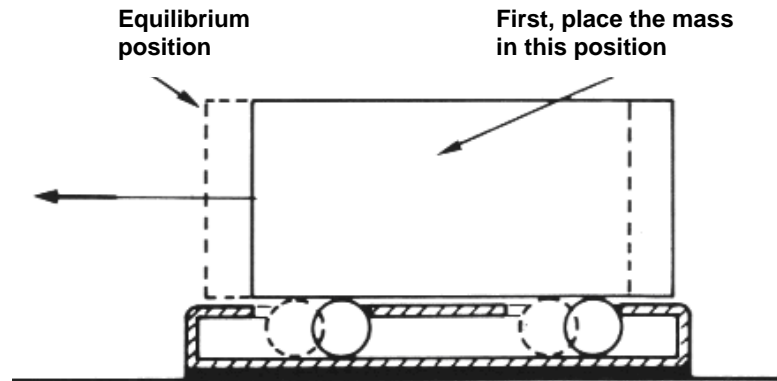

Let α the angle of the rail relative to the horizontal plane (as read on the graduated dial), F the mass hung on the wire (the hook mass shall be taken into account as required) and P the mass of the parallelepiped (weighed beforehand).

Then the following relationship is verified to within the error margin:

$$F = P \cdot \sin \alpha$$

5.2 Measurement of the normal plane reaction


The principle consists in replacing the plane reaction by an equal force applied through a wire tied onto the upper side of the parallelepiped.

At equilibrium, the action of the rail on the mass is zero; the removal of the rail must not alter the mass position, which can be achieved with this apparatus by loosening the screw used to unite the pulley and the incline (without contacting the fastening fixture) then by lowering the rail.

To measure the force R , it is first necessary to mount the second pulley on a second universal rod stand by using a fastening fixture.

The exact mass is determined by adding masses until the first ball starts to roll (it is heard meeting the extremity of its housing).

Then the following relationship is verified: $R = P \cos \alpha$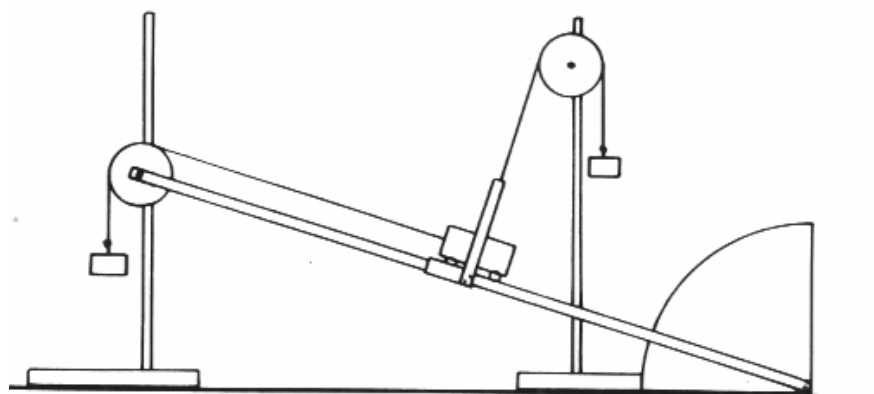


5.3 Set up with dynamometers

The experiment principle is the same. The forces F and R are measured by using either “teaching” dynamometers or “lab work” dynamometers.

In this case, the pulleys are no longer needed but their spindle will be used as a mounting means (for that, the pulleys should be reversed, i.e. mounted at the rear of the fastening fixture).

The dynamometers are installed between the parallelepiped hooks and the two spindles. The force is adjusted by moving the fastening fixture and the parallelepiped farther away or closer, respectively.

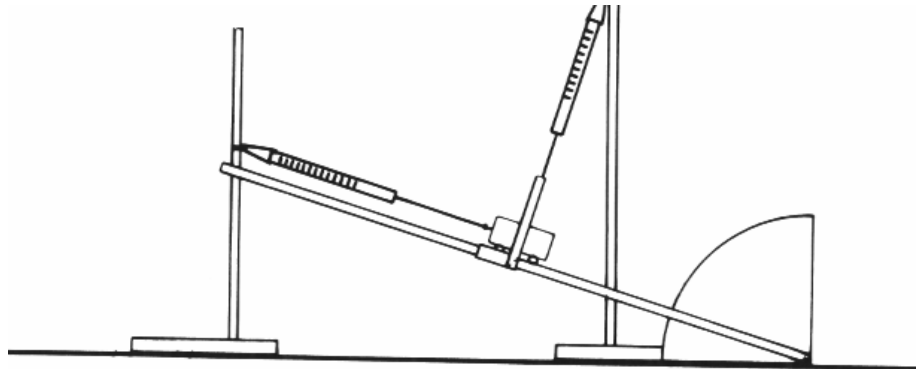6 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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Service gratuit * :
+ 33 (0)2 32 29 40 50

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Rubrique FAQ

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

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