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## *Mechanical and Optic*

Réf :  
**323 050 – 323 060**

Français – p 1

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**0 825 563 563\***

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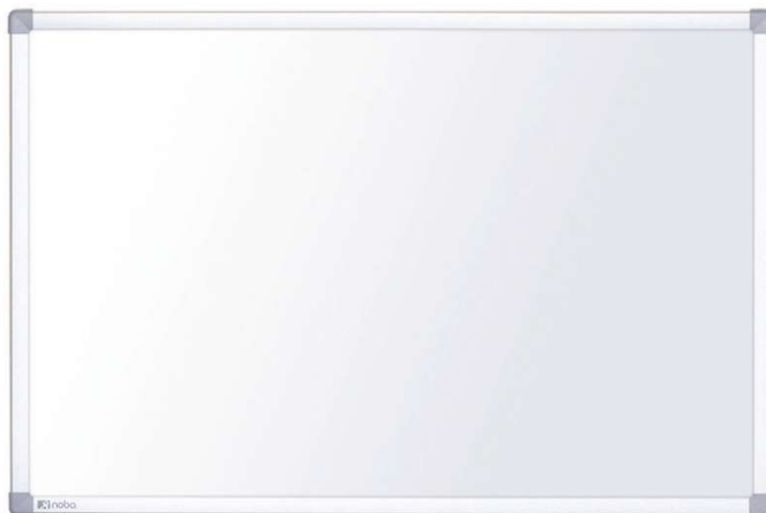
## 1. Description

This metal panel allows you to carry out your mechanical or optical experiments which may require fixing elements with a magnetic base. It can be positioned horizontally or vertically depending on your need. The rewritable surface offers the possibility of writing with markers and erasing with the supplied wipe. The board is also supplied with a wall mounting system and all its accessories for more conventional use.

## 2. Composition

1 rewritable metal board 600 x 400 mm or 900 x 600 mm :

Accessories supplied (wall brackets, holder for felt-tip pens, felt-tip pen, microfiber cloth):

2 Modumontage feet :

2 Polynux tightening nuts :

2 Modumontage rods length 400 mm :


### 3. Mounting

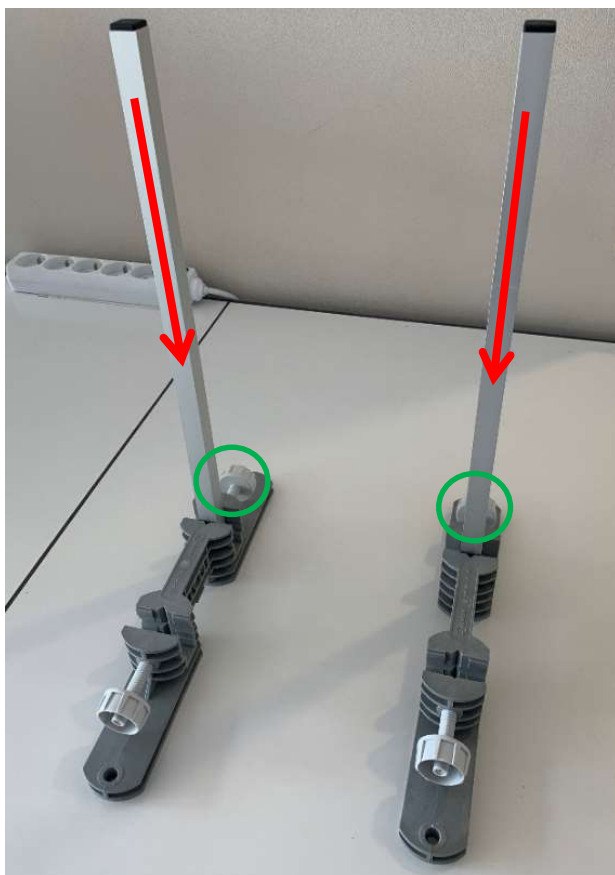In order to install the table on its support, please follow the following assembly steps:  
**Step 1:** place the 2 Modumontage feet on a flat surface

**Step 2:** Place the 2 Modumontage rods in the 2 notches provided, having previously unscrewed the 2 knurled screws circled on the visual:

**Step 3:** Screw the 2 knurled screws on each foot in order to secure the rod with the foot

**Step 4:** Position the 2 Polynux nuts in this way on each foot respecting the directions of insertion indicated by the arrows and allowing a height of around 15 cm between the upper part of the foot and the nut

**Step 5:** Tighten the 2 Modumontage screws on each rod to make the whole united

**Step 6:** Place the table as below

**Step 7:** Tighten the 2 Modumontage screws on the panel face

## **4. After-sales service**

The device is under a 2-year guarantee, it must be sent back to our workshops.  
For any repairs, adjustments or spare parts please contact:

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## Direct connection for technical support

A team of experts  
at your disposal  
from Monday to Friday  
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request  
immediatly to provide you  
with the right answers regarding  
your activity field : Biology, Physics,  
Chemistry, Technology.

### Free service\*

**+33 2 32 29 40 50\*\***

\* Call cost not included.

\*\* Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux