Mécanique

Dynamique

Mechanics

Dynamics

Réf :
332 056

Français – p 1

English – p 7

Version : 7112

Mobile jet d'encre – Pack rénovation

Mobile Inkjet – Refit

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0 825 563 563*

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1 Description

This assembly is adapted for the study of two-dimensional movements.
A horizontal plane support is necessary for the study of movements.

If you already have a Jeulin glass table (references 332010 and 332021) and accessories such as accessories for shocks, launcher, etc. then this offer is ideal for you. Accessories for old spark versions are compatible with your refit.

1.1 Composition and installation

This pack is composed of:

- 2 air cushion pucks
- 2 accumulators 9 V-300 mAh
- 1 power supply 12 V-500 mA
- 1 "Y" cable
- 2 humidifying bases
- 1 remote control
- 1 edge marking accessory
- 1 accessory to search for the centre of inertia of a system composed of 2 connected pucks


1.1.1 Air cushion pucks

Ink cartridges and batteries supplied are not installed in pucks at the time of delivery.

The user should install these elements in pucks as described below:

1/ Open the puck (2 screws)

2/ Open the sealing cover concealing the ink cartridge and the pump

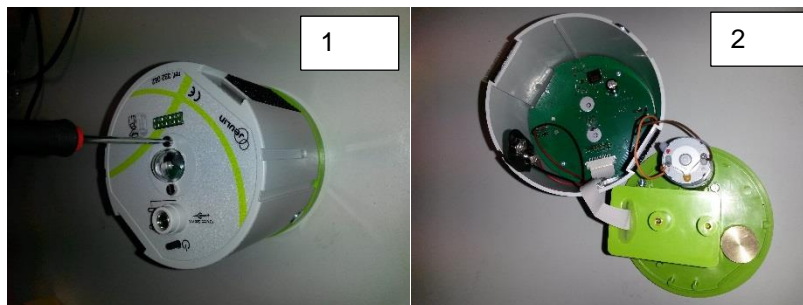Warning: the removal of the hood should be no sudden move to avoid breaking flounder

3/ Put the ink cartridge into place (see picture): take care that that it click fits properly in its housing

Close the sealing cover

Caution: Position the seal in its notch before closing the cover to maintain a perfect seal. In the case where the seal would be poorly positioned, the air cushion would not be strong enough to ensure satisfactory lift.

4/ Connect the battery and put it into its housing (between the 3 studs)

5/ Close the cover taking care not to wind the ink cartridge ribbon cable. Make use of the guide slots when closing the puck.

6/ Press the « On/Off » button and check that the LED comes on. To change LED color, press the “on/off” button during 3 seconds.

1.1.2 Humidifier base

2 humidifier bases are provided to protect ink cartridges installed in pucks.

Air cushion puck on humidifier base

It is **ESSENTIAL** that the puck should be put back on this base after each manipulation. This will prevent the ink cartridge from drying out.
 The sponge included in the base must be kept damp.

Do not touch the print head unless it is dry.
 In this case, apply isopropyl alcohol onto the print head carefully using a cotton swab.

If marking contains 2 lines for a puck, then the print head has been damaged. The print cartridge must be replaced.

1.1.3 Remote control

The time Δt between two successive markings is always constant so that the rate variation can be observed from the distance variation.

A display shows this Δt .

This remote control includes 3 buttons.

The central button switches the remote control on. For a long press, this button triggers printing of the Δt shown.

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The remote control switches off automatically after 5 minutes.

1.1.4 Markings external to the puck

1.1.4.1 Marking at the edge of the puck

Edge marking accessory

The edge marking accessory is composed of an ink cartridge support that is marked synchronously with the puck. The second element is a counterweight to balance the puck.

The various steps in assembly of this accessory are described below:

1/ Place an ink cartridge (to be ordered separately or to be taken from one of your pucks) in the ink cartridge support.

2/ Fix and connect the cartridge support onto the puck


3/ Place the counterweight on the puck facing the ink cartridge support.

Your puck is ready for synchronised marking at its centre and on its edge.

1.1.4.2 Marking at the centre of two connected pucks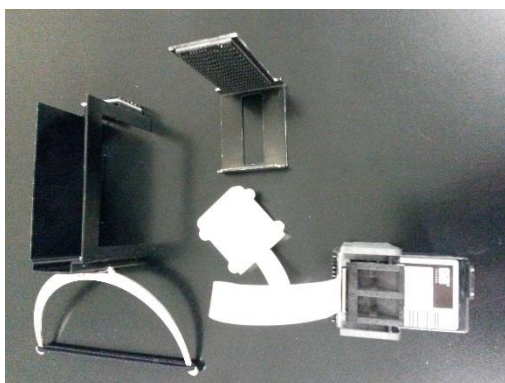


1 - Centre of inertia marking assembly (the ink cartridge should be taken from one of the pucks)

Step 1: Place your ink cartridge in the cartridge support


Step 2: Pass the connector into the window of the L with velcro tape then scratch the cartridge support in the L.

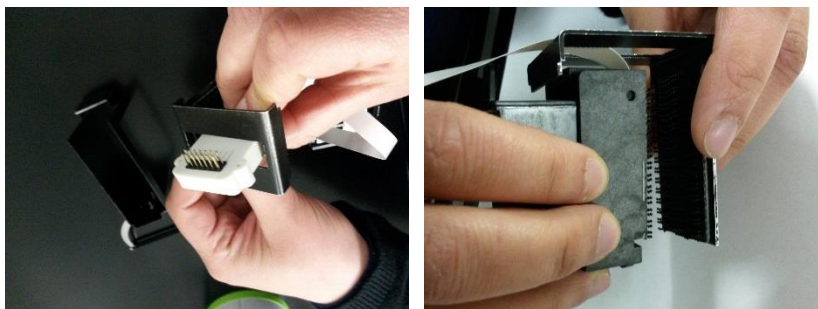
Step 3: Use velcro tape to fix the rectangle onto a puck and then connect the puck using the ring and the elastic provided on the diametrically opposite side.

Step 4: Slide the assembly with the ink cartridge into the rectangle.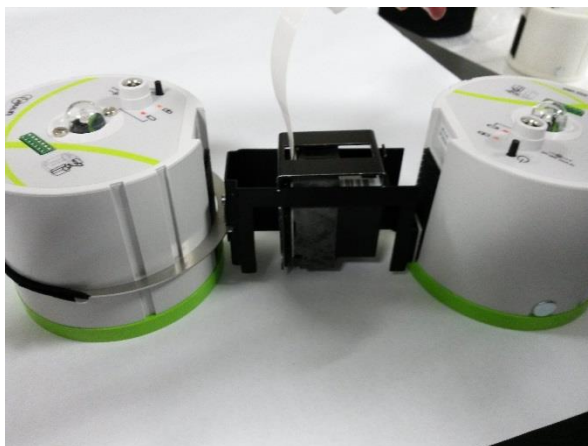

Step 5: Connect this marking accessory on top of the puck containing an ink cartridge (see picture below).

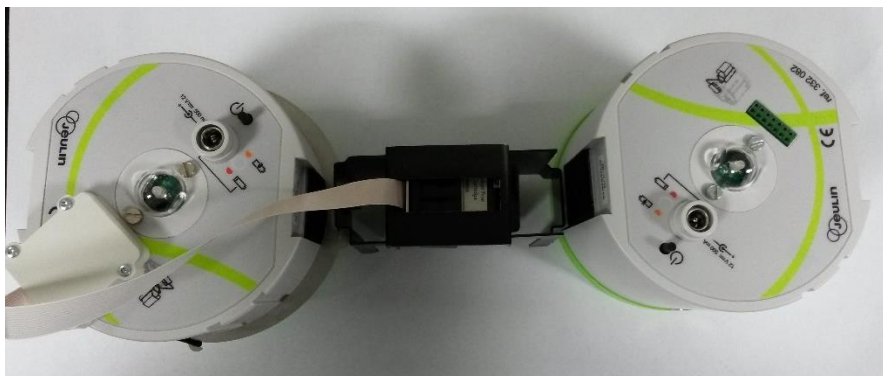Usage precautions: Take care to position the two pucks symmetrically about the central point of the set of two connected pucks. See picture above.

2 After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.
For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
468 rue Jacques Monod
CS 21900
27019 EVREUX CEDEX France

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Direct connection for technical support

A team of experts
at your disposal
from Monday to Friday
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request
immediatly to provide you
with the right answers regarding
your activity field : Biology, Physics,
Chemistry, Technology.

Free service*

+33 2 32 29 40 50**

* Call cost not included.
** Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux