Méthode ESAO®

ESAO® method

**Ref :
452 122**

Français – p 1

English – p 5

Version : 6005

Adaptateur Ethanol ESAO® 4

Ethanol adaptator ESAO® 4

1 Généralités

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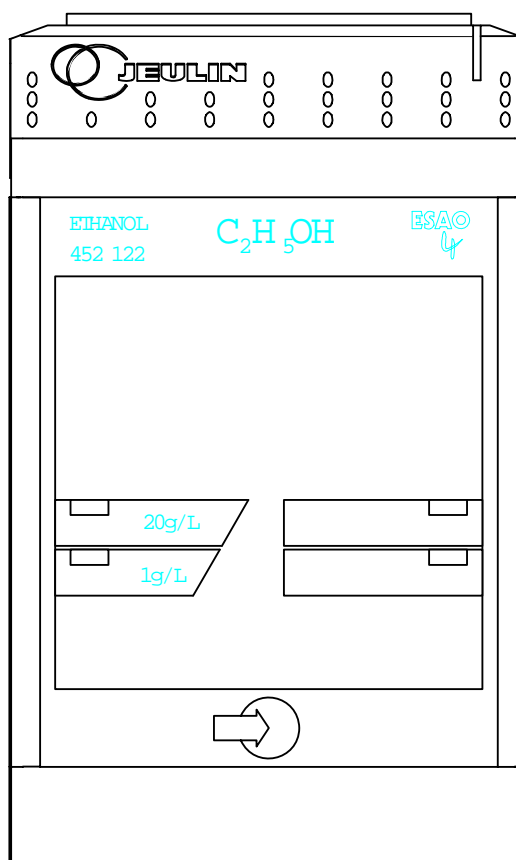
1 General information

This adaptor is designed to measure ethanol concentration using the ESAO method.

It is used with the ethanol probe: Part no.: 453 084.

Its function is to amplify and linearise the probe signal to adapt it to the input of the ESAO interface.

It enables the user to select the type of measurement (in air or in water) and to calibrate the probe using a potentiometer and a push-button.

2 Use

2.1 Connecting the adaptor

The ethanol adaptor is designed to operate with the ESAO system.

With ESAO 4:

- ESAO 4 console Part no. 451 101
- ESAO 4 interface card Part no. 451 300

Or

With ESAO 4+ console Part no. 451 401

The adaptor is connected using the connector located at the back of the box.
Follow the guide slots to connect it.
Engage it fully.

2.2 Connecting the ethanol probe

This adaptor is used with the ethanol probe part no. 453 084.
The probe is connected using the 5-pin DIN header located in front of the adaptor.

2.3 Choice of the type of measurement

On the adaptor:

symbolises measurement in air

symbolises measurement in water.

The choice of the measurement type is done by the software.

2.4 Commissioning

- Check that the ESAO card is correctly installed in the computer.
- Check that the console is connected to the ESAO card
- Connect the adaptor
- Connect the ETHANOL probe.
 - Switch on the computer.

2.5 Software packages

In addition to the software packages dedicated fully or partially to the measurement of ETHANOL concentrations and which therefore directly recognise it, this adaptor may be used with the general-purpose software:

Généris

Part no. 000 720

Configuring the software with the adaptor, configure according to the following model:

Adapters file edit

Nr: 38 Name: Ethanol
type: Esao 4 Quantity: Ethanol_rate

Type: SINGLE SENSOR ☒ Double function gas
liquid

Unit channel 1: %éthanol_air
Unit channel 2: g/L Option: User's calibration

Number of calibers: 2

Min: 0 0
Max: 2.5 0.125
0/2.5 % 0/0.125 %

Function 2
Min: 0 0
Max: 20 1
0/20 g/L 0/1 g/L

Save Current file: Physics secondary school Exit

2.6 Calibration of the ETHANOL probe

The calibration consists in adjusting the gain of the adaptor, the measuring system showing 0 at zero concentration.
The calibration solution may be prepared by diluting denaturated alcohol (or pure alcohol for better precision) in water.

Operating procedure:

It is mandatory to calibrate at the same temperature than the planned measurement and stir during this period.

With a general-purpose software (Generis or Serenis)

- Put the software in the adaptor test position on the ethanol adaptor channel
- Place the selector on the gas or liquid position depending on the planned experiment environment
- Immerse the electrode in the calibration solution
- Wait for 10 min. Stabilisation of the sensor
- Using the button, adjust the value read with the value in the table below.

1 ml of alcohol /l of water	0.8 g/l	0,1%
20 ml of alcohol /l of water	16 g/l	2%

2.7 Technical characteristics

Measuring ranges	in gases:	from 0 to 0.125 %
		from 0 to 2. 5 %
	in liquids	from 0 to 1 g/l
		from 0 to 20 g/l
Response time		<1 sec
Accuracy		10% of the entire scale
Mass		260 g
Dimensions		105 x 66 x 49 mm

3 After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.

For any repairs, adjustments or spare parts, please contact:

JEULIN – TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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Service gratuit * :
+ 33 (0)2 32 29 40 50

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Rubrique FAQ

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

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