Ex.A.O.

Accessoires respiration animale

Data logging

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0 825 563 563*

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1. General Information

This kit consists of two chambers 40 and 250 ml in volume, respectively, for performing experiments related to respiratory gas exchange.

The chambers are transparent and their volumes are small, thus they are ideally suited for studying respiration in small animals (mealworms, crickets, stick insects, etc.), plants, mushrooms,

Ideal for the study of animal respiration, they allow optimization of the result by amplifying the visualization of the magnitude studied, such as the drop in the O₂ concentration for example.

The various adapter rings make it possible to use the various probes available from Jeulin.

2. Contents

- One 250 ml chamber with a lid perforated with 3 holes:
 - 2 for gas exchanges with a diameter of 27 mm
 - 1 for the temperature of diameter 3 mm
 -
- One 40 ml chamber with a lid
- 2 adapter plugs
- 2 stoppers 12 and 3 mm in diameter

3. Instructions for use

3.1 The 250 ml chamber

- Unscrew the lid and then insert the O₂ and CO₂ probes in the holes provided for this purpose. The chambers are supplied with adapter plugs for adjusting the size of the holes to the diameter of the probe you are using (§ 3.3)
- Plug the holes not being used with the stoppers.
- Place the animal or animals in the chamber.
- Screw the lid back in place and start taking the readings immediately.

We recommend using the 40 ml chamber for experiments with plant material.

3.2 The 40 ml chamber

- Place the animal in the chamber
- Insert the adapter ring and place the probe therein
- Start taking the readings

This chamber can also be used to study gas exchange in plants. For this application, we recommend using two light sources in order to avoid shadow zones.

3.3 The adapter plug

The adapter cap is used to hold any type of probe with a diameter between 12 mm and 22 mm

The 12 mm diameter probes can be inserted directly.

For probes with a diameter greater than 12 mm, the cap must be cut to the desired diameter.

Proceed as follows : (The example below is provided for a 15.1 mm diameter probe.)

- 1) Insert the probe into the adapter plug

- 2) Locate the stop zone using your fingers

- 3) Cut out, using scissors, the stopper at the mark

- 4) Insert the probe into the cap to check that it is properly maintained

4. After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.
For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
468 rue Jacques Monod
CS 21900
27019 EVREUX CEDEX FRANCE

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Direct connection for technical support

A team of experts
at your disposal
from Monday to Friday
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request
immediatly to provide you
with the right answers regarding
your activity field : Biology, Physics,
Chemistry, Technology.

Free service*

+33 2 32 29 40 50**

* Call cost not included.

** Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux