# Méthode ESAO®

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## **ESAO® method**

*Stretch reflex study*

Ref :  
453 068

Français – p 1

English – p 4

Version : 9010

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## *Electronic Martoreflex*

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## 1. Overview

The Electronic Martoreflex is designed to function with all ESAO® JEULIN's interface.

It allows to :

- Cause a stretch reflex by hitting tendon (ankle or patellar reflex),
- Synchronize recording of the muscular response by the ESAO® system.

## 2. Description


- Steel hammer
  - Length : 230 mm (without cord)
  - Diameter of the head : 32 mm
  - Weight : 202 g
- Cable connexion : longueur 1,5 m. The cord is integrated to the Martoreflex and ends with 2 securized plugs, with a 4mm diameter use the interface connection.
- Piezoelectric integrated system trigerring signal.

## 3. Use

The electronic Martoreflex serves to synchronize the acquisition trigerring action when it hits the tendon, typically the achille's tendon.

To do this :

Plug the cord of the Electronic Martoreflex to the sync input of the ESAO® interface : the black plug into the black socket (with the mass symbol) the red plug on the sync socket with red brand.

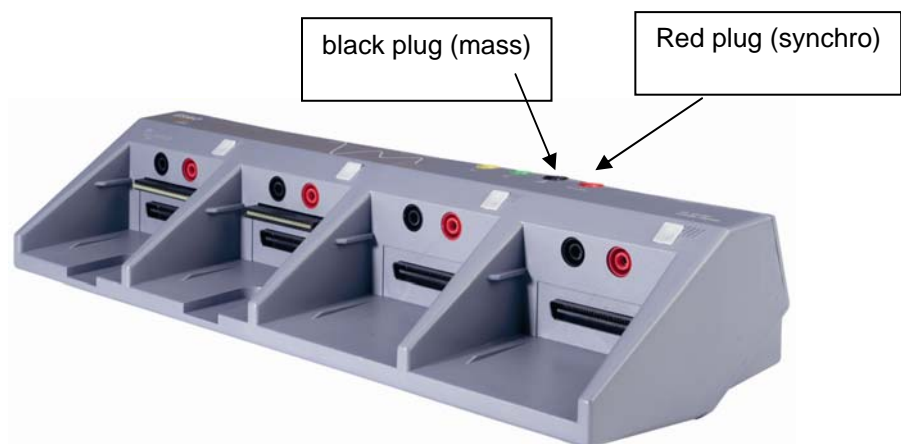
*Example : Terminal branching of electronic Martoreflex for ESAO® interface VISIO (réf. 451460) and Visio+ (réf. 451461).*

The response of muscle is typically recorded using an electrophy sensor.

## Software parameter settings


In the setup menu of the Scientific Workshop, external sync function must be activated to trigger the acquisition when the electronic Martoreflex hits the tendon.

To do this in the settings tab :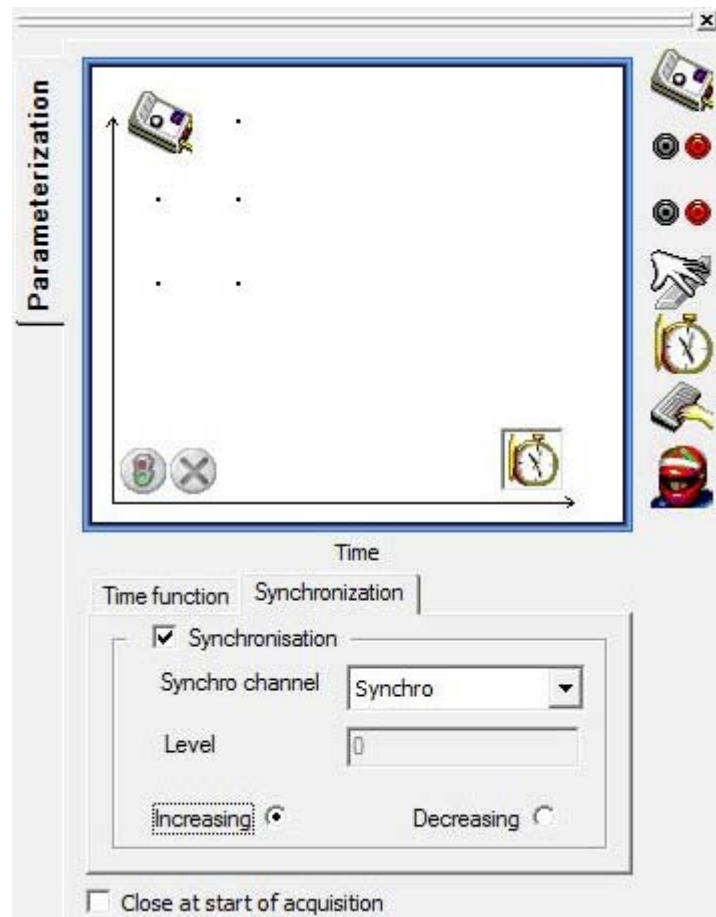

- Drag the icon of the electrophysiology sensor on vertical axis.
- Drag the icon of the timer in horizontal axis.
- Then click on sync and check the box to activate the timing synchronization mode.
- Then in the dropdown menu, select the "synchro" with a left click.

To start the acquisition, click on the following icon   
 Having given a name to the experience, click on « Start », the message  
 « pending synchro » appear then the window settings.

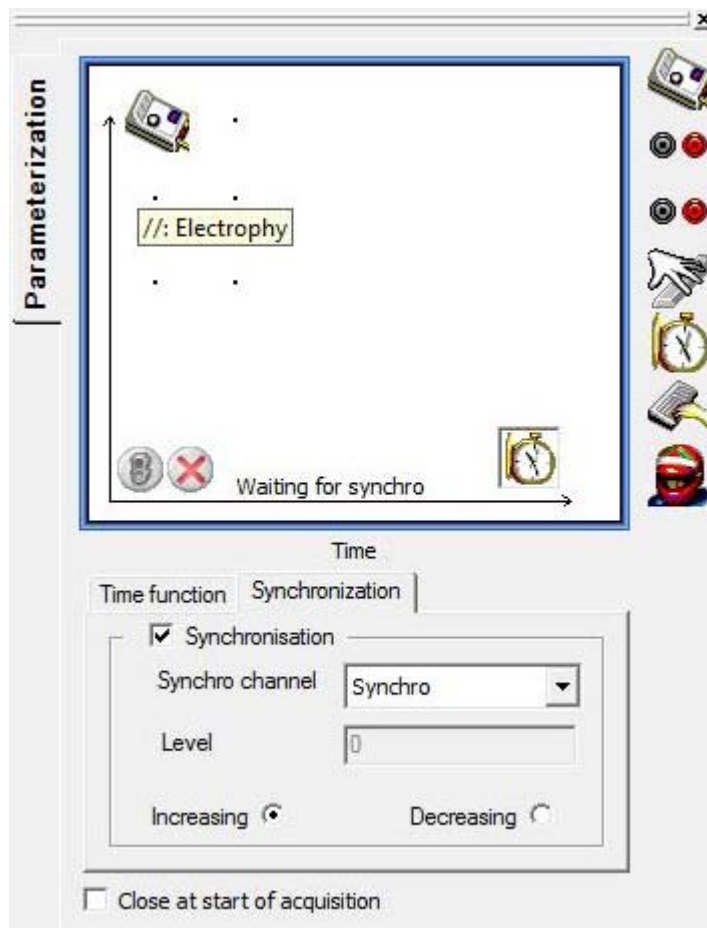The triggering of the acquisition begins when the electronic Martoreflex will be in use.

#### 4. Care and maintenance

For any repairs, adjustments or parts, please contact :

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

## Direct connection for technical support

A team of experts at your  
disposal from Monday  
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

### Free service \*

**+ 33 (0)2 32 29 40 50\*\***

\* Call cost not included

\*\* Only for call from foreign countries