Méthode ESAO[®]

ESAO[®] method

**Réf :
453 084**

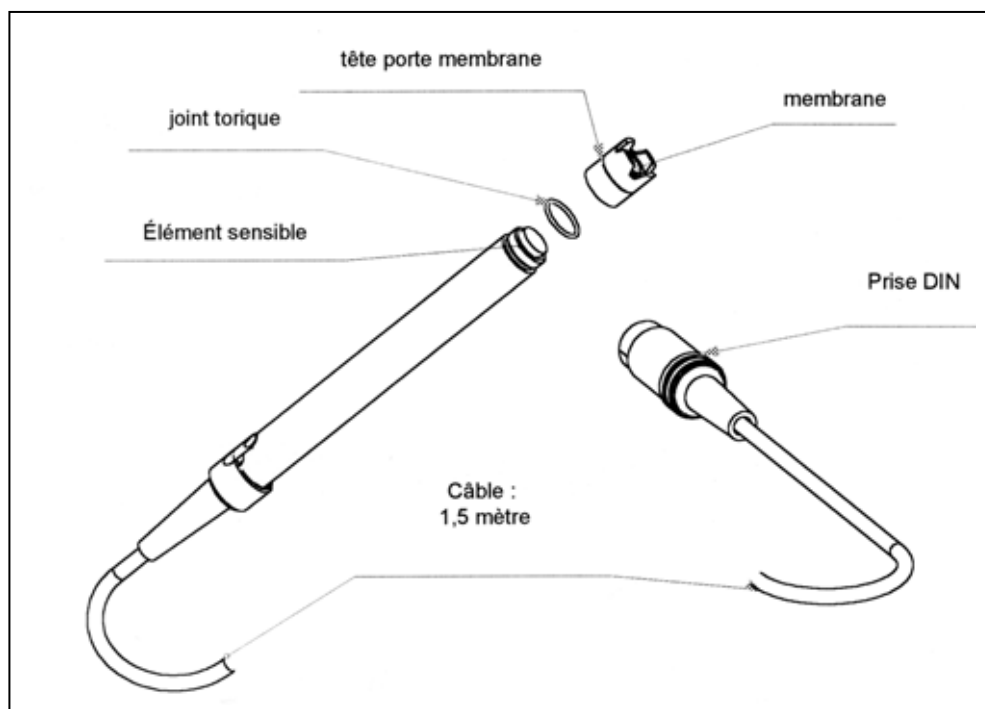
Français – p 1

English – p 5

Version : 1102

**Sonde éthanol
*Ethanol probe***

1 Description

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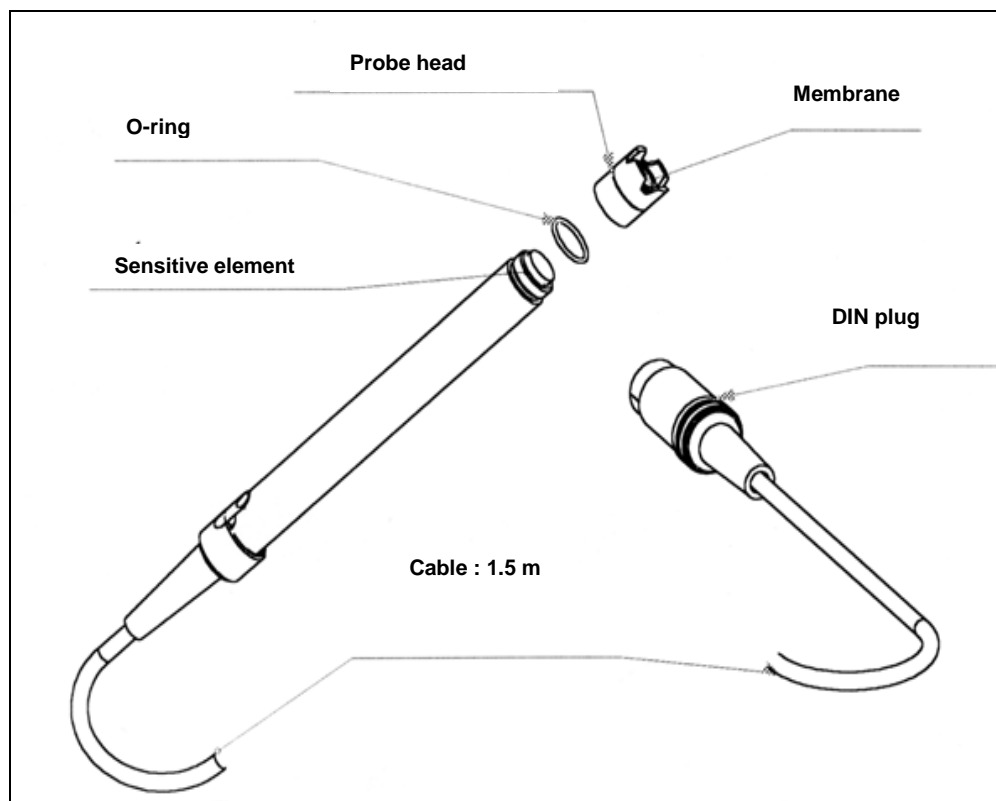
0825 563 563*

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1 Description

The probe is made up of an electronic component that measures the ethanol level. This is a semi-conductor **that is sensitive to water**. It is physically protected by the membrane of the probe head. The probe does not contain any electrolyte.

To increase the service life of the component: do not expose it to very high levels of ethanol. Its maximum measuring capacity is **20 g/l or 2.5 %**.

2 Commissioning

2.1 Connection

Connect the probe to the ESAO® 4 (Part no.: 452 122) or ESAO® 3.1 (Réf : 452 040).

Connect the probe 5 min before use; its temperature must stabilise before it is calibrated.

2.2 Calibration

There is no zero adjustment. Select the rated capacity with which you wish to use the adaptor (in case of any doubt, use the rated capacity 20 g/l).

Adjust the gradient:

1st rated capacity: from 0 to 20g/l (i.e. a maximum level in air and liquids of 2.5 % or 2.5° of alcohol).

2nd rated capacity: from 0 to 1g/l (i.e. 0.125% in air and in liquids).

Quantity	Concentration	Level
20 ml of alcohol /l of water	16 g/l	2 %
1 ml of alcohol /l of water	0.8 g/l	0.1 %

Prepare the calibration solutions by diluting denaturated alcohol (or pure alcohol for better precision) in water.

3 Precautions for use

The sensor does not get spent unless it is exposed to concentrations greater than 25000 ppm of alcohol (20g/l or 2.5%). Do not dip in alcoholic beverages.

Only the end with the sensitive element (equipped with the probe head) may be immersed for measurement.

Do not immerse the probe for more than 1h. The pressure on the membrane should not be too high (avoid placing the probe in a large volume of water).

4 Maintenance

4.1 Cleaning

Clean with clear water, avoid violent jets of water on the membrane.

Check regularly the integrity of the membrane, which ensures a proper discrimination of ethanol and protects the sensor from water.

Changing the membrane: unscrew the probe head and replace it with a new one.

The active component is sensitive to water, it must not be wetted. If you have any doubt about the water tightness of the probe head or the membrane, change the head and the O-ring if necessary (rubber ring).

4.2 Storage

Short-term storage (in between 2 experiments):

Disconnect the probe between every experiment, rinse it if necessary. Store it dry.

Long-term storage (15 days or more):

Store the dry probe after rinsing.

Test:

In case of a doubt regarding the operation of the probe, unscrew the probe head and place the end of the probe on the back of your hand. You will get a slight sensation of heat if the probe is operating correctly.

5 Problems and Solutions

Problems	Causes	Solutions
Not detected by the software or aberrant scale.	Adaptor not configured or improperly configured.	Reconfigure the software, refer to the adaptor instructions.
The value read does not move off zero until the probe is in presence of alcohol.	The DIN plug is not fully plugged in.	Plug in fully.
The curve suddenly disappears OR the value is blocked on full scale, the probe is placed in ambient air (very little alcohol vapour).	The sensitive element is damp, the membrane has a hole in it or the probe head is not screwed on properly.	<ul style="list-style-type: none"> - Remove the probe from the preparation if necessary. - Dismantle the membrane holder head. - If necessary, wipe the sensitive element. - Wait till the value indicated returns close to zero (probe placed in ambient air). - Change the head if necessary. - Recalibrate.
The response time increases with time.	The membrane is partially blocked by deposits.	<ul style="list-style-type: none"> - Replace the probe head with a new one. - You may clean the old head by soaking it in distilled water.

6 Technical specifications

Principle: semi-conductor component-type probe sensitive to dissolved alcohol or alcohol in the form of vapour.

Operating temperature: from 5 °C to 40 °C.

Influence of other gases: carbon monoxide, isobutane, hydrogen and methane to a lesser extent.

Detection level: 0 to 25000 ppm.

Water repellent membrane permeable to ethanol and resistant to diluted organic solvents.

Diameter: 12 mm (adapted to bioreactor 2), length 120 mm.

7 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries