Mécanique

Dynamique

Mechanics

Dynamics

Réf :
453 142

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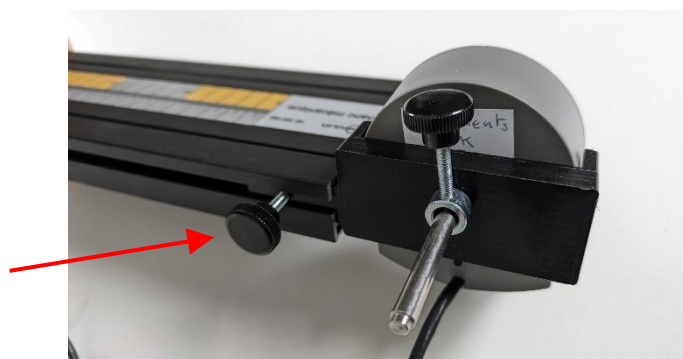
Serrer le bouton de serrage

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Table des matières

1. Content	6
2. Preparing the electronic stepped pulley holder	6
3. Fixing the electronic stepped pulley on the holder	7
4. Fixing the holder on the mechanics bench	7

1. Content

- 1** - 1 electronic stepped pulley holder
- 2** - 1 tightening ring
- 3** 2 tightening knobs

2. Preparing the electronic stepped pulley holder

Screw one of the tightening knobs onto the tightening ring.

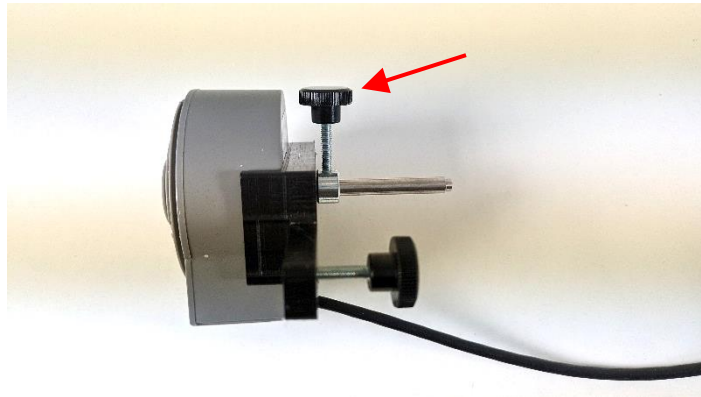
Screw the 2nd tightening knob onto the support.

3. Fixing the electronic stepped pulley on the holder

Insert the pulley onto the holder. Position the tightening ring on the rod of the electronic pulley.


Tighten the tightening knob.

4. Fixing the holder on the mechanics bench

Insert the pulley holder into the mechanics aluminium track up to the stop.

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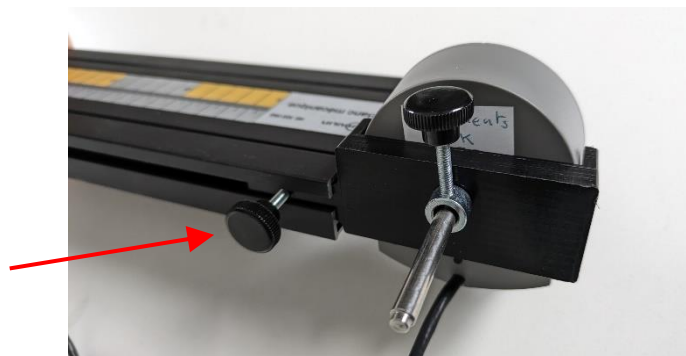Direct connection for technical support

A team of experts
at your disposal
from Monday to Friday
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request
immediatly to provide you
with the right answers regarding
your activity field : Biology, Physics,
Chemistry, Technology.

Free service*

+33 2 32 29 40 50**

* Call cost not included.

** Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux