Foxy
Capteur

Foxy
Sensor

Réf :
482 035

Français – p 1

English – p 3

Version : 2101

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27 019 EVREUX CEDEX FRANCE
0 825 563 563 *

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1 General

This sensor measures the electrical voltage as an instantaneous or effective (TRMS - True Root Mean Square) value.

2 Use

This sensor can be used with any console in the Foxy range.

3 Connection to the Foxy console

- Place the sensor opposite the console connector
- Slide the connector following the console guide
- Push home until it clips in.

3.1 Commissioning

3.2 Use in stand-alone or connected mode

3.2.1 In stand-alone mode

Select the type of measurement chosen: instantaneous or effective (TRMS) using the orange button. A triangle indicates the selected mode.
Select the calibration, a triangle appears in front of the calibration value.
The display value is the value measured continuously.
If the value measures exceeds the selected calibration, dashes appear on the display.

3.2.2 When connecting to the computer

The sensor display switches off. The values measured are visible only on the computer and the buttons become inactive.

4 Additional equipment

No additional equipment is needed.

5 Technical characteristics

- Calibrations: $\pm 50V$, $\pm 15V$, $\pm 2.5V$, $\pm 0,1V$ (display to four significant figures)
- Accuracy: 1% of measurement
- Bandwidth : ≥ 100 kHz at -1%

6 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries