Foxy
Capteur

Foxy
Sensor

Réf :
482 036

Français – p 1

English – p 3

Version : 2101

Capteur thermocouple Foxy

Foxy thermocouple sensor

1 Généralités

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0 825 563 563 *

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1 General

This sensor measures the temperature when combined with a type K thermocouple temperature probe.

2 Use

This sensor can be used with any console in the Foxy range.

3 Connection to the Foxy console

- Place the sensor opposite the console connector
- Slide the connector following the console guide
- Push home until it clips in.

3.1 Commissioning

Adjustment/Calibration: No adjustment is required.

3.2 Use in stand-alone or connected mode

Stand-alone mode offers all the functions of a multimeter.
The instantaneous temperature value is visible on the display.
Changing the calibration is achieved by briefly pressing the calibration button.

A triangle on the display indicates the selected calibration.

When connected to the computer, the sensor display goes out. The measured values are visible only on the computer and the button becomes inactive.

4 Additional equipment

The sensor should be used with a type K probe with flat pins.

5 Technical characteristics

- Calibrations: 0/1000 °C and – 50/+150°C
- Accuracy: 1% ± 2 LSB of value.
- Measuring noise: 0.05% of full scale
- Response time (bare probe and water at 100°C) < 1 s.

6 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries