**Foxy**  
Capteur

**Foxy**  
Sensor

Réf :  
482 037

Français – p 1

English – p 3

Version : 6109

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27 019 EVREUX CEDEX France  
0 825 563 563\*

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## 1 General

This sensor measures luminous lighting up to 50 kLux using the probe supplied.

## 2 Use

This sensor operates with the Foxy console and can be used:

- in stand-alone mode, showing the results on the sensor display
- connected to a computer, displaying the data on a computer

## 3 Connection to the Foxy console

The sensor is connected using the connector located on the unit:

- Align with the chosen measuring channel
- Engage the unit on the console guide
- Push well home.

For a use with the bioreactor, you have to remove the rubber protection of the probe.

## 4 Operation in stand-alone mode

### Commissioning:

- Connect the sensor to the console
- Power-on the console

**Display:**

The measurement is directly on the luxmeter sensor display.

## 5 Operation in connected mode

**Commissioning:**

- Connect the sensor to the console
- Switch on the console
- Connect the console to the computer

**Display:**

The luxmeter sensor display is inactive, the measurements are displayed directly on the computer screen, via the acquisition software.

## 6 Technical characteristics

<b>Measuring range</b>	From 0 to 50,000 Lux
<b>Resolution</b>	12.5 Lux
<b>Accuracy</b>	$\pm (1\% + 2 \text{ LSB})$ of the value
<b>Adjustment/Calibration</b>	None

## 7 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

**JEULIN - TECHNICAL SUPPORT**  
468 rue Jacques Monod  
CS 21900  
27 019 EVREUX CEDEX FRANCE  
+33 (0) 2 32 29 40 50

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## Direct connection for technical support

A team of experts  
at your disposal  
from Monday to Friday  
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request  
immediatly to provide you  
with the right answers regarding  
your activity field : Biology, Physics,  
Chemistry, Technology.

### Free service\*

**+33 2 32 29 40 50\*\***

\* Call cost not included.  
\*\* Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux