Foxy
Capteur

Foxy
Sensor

Réf :
482 051

Français – p 1

English – p 4

Version : 2106

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Foxy blood pressure meter sensor

1 Généralités

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0 825 563 563 *

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1 General

This sensor measures arterial pressure and allows calculating systolic and diastolic pressure. It is used with a pressure measuring armband.

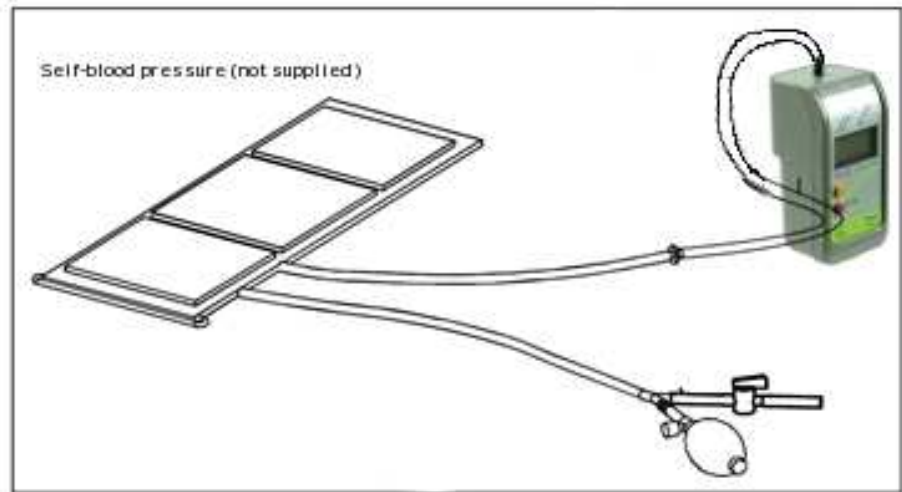
2 Use

This sensor can be used with any console in the Foxy range.

2.1 Connecting the sensor to the console

- Place the sensor opposite the console connector
- Slide the connector following the console guide
- Push home until it clips in.


Connecting the sensor to self-blood pressure

2.2 Commissioning

Adjustment/Calibration: No adjustment is necessary.

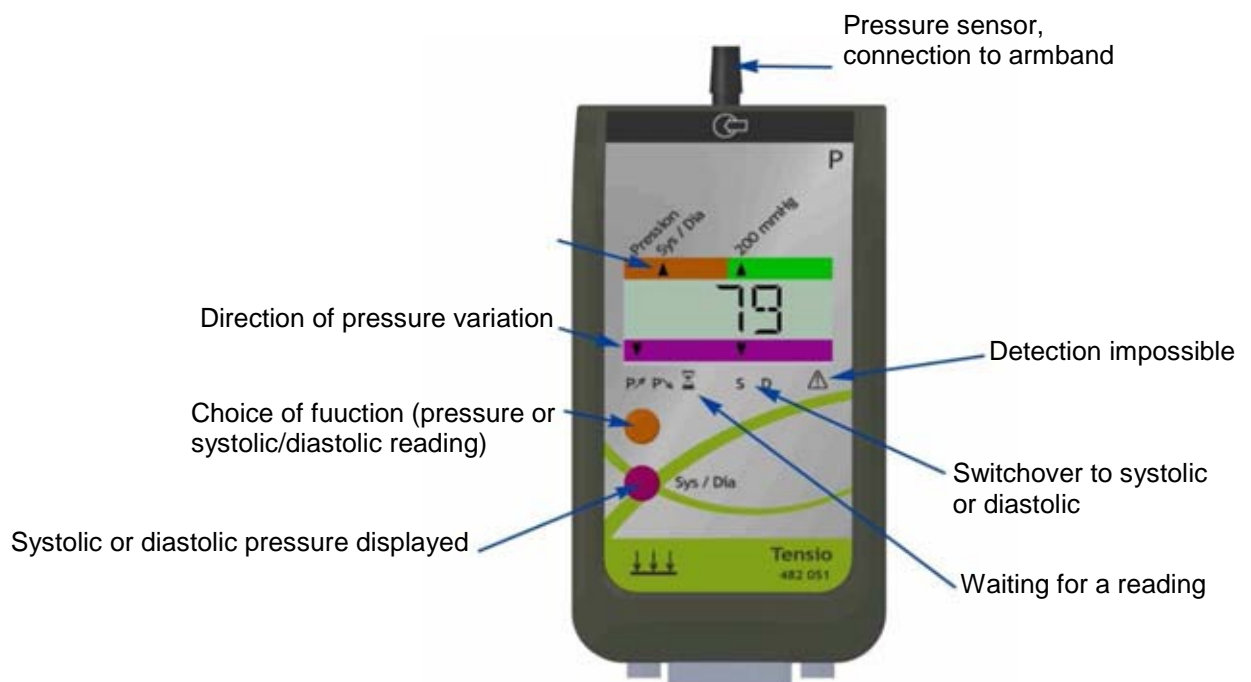2.3 Use in stand-alone or connected mode

2.3.1 In stand-alone mode

The instantaneous pressure is visible on the display.

The sensor offers all the functions of a blood pressure meter.

To determine the systolic and diastolic pressure,

- Select the function (triangle opposite sys/dia) by pressing the orange button.
- Perform a long press on the orange button : the arrow is over the sign **P** 
- Inflate the armband to a value close to 200 mm Hg.
- Open the armband release valve, the arrow indicates the phase of descending pressure then the egg timer indicates the search for the systolic and diastolic values.

After the measurement, press the display switch to display systolic pressure, press again to display the diastolic pressure.

2.3.2 Connecting to a computer

The sensor display goes out. The values measured are visible only on the computer and the buttons become inactive.

The measuring principle is the same as in stand-alone mode, simply follow the instructions on the computer screen.

3 Additional equipment

This sensor is used with a armband inflated using a pump.

4 Technical characteristics

- Calibration: 200 mm Hg, differential sensor
- Accuracy: 1%

5 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
468, Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0) 2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries