Foxy
Capteur

Foxy
Sensor

Réf :
482 053

Français – p 1

English – p 4

Version : 2101

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Foxy ethanol sensor

1 Généralités

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27 019 EVREUX CEDEX FRANCE
0 825 563 563 *

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1 General

The Ethanol sensor combined with the ethanol probe (ref. 453084 not supplied) measures the concentration of ethanol in air (%) or in water (in mg/L). It is particularly suited to experiments involving cellular metabolism, such as alcoholic fermentation of yeasts.

2 Use

The Ethanol sensor can be used in stand alone or connected mode.

2.1 Connection to the console

- Align the unit with the chosen measuring channel.
- Engage the unit on the console guide.
- Push well home.

2.2 Connection of the probe

The sensor is used with the ethanol probe ref. 453084. The probe is connected using the 5 pin DIN socket located on top of the sensor.

2.3 Stand-alone mode

Connect the probe to the sensor 5 minutes before use: the temperature must be stabilised prior to calibration. During this period, an arrow flashes under the egg timer in the display.

Select the required experiment medium (air or water) by pressing the orange button.

Select the experiment calibration by pressing the green button. There are 2 calibrations in air mode (0 to 0.125% and 0 to 2.5%) and 2 calibrations in water mode (0 to 1 g/L or 0 to 20 g/L).

In the case of doubt on the choice of calibration, use the largest one.

2.4 Calibration

Calibration of the ethanol probe is made at a single point.

Depending on the calibration selected, prepare an ethanol solution:

- Calibration 0 to 20 g/L:

Prepare a solution with 20 mL alcohol/L of water and adjust the slope to 16 g/L that is 2% by content.

- Calibration 0 to 1 g/L:

Prepare a solution with 1 mL alcohol/L water and adjust the slope to 0.8 g/L that is 0.1 % content.

Remark: Prepare calibration solutions by diluting denatured alcohol.

To adjust the slope:

- Soak the probe in the chosen calibration solution.
- Press the small violet button for a few seconds. A flashing arrow appears on the violet band of the slope sign.
- Adjust the value displayed using the centre button to display the value closest to the calibration value.
- Validate by briefly pressing the violet button.
- The arrow then appears above OK.

If the arrow appears above the "Caution" sign, check the ethanol probe is operating correctly. For more details consult the instructions for the ethanol probe ref. 453084.

2.5 Connected mode

Start the onboard software. The sensor is recognised automatically.

Select the experiment medium and the calibration using the corresponding tabs in the software.

Calibrate the probe as indicated in paragraph 2.4 "Calibration"*. Use only the central adjustment button displaying the value on the "Measuring" tab.

Remark: The sensor display is inactive when used in connected mode.

* Remark: Of course it is not necessary to perform the calibration sequence each time the software is launched.

3 Technical characteristics

- Calibration at 1 point
- Operating media: 2
 - Air
 - Water
- Calibrations: 2 in Air mode
 - 0 to 0.125 %
 - 0 to 2.5 %
- Calibrations: 2 in Water mode
 - 0 to 1 g/L
 - 0 to 20 g/L
- Response time at 90 %: ≤ 1 second
- Accuracy: 10 % of full scale

4 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries