# Capteur Foxy / Tooxy / Airnext

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## ***Foxy / Tooxy / Airnext sensor***

*Distance measurement*

Réf :  
482 058

Français – p 1

English – p 4

Version : 4109

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***Motion sensor***

## 1 Généralités

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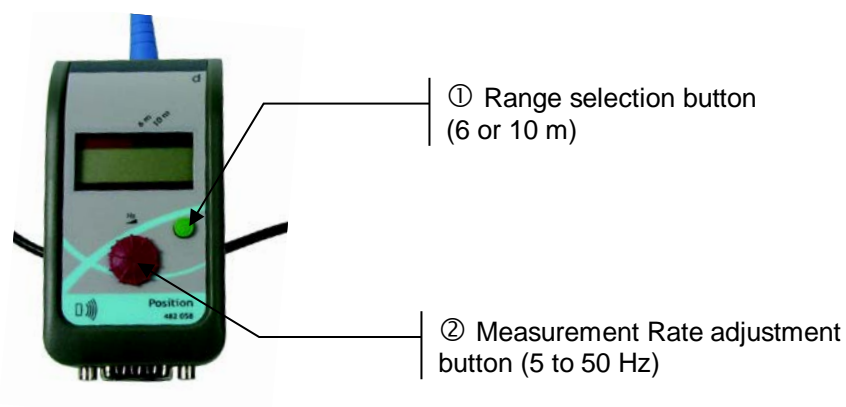
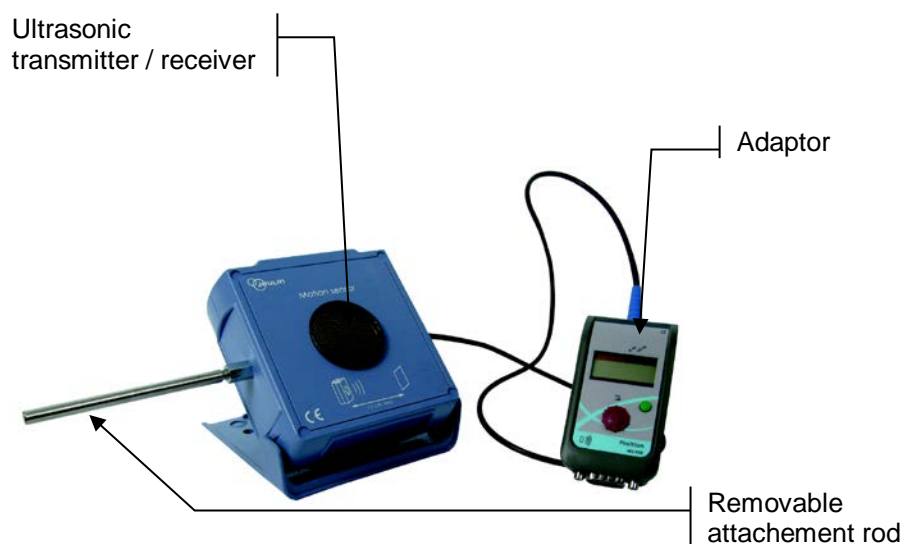
0 825 563 563 \*

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## 1 General description

This sensor, when connected to a Tooxy, Foxy or AirNext, measures the position of fixed or moving objects lying in the range 0.15 m to 10 m along the axis of its transmitter / receiver. Measurements are achieved by the ultrasonic technique.

## 2 Use

### 2.2 Foxy<sup>®</sup>, Tooxy<sup>®</sup> and AirNext<sup>®</sup> systems

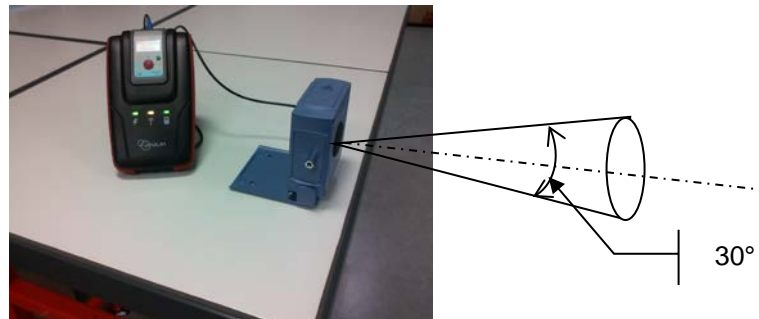
This sensor operates with Foxy<sup>®</sup>, Tooxy<sup>®</sup> and AirNext<sup>®</sup> systems

### 2.3 Initial operations

The objects to be detected must evolve in a cone of about 15° angle around the transmitter / receiver axis. The sensor will detect the closest object in this area.

Make sure to remove all non desired objects from the above defined measurement area.

### 2.4 Special case: object moving on a plane

Aim the motion sensor slightly up to avoid detecting the tabletop or the track on which the object is moving.

Note that this precaution is true for a distance greater than about 80cm. It depends on the inclination of the sensor with respect to the working plane

### 2.5 Measurement

- Connect the adaptor to the console or Pod (in AirNext case)
- Power up the Pod (in AirNext case) and the console
- Using button ① set the suitable range for the planned experiment
- Using button ② set the suitable measurement rate for the planned experiment:

This setting can be operated while measuring in order to monitor its effect and thus adapt it to obtain the best possible result

The fly time technique used to measure the distance implies that:

$$F(\text{Hz}) < \frac{\sim 340}{2 * d}$$

Where F is the measurement rate and d the measured distance

Consider the below table when choosing the suitable rate versus distance

d (m)	≤ 3,4	4	5	6	7	8	9	10
F <sub>max</sub> (Hz)	50	42	34	28	24	21	18	17

### 3 Technical specifications

Range	0.15 m to 10 m
Measurement rate	Adjustable from 5Hz to 50Hz

### 4 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts please contact:

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## Direct connection for technical support

A team of experts  
at your disposal  
from Monday to Friday  
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request  
immediatly to provide you  
with the right answers regarding  
your activity field : Biology, Physics,  
Chemistry, Technology.

### Free service\*

**+33 2 32 29 40 50\*\***

\* Call cost not included.  
\*\* Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux