**Ex. A. O.**

Capteur

***Data logging***

Sensor

Réf :  
**482 083**

Français – p 1

English – p 4

Version : 7112

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***Colorimeter sensor***

## 1. Description

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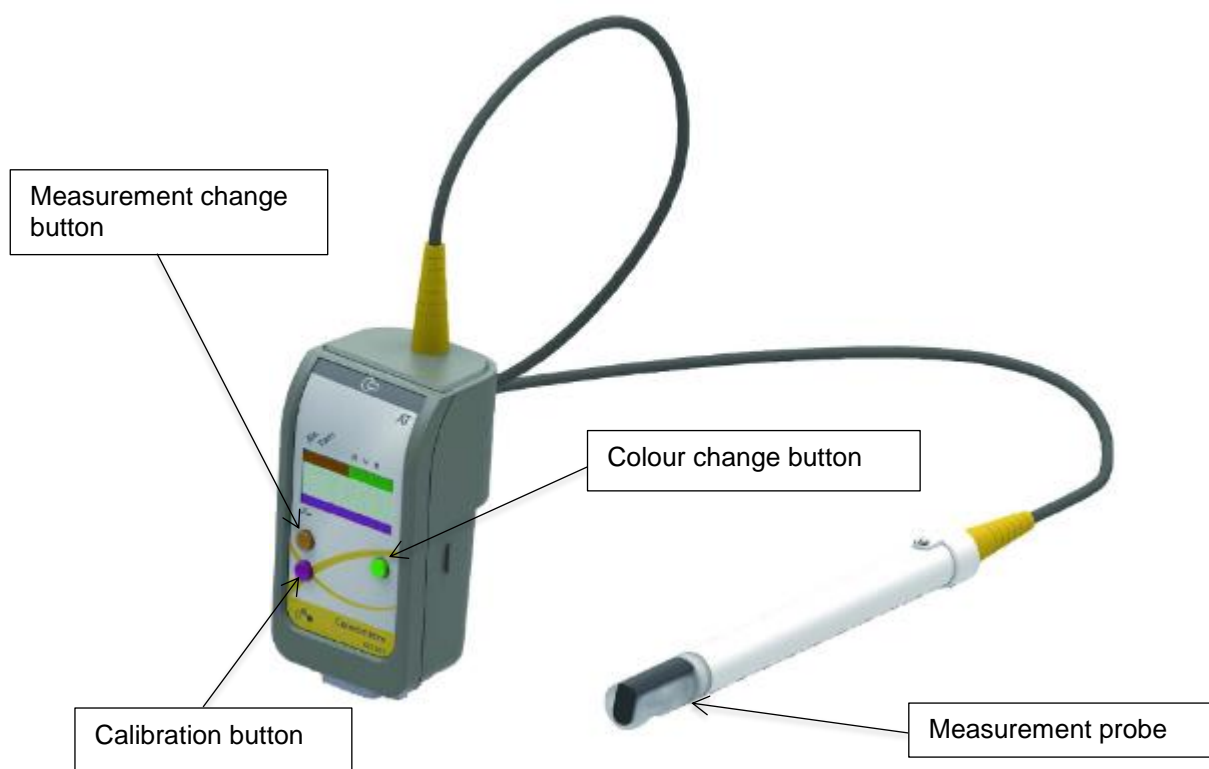
**0 825 563 563\***

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## 1. Description

The colorimeter sensor measures the absorbency and transmittance of a solution. It can be used for data capture or a direct reading via its display.

Three colours are available:

- Red
- Green
- Blue

## 2. Use

Thanks to this probe, the sensor can be used directly in the solution being analysed, without using a special tank. Measurements can therefore be taken in a beaker, the Jeulin bio-reactor or in any other container which can hold a 12mm diameter probe.

For more precise measurements, must be kept dark the enclosure where the measurement is taken.

### 2.1 Datalogging

This sensor is used with Foxy®, AirNeXT® and Tooxy® interfaces

### 2.2 Inserting the sensor

- Place the sensor facing the interface connector
- Slide the sensor following the interface guide
- Push until it clicks

## 2.3 Use in measurement mode

- Select the desired colour using the green button
- Select the type of measurement from absorbency and transmittance using the orange button
- Place the probe in the recipient used for measurement just with the solvent
- Press and hold the purple button, a flashing arrow will appear below the display
- Wait 10 seconds and press the purple button again

The probe is calibrated and ready to use

**Note 1:** It is important to calibrate the probe in the same conditions as the measurements to avoid creating a bias

**Note 2:** For absorbency calibration, the value after calibration is 0. For transmittance calibration, the value after calibration is 100.

## 2.4 Use with an interface

The buttons are inactive when the sensor is connected to an interface. Adjustments are made via the software.

First, click on the  sensor icon and place it on the Y axis.

To adjust measurement and colour, go to the Calibre tab.

For calibration:

- Place the probe in the recipient used for measurement just filled in with the solvent and go to the Adjustment tab
- Click on Start
- For absorbency, enter 0
- For transmittance, enter 100
- To finish, click on adjust

The probe is calibrated and ready to acquire data.

## 3. Technical specifications

### Absorbency

Range	0 to 3 Abs
Accuracy	$\pm 2\%$

### Transmittance

Range	0 to 100%
Accuracy	$\pm 2\%$

### Available colours

Blue	460 nm
Green	520 nm
Red	630 nm

Calibration	At one point
-------------	--------------

## 4. After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.  
For any repairs, adjustments or spare parts please contact:

**JEULIN – TECHNICAL SUPPORT**  
468 rue Jacques Monod  
CS 21900  
27019 EVREUX CEDEX FRANCE

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## Direct connection for technical support

A team of experts  
at your disposal  
from Monday to Friday  
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request  
immediatly to provide you  
with the right answers regarding  
your activity field : Biology, Physics,  
Chemistry, Technology.

### Free service\*

**+33 2 32 29 40 50\*\***

\* Call cost not included.

\*\* Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux