Ex.A.O.

Capteur Foxy/Tooxy/AirNeXT

Data logging

Sensor

Réf :
482 101

Français – p 1

English – p 4

Version : 8110

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0 825 563 563*

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1. Description

This sensor makes temperature measurements within the range from -50 °C to +150 °C.

Measurements are made using a standard 6 mm diameter Pt 1000 probe.

2. Use

This sensor operates with a console in standalone mode (without computer), or in connected mode (with computer) after being connected to the console.

2.1 Connection to the console

The sensor is connected using the connector on the back of the housing:

- bring the sensor in front of the selected measurement channel
- insert the sensor into the console guide system
- follow the console guide system.
- push all the way in

2.2 Measurement type

The sensor measures the temperature on a scale from -50 °C to +150 °C.

2.3 Operation in standalone mode

a. Start up

- connect the sensor
- switch the console on

b. Display

An instantaneous temperature value is displayed directly on the display screen present on the housing.

2.4 In connected mode

a. Start up

- connect the sensor to the console
- switch the console on
- connect the console to the computer

b. Display

Measurements are displayed directly on the computer screen, either during an acquisition or using the “display” tab

3. Technical characteristics

Measurement range	-50/+150 °C
Resolution	0,05 °C
Precision	± (1 % + 2 LSB) of the value
Response time at 90%	≤ 1 minute

4. After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.
For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
468 rue Jacques Monod
CS 21900
27019 EVREUX CEDEX FRANCE

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Direct connection for technical support

A team of experts
at your disposal
from Monday to Friday
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request
immediatly to provide you
with the right answers regarding
your activity field : Biology, Physics,
Chemistry, Technology.

Free service*

+33 2 32 29 40 50**

* Call cost not included.

** Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux