ExAO-NG

Capteur

ExAO-NG

Sensor

Réf :
482 103

Français – p 1

English – p 3

Version : 3106

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0 825 563 563 *

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1 General

This sensor measures the relative humidity of air. The \varnothing 12 mm probe cannot be removed.

Cannot be used in a liquid medium.

2 Use

This sensor operates with the interfaces Tooxy, Foxy and AirNeXT in standalone mode (without computer) or in connected mode (with computer).

2.1 Connecting the sensor to the console

- Bring the sensor in front of the console connector
- Slide the connector along the console guide rail.
- Insert until it clicks into place.

2.2 Start up

Adjustment / calibration: this sensor does not require any adjustment.

2.3 Use in standalone mode and in connected mode

In standalone mode, the sensor displays the value of the relative humidity with a display precision of 1:10.

When the sensor is connected to the computer, the sensor display screen goes off. Measured values are only visible on the computer.

3 Additional equipment

The sensor is supplied with its probe.

4 Technical characteristics

Range	0 to 100 % RH
Resolution	0.25 % RH
Precision	$\pm (1 \% + 2\text{LSB})$ of the value
Measurement time with onboard software	1s

5 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
468, Rue Jacques Monod
CS 21900
27019 EVREUX CEDEX France

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries