ExAO-NG

Capteur

ExAO-NG

Sensor

Réf :
482 104

Français – p 1

English – p 3

Version : 3106

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Pressure sensor

1 Généralités

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0 825 563 563 *

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1 General

This sensor measures the pressure in non-corrosive gases.

2 Use

This sensor operates with the Tooxy console in standalone mode (without computer) or in connected mode (with computer).

2.1 Connecting the sensor to the console

- Bring the sensor in line with the console connector
- Insert the connector along the console guide rail.
- Insert until it clicks into place.

2.2 Start up

Adjustment / calibration: no adjustment is necessary.

2.3 Use in standalone mode and in connected mode

In standalone mode, the absolute pressure value is displayed on the display screen.

When connected to the computer, the sensor display screen goes off. The measured values can only be seen on the computer.

3 Additional equipment

This sensor is used with a 4 mm diameter PVC tube (supplied).

4 Technical characteristics

| | |
|--|---|
| Range | 0 to 2000 hPa |
| Resolution | 0.5 hPa |
| Precision | $\pm (1 \% + 2\text{LSB})$ of the value |
| Measurement time with onboard software | 1 s |

5 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
468, Rue Jacques Monod
CS 21900
27019 EVREUX CEDEX FRANCE

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries