Ex.A.O.

Capteur Foxy / Tooxy / AirNeXT

Data logging

Sensor

Réf :
482 105

Français – p 1

English – p 4

Version : 8106

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Heart rate sensor

1. Généralités

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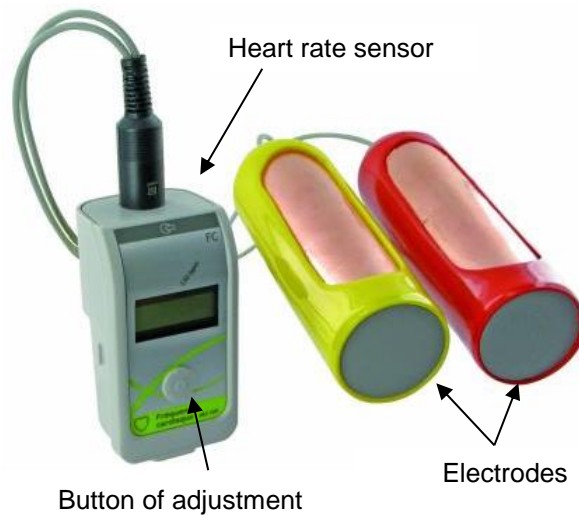
1. General

The heart rate sensor detects and records heart electrical phenomena.

It measures the heart rate in real time.

It is composed of the sensor itself connected to a Tooxy console and two electrodes connected to the sensor through a DIN connection cable.

2. Use

The sensor may be used:

- In standalone mode, displaying data on the sensor display screen.
- In connected mode, connected to a Tooxy console and displaying data on a computer.

2.1 Connection to the console

- Bring the sensor in line with the selected measurement channel.
- Insert the sensor along the console guide rail.
- Insert until it clicks into place.

2.2 Start up

In standalone mode: Display the heart rate on the display screen. Take an electrode in each hand. There is no need to squeeze them too tightly. Preferably take the red electrode in your right hand and the yellow electrode in your left hand.

The heart rate will be displayed after a few seconds. Several arrows are displayed flashing under this value; they should be between 4 and 5. If the displayed number is lower or higher, you should use the adjustment button to adjust the amplification.

In connected mode: Adjust again if necessary (see adjustment in standalone mode), then run the onboard software and the heart rate sensor will be recognised. Start the acquisition.

3. Technical characteristics

Measurement range	0 to 220 bpm (beats per minute)
Measurement precision	± 1 bpm of the maximum value
Number of decimals displayed in standalone mode	1

4. After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.
For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
468 rue Jacques Monod
CS 21900
27019 EVREUX CEDEX FRANCE

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Direct connection for technical support

A team of experts
at your disposal
from Monday to Friday
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request
immediatly to provide you
with the right answers regarding
your activity field : Biology, Physics,
Chemistry, Technology.

Free service*

+33 2 32 29 40 50**

* Call cost not included.

** Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux