**Tooxy**

Capteur

**Tooxy**

Sensor

Ref :  
482 202

Français – p 1

English – p 4

Version : 0112

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***Ammeter sensor***

## 1 Généralités

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0 825 563 563 \*

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## 1 General

This sensor measures instantaneous current intensities of up to  $\pm 1$  A. The measurements are made by a differential input with safety banana jacks (4mm diameter).

## 2 Use

This sensor operates with the Tooxy console in standalone mode (without computer), or in connected mode (with computer) after being connected to the Tooxy console.

### 2.1 Connection to the Tooxy console

The sensor is connected using the connector on the back of the housing:

- Bring the sensor in line with the selected measurement channel
- Insert the sensor along the console guide rail
- Follow the Tooxy console guide system
- Push all the way in

### 2.2 Type of measurement

The sensor measures the current intensity that passes through it.

## 2.3 Choice of the rate

The two available rates are  $\pm 0.1$  A and  $\pm 1$  A.

A button available on the sensor housing is used to modify the sensor range.

To modify the sensor range

## 2.4 Operation in standalone mode

### a. Start up

- Connect the sensor
- Switch the console on

### b. Display

An instantaneous current value is displayed directly on the display screen present on the housing.

## 2.5 In connected mode

### a. Start up

- connect the sensor to the console
- Switch the console on
- Connect the console to the computer

### b. Display

Measurements are displayed directly on the computer screen, either during an acquisition or using the "display" tab.

### 3 Technical characteristics

Measurement range:	- range $\pm 0.1$ A - range $\pm 1$ A.	from $-0.1$ A to $0.1$ A from $-1$ A to $1$ A
Resolution	- range $\pm 0.1$ A - range $\pm 1$ A.	$50\mu\text{A}$ $500\mu\text{A}$
Precision	- range $\pm 0.1$ A - range $\pm 1$ A.	$\pm (1\% + 2 \text{ LSB})$ of the value $\pm (1\% + 2 \text{ LSB})$ of the value
Passband	$\geq 10$ kHz	
Protection	Up to $10$ A per resettable fuse	

### 4 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

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## Direct connection for technical support

A team of experts  
at your disposal  
from Monday to Friday  
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request  
immediatly to provide you  
with the right answers regarding  
your activity field : Biology, Physics,  
Chemistry, Technology.

### Free service\*

**+33 2 32 29 40 50\*\***

\* Call cost not included.

\*\* Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux