Anatomie

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Anatomy

Model - dissection

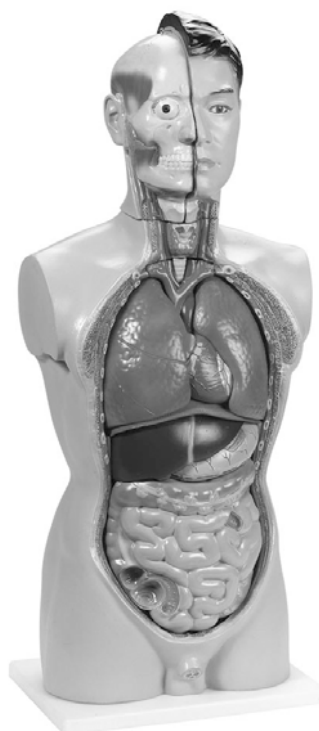
**Ref :
512 062**

Français – p 1

English – p 7

Version : 6010

Tronc humain asexué
Asexual human trunk

NOMENCLATURE

LA TÊTE

Le crâne

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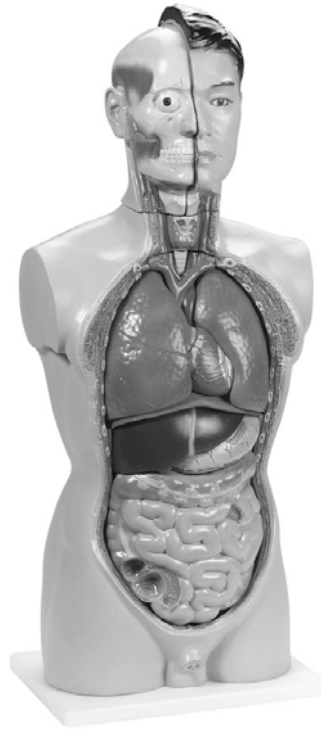
The brain

1. Dura mater
2. Braincase
3. Frontal lobe
4. Parietal lobe

5. Occipital lobe
6. Temporal lobe
7. Cerebellum
8. Corpus callosum
 - a. face
 - b. knee
9. Transparent wall
10. Thalamus. The third ventricle of the brain is found between the two thalamus
11. Interventricular hole (Hole of Monroe, allowing the third ventricle to communicate with the two lateral ventricles).
12. Aqueduct of Sylvius, joining the third and the fourth ventricle
13. Epiphysis or pineal gland
14. Dome of the arch
15. Medullar body
16. Pons Varoli
17. Medulla oblongata
18. Oliva medulla oblongata
19. Mamillary bodies
20. Optic chiasma
21. Hypophysis
22. Optical nerve
23. Olfactory nerve
24. Trigeminal nerve
25. External ocular motor nerve
26. Facial nerve
27. Auditory nerve

THE TRUNK

Region of neck

1. Hyoid bone
2. Thyroid median cartilage
3. Thyroid lateral cartilage
4. Thyroid cartilage
5. Crico-thyroid muscle
6. Thyroid gland
7. Tracheal cartilage
8. Ligament between the tracheal cartilages

The heart

1. Right ventricle
2. Aorta
3. Cross of the aorta
4. Thoracic aorta
5. Left auricle
6. Left ventricle
7. Pulmonary artery
8. Right auricle
9. Superior vena cava

10. Inferior vena cava
11. Pulmonary vein

The trachea

12. Bifurcation of the trachea
13. Right lung
14. Left lung
15. Bronchial branches

"A" Right lung

16. Superior lobe
17. Median lobe
18. Inferior lobe

"B" Left lung

19. Superior lobe
20. Inferior lobe
21. Oesophagus
22. Diaphragm

Abdominal cavity

1. Spleen
2. Right kidney
3. Renal cortex
4. Medulla
5. Renal pyramids
6. Right renal artery
7. Left adrenal
8. Left kidney
9. Left ureter
10. Pelvis.
11. Abdominal aorta

Liver

12. Liver
13. Right lobe
14. Left lobe
15. Quadrate lobe
16. Caudate lobe
17. Hepatic artery
18. Portal vein
19. Hepatic duct (excretion of liver)
20. Cystic duct (excretion of gallbladder)
21. Common bile duct (joining the above two)
22. Gallbladder

Stomach

- 23. Stomach
- 24. Cardia
- 25. Fundus of the stomach
- 26. Pylorus
- 27. Small curve
- 28. Large curve
- 29. Pancreas

Small intestine

- 30. Duodenum
- 31. Jejunum
- 32. Ileum

Large intestine

- 33. Caecum
- 34. Appendix
- 35. Colon :
 - a. Ascending colon
 - b. Transversal colon
 - c. Descending colon
- 36. Sigmoid loop
- 37. Rectum

Urogenital system

- 1. Pubis
- 2. Right ureter
- 3. Bladder
- 4. Prostate
- 5. Urethra: prostatic section
- 6. Seminal vesicle
- 7. Vas deferens

Mounting of the different organs

During the mounting, the organs are placed from the bottom of the abdominal cavity towards the head.

The notions of right and left refer to the trunk and not to the user.

- 1. Place the bladder (**3**), the number should be visible, the prostate (**4**) behind and oriented downwards.
- 2. Place the right half-kidney (**2**), press the pin in position.
- 3. Place the oesophagus (**21**)/thoracic aorta (**4**), press the pin in position.
- 4. Place the trachea on the oesophagus (**7**), press the pin in position.

5. Place the left half-lungs not numbered in the dorsal part, place the right half-lung not numbered in dorsal part.
6. Place the spleen (1)/pancreas(29)/duodenum(30): the spleen is inserted under the dorsal left half-lung, then insert the pin in position.
7. Place the intestinal organs (30-31-32).
8. To be mounted separately: place the liver (12) in the diaphragm (22). Insert the stomach (23) under these organs; the cardia of the stomach (24) fixed in at the level of the small opening of the diaphragm.
9. Fit this mounting diaphragm/liver/stomach on the transversal colon (35 b). The stomach is placed in front of the spleen (1) and the diaphragm under the lungs. Lift them slightly; pull the intestine to the front if necessary for holding all the organs.
10. Place the heart: the pins in top position enter the slots provided for this purpose on each side of the thyroid gland (6). The arteries (red) and the veins (blue) should be continuous.
11. Place the right and left half-lungs. Fix them under the ribs.
12. Insert the brain into the skull.
13. Place the left side of the head in the pins provided on the trunk, then assemble the right half, hold the model in place with the hook.

After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediately to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service * :
+ 33 (0)2 32 29 40 50

** Call cost not included*