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Faune terrestre

Environment

Fauna

Ref :
525 007

Français – p 1

English – p 3

Version : 6010

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
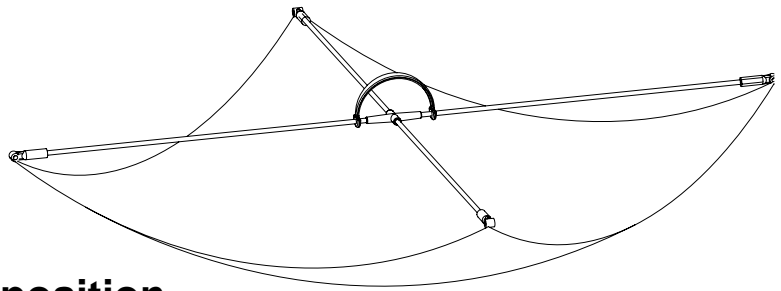
1 Composition

- 1 rod of 100 cm (1)
- 2 rods of 50 cm (2) and (2')
- 1 cross with hook (3)
- 1 handle (4)
- 1 washable cloth net (75 x 75 cm) with 4 junctions (5)

2 Procedure

2.1 Set-up:

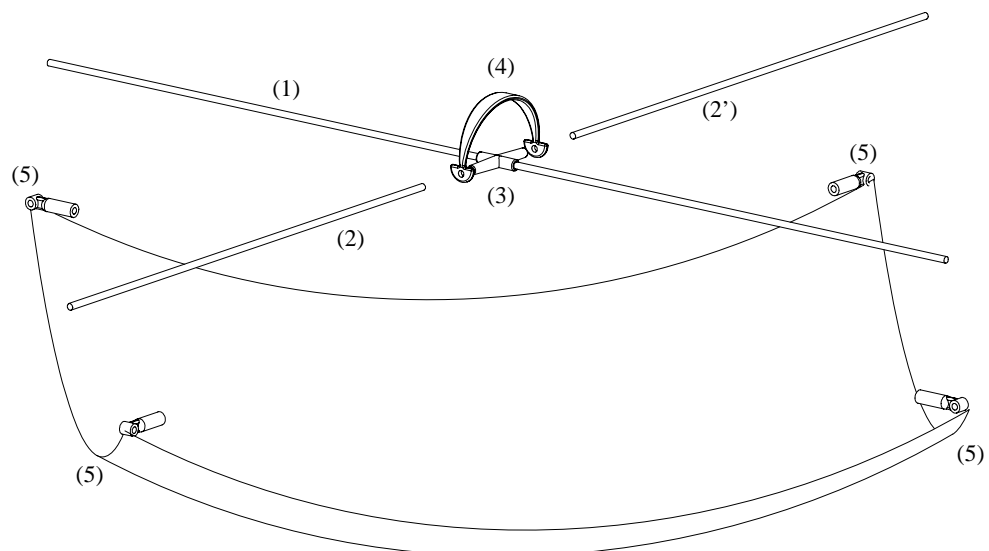- Fit and slide the cross with hook (3) on the large rod (1).
- Position the handle (4) on each side of the cross and slide in the two small rods
- (2) and (2').
- Fit each of the four junctions (located on the corners of the cloth) (5) on the frame mounted previously.

2.2 Collection of fauna

- Place the net under the branches to be studied.
- Hit the branch with a rod or a stick.

- Let the insects drop into the umbrella, collect them carefully without damaging them (using an insect collector, for example) and transfer them into a ventilated flask for observing them.
- Label the flasks (place of collection, tree no., date, etc).
- For observing the collection extemporaneously, fix the animals for preserving them (ethanol 70°).

3 Suggestions for usage

- Discovery of the fauna in trees.
- Discovery of the fauna in bushes.
- Comparison between trees or bushes of different species.
- Estimation of the biomass of invertebrates in the fauna.
- Comparison of the fauna before and after the haying season, during different seasons, between two years, before or after chemical treatment, etc.
- Reflection on the process employed and the strategies used
- Reflection on the preservation of individual collected.

4 Storage

Dismantle the net and the frame before storing it in its original packing.

Shake the net vigorously for cleaning it before folding. If it is wet, allow it to dry in order to prevent fungus growth.

It is possible to wash the net if necessary (hand wash, temperature less than 30°C).

5 Additional equipment required

- Insect collector Ref. **525 001**
- Ventilated flask for collection of fauna.
- Sample labels.
- Ethanol 70°.
- Stick or rod.

6 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediately to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service * :
+ 33 (0)2 32 29 40 50

** Call cost not included*