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Faune terrestre

Environment

Fauna

Ref :
525 023

Français – p 1

English – p 5

Version : 7004

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Berlese funnel student

1 Fonction

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NOTES

1 Function

The material is intended for extraction of living beings from the soil: the students thus discover that this environment often shelters a wide variety of fauna.

2 Composition

The product is made up of four parts:

- A black shockproof polystyrene container, for receiving the soil sample (manure or other material), Ø external 94 x height 55 mm
- A sieve Ø 86 mm with plastified grids retaining the soil but allowing the organisms to pass through
- A support/funnel unit in shockproof black polystyrene whose surface and the slope are designed for guiding the organisms downwards; Ø max. 100 x height 80 mm
- A collection pot in transparent polyethylene Ø 70 (base) and 88 (top) x h 70 mm.

It is delivered with two accessories:

- A brush held in place by a clip for removing the organism from the collector without damaging them and placing them in position for observation (mounting glass, Petri dish, cavity slide...)
- A small magnifying glass whose handle fits into a ring for a primary observation of the animal forms collected.

3 Installation

If the equipment was stored dismantled:

- Fit the funnel on the collection pot (or if you prefer, place it on other recipient) where some water or preserving liquid (water + alcohol) should be placed.
- Place the sieve on the funnel.
- Place the container above the sieve so that it is correctly fixed.

❖ **Equipment necessary:**

- An articulated lamp with reflector,
- Alcohol for preparing the preserving liquid, for example ethanol denatured.

❖ **Accessories recommended:**

- A binocular lens,
- Cavity slides,
- Or for sorting with the forceps, 1 slide with two cavities,
- Or for allowing the animals to move without losing them opaque slide with sliding cover slip.

4 Usage

- Fill the container with the sample. Record its condition: compact/gritty/sandy, dry/wet, rich/poor in plant debris, etc. Adjust the reflector of the lamp very close to the manure (about 1 cm) and light it during about 15 minutes: the organisms are driven downwards by the strong light and heat, pass through the sieve and fall on the smooth surface of the funnel that guides them into the collector.
- Turn off the lamp and remove it. Remove from their position the magnifying glass and the brush carefully (without shaking), lift the apparatus vertically and separate the collector from the funnel + sieve + container + sample assembly by tilting in the direction of a window of the funnel.
- Collect the organisms with the brush and place them on the slide or watch glass; observe them with the magnifying glass, then with the binocular lens, sort them according to their morphology: this may be a first approach to their classification. Correlate their number with the condition of the sample collected initially.

5 Maintenance

Clean all the components for running water without rubbing and allow to dry. Avoid the use of any abrasive material that may damage the surface conditions, affecting the efficiency of the funnel or the transparency of the collector.

6 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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Service gratuit * :
+ 33 (0)2 32 29 40 50

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Rubrique FAQ

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

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