Loupes

Magnifiers

**Ref :
571 117**

Français – p 1

English – p 5

Version : 7004

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NOTES

1 Purpose

The binocular magnifier is meant for the stereoscopic observation of any biological or geological material, which cannot be observed under the microscope due to its thickness and/or its opacity.

2 Description

- Fixing screw with lock mechanism: to prevent the body of the magnifier from falling in case of uncontrolled releasing.
- Reversible disc (black and white)
- Base forming the stage plate: 170 x 115 mm.
- Disc centred on the optical axis:
 - . diameter 60 mm.
 - . in unbreakable altuglass.
 - . removable
 - . reversible (black side, white side).
- Magnifier body support: Metal column: $h = 150$ mm, diameter 14 mm. Wide base (diameter 20 mm) forming a stop.
- Body of the magnifier consisting of the optical block and an arm.
- Fixing of body on the column: by fixing screw (knurled knob).
- Up and down travel stops.
- Body capable of rotating in all directions.
- Straight binocular head.
- Eyepiece separation adjustment.
- Independent adjustment of one of the eyepieces

- Rapid adjustment: positioning of the magnifier body by the lock screw.
- Locking device at the level of the fixing screw of the body on the column (ensures that the body of the magnifier does not fall in case of uncontrolled releasing).
- Focussing through dual knob rack and an adjustable slip clutch (prevents damage of mechanism at end of travel).
- A pair of wide field eyepieces with large aperture (diameter of front lens 15 mm) 10 X.
- A pair of interchangeable fixed achromatic objectives 2 X.
- A pair of collapsible eye cups in rubber fitted on the eyepieces.
- Anti-reflection coating on eyepieces and objectives.
- Total magnification: x 20.
- Optical field: diameter 10 mm.
- Maximum height of the sample for observation: 90 mm.
- Maximum amplitude of fine adjustment: 40 mm.
- Total amplitude (all adjustments taken into account) : 135 mm.
- Delivered in a polystyrene box with dust cover.
- Size, weight of magnifier:
 - max.: 330 X 170 x 115 mm – 1.48 kg.
 - minimum: 210 x 170 x 115 mm.
- Size, weight (magnifier + box):
 - 310 x 220 x 140 mm – 1.650 Kg.

3 Operating instructions

3.1 Preparation

- Take out the magnifier from its packing.
- Place the eyecups on each eyepiece: these facilitate observation by shielding stray light.
- Choose the white or black side of the central disk of the stage plate depending on the object to be observed.
- Place the object to be observed on the stage plate.

3.2 Focussing

❖ Rapid pre-adjustment:

- Loosen the fixing screw of the magnifier body on the column (a locking mechanism prevents its accidental fall.)
- Slide the body so that it is placed at around 7 or 8 cm above the object.
- Tighten the fixing screw.

❖ **Focussing:**

- Adjust the eyepiece separation.
- Focus by using the dual knob of the rack.
- Adjust vision by turning the knurled ring of the adjustable eyepiece.

3.3 Special cases : dissection

For objects that are particularly bulky, which cannot be placed on the stage plate, just rotate the body of the magnifier on its support after having unscrewed the fixing screw (possible rotation 360 °) around the axis.

4 Maintenance

The binocular magnifier is delivered in a polystyrene box. A cover is provided to protect the magnifier from dust when not in use.

The tightness of the focussing mechanism can be adjusted by a special key delivered along with each instrument

All mechanical parts of the magnifier are in metal. These must be cleaned regularly with a soft cloth soaked in acid free Vaseline. If required clean with spirit before lubricating.

During cleaning and lubricating, protect the optical parts of the magnifier.

In order to ensure dust free interior, never remove the eyepieces.

Never touch the lenses. Clean them as and when required with an optical cleaning paper Reference 803 092

Never dismantle the eyepieces and objectives for cleaning. If necessity arises, return them to us.

Transport: while transporting, carefully pack the magnifier in its original box to protect it from possible shocks.

5 After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.

For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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Service gratuit * :
+ 33 (0)2 32 29 40 50

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Rubrique FAQ

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

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