Observation

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Observation

Magnifiers

Réf :
571 120

Français – p 1

English – p 5

Version : 7001

**Loupe binoculaire série
"T" X20 X40**

***Binocular magnifier serie
« T » X20 X40***

1 Objectifs

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NOTES

1 Purpose

The binocular magnifier is meant for the stereoscopic observation of any biological or geological material, which cannot be observed under the microscope due to its thickness and/or its opacity.

2 General features

- Ring mounted optical head, 360° rotatable
- Adjustable column height
- Base fitted with two clamps fixed by screws to prevent its loss.
- Focussing by smooth and reliable bilateral control.
- Eyepiece separation adjustment and dioptré adjustment on the left eyepiece holder.
- Wide field eyepieces inclined for your comfort, fixed by screws, and flexible rubber eyecups.
- Good field depth; high optical resolution.
- Dust cover.

12 V 10 W low voltage top and bottom lights – Top light used for opaque objects (reversible opaque stage plate) – Bottom light used for translucent objects (reversible opaque stage plate).

3 Optical features

Pair of eyepiece	Pair of parfocal objectives on turret	Resultant magnification	Visual field	Working distance
x 10	x 2 et x 4	x 20 et x 40	10 mm and 5 mm	48 mm

4 Utilisation and adjustment

4.1 Utilisation and adjustment

This can be adjusted to suit the thickness of the object and the selected magnification (option, see accessories at the end of the manual), with the help of the rear knurled knob, to be unscrewed while holding the body of the magnifier.

4.2 Objectives and eyepieces

The wide field eyepieces x 10 are maintained in position by screws, in order to prevent loss or accidental fall. Their range of separation is adjustable from 51 to 75 mm.

The left eyepiece holder is fitted with a dioptré setting ring(± 5 mm) : set the focus for the right eye (see below), then without disturbing the focus, observe with the left eye and turn the ring until a clear image is obtained.

4.3 Focussing

The instrument can be adjusted either with the right or the left hand by the dual control for the rack and pinion system which is located on both sides of the column. At the end of travel, the ring of each control knob slips without driving the pinion, thus preventing any deterioration of the rack.

4.4 Change in magnification

By a simple rotation of the turret, we can move from the objectives x 2 to the objectives x 4 without altering the focus.

A screwdriver is needed to change the eyepiece. Note: Do not lose the fixing screws. Various pairs of eyepieces x 5, x 15, x 20 form part of the list of accessories.

5 Maintenance

5.1 Protection from dust

After use, immediately place the dust cover on the instrument. If dust has fallen on the lenses, use a syringe or a rubber bulb blower or a sable brush: the lenses are fragile and dust can cause scratches.

5.2 Finger marks

Do not touch the lenses. In case of fingerprints remove them with the help of a soft cloth soaked in xylene.

5.3 Humidity

Humidity, especially when combined with high temperatures is particularly harmful to optical elements. It is preferable to store the instruments in a dry place and in case of very exposed climates, use the polystyrene boxes with desiccative sachets.

5.4 Dismantling

In case the binocular magnifier needs to be dismantled (for maintenance), the instrument must be entrusted to the supplier or to a qualified workshop.

5.5 Bulb replacement

Unplug the 3 pin lead :

- Bottom lighting: open the clamps and lift the stage plate to easily access the bulb.
- Top lighting: unscrew the knurled knob located on the right of the illuminating block. The bulb can be seen after removing the lamp body.

6 Accessories

Spare bulb Reference 574 152.
Fuse-shaped bulb 12 V and 10 W.

7 110 Volts power supply

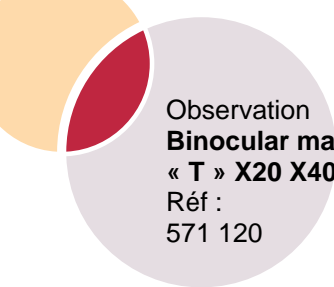
- 1- Unplug the instrument from the power supply
- 2 – Unscrew the backplate (4 Phillips star head screws).
- 3 – Identify the 3 transformer primary wires: black = neuter, yellow = 117 V, red = 230 V. A brown wire is connected at the end of the red wire. Cut the end of the brown wire and bare the wire for 0.5 cm . Also cut the end of the yellow wire and bare the wire for 0.5 cm. Connect the yellow and brown wires with a connecting block.
- 4 – Replace the blackplate
- 5 – Stick a label 110 V at the back of the instrument above the main lead.

8 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service * :
+ 33 (0)2 32 29 40 50

** Call cost not included*