Microscopie

Microscopy

**Ref :
571 234**

Français – p 1

English – p 5

Version : 4101

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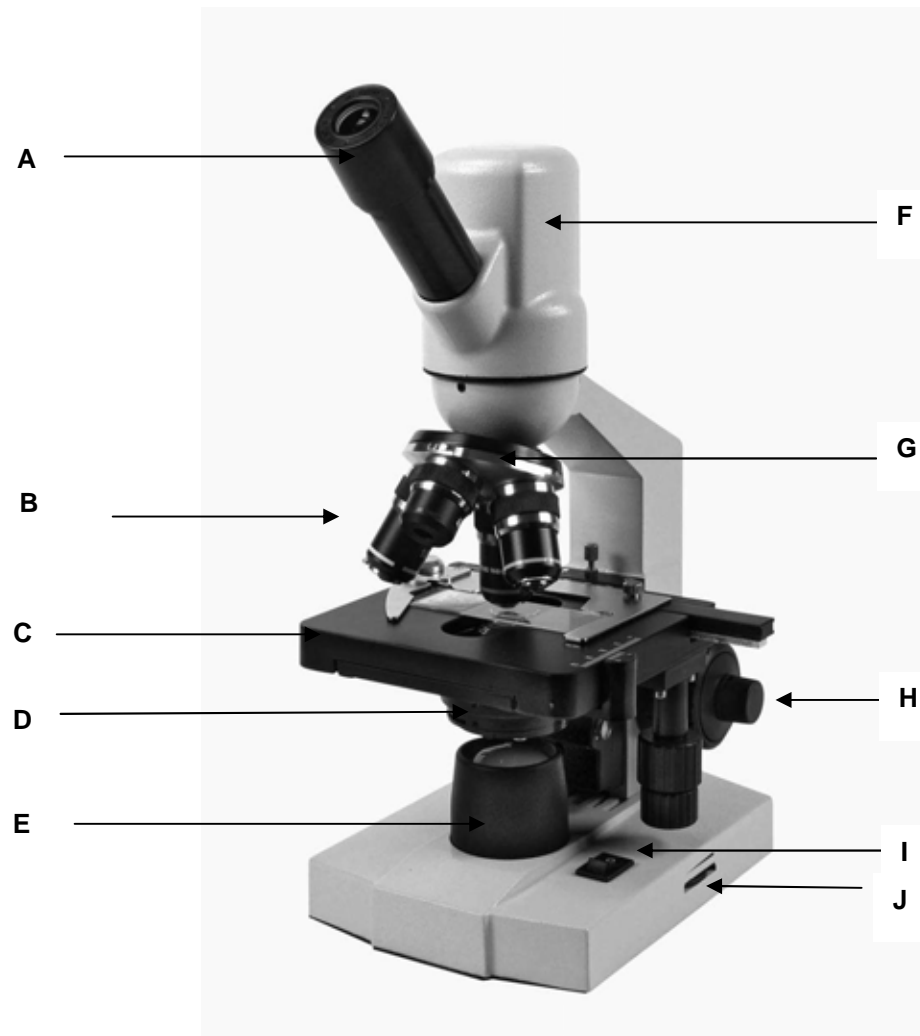
A Wide field ocular

B Objectives

C Rectangular stage with over-stage

D Condenser, iris diaphragm

E LED illumination

F Monocular head with camera

G Triple nosepiece

H Coarse and fine coaxial adjustment

I Dimmer switch

J On/Off

2 Getting started with the microscope

Take the microscope out of its packaging, remove the various protections.
Keeps the package for extended storage or possible future shipping.

Place the eye piece in the eyepiece holder tube.

3 Getting started with the camera

1. Insert the CD delivered with a microscope in your drive.
2. Link the microscope to your computer using the USB cable supplied.
3. The camera is detected. The classic Windows installation process starts. Wait until it is completed.

The camera can now be used with your image capture and processing software, e.g. Picturio.
Scopelimage software is supplied with the camera.

4 Installing the software supplied

When you insert the CD, a window is displayed :

Click on software and the installation process will start up.
Follow up the process until it is completed.
Reboot your computer.

Your video microscope is now ready for use.

5 Use and adjustment

- To move, hold the microscope exclusively by the inclined part of the stand.
- Switch on and adjust the brightness using the switch on the right-hand side of the base to obtain the desired lighting.
- Insert the preparation.
- The 360° adjustable observation head and the tube inclined at 45° guarantee excellent ergonomics and allow for extended use.
- The para-focal lenses facilitate focusing when changing the enlargement. The 40X and 100X lens have a retractable mount which avoids damaging the preparation in the event of accidental contact.
- Focus by vertical movement of the plate using a double coaxial macro and micro knurled knob for fine adjustment.
- A light condenser and an adjustable iris diaphragm are located under the plate.
- Adjust the lighting by closing the diaphragm to increase the contrast and depth of field but without exaggerating so as to affect the resolution.

6 Technical features

- Stable and robust metal stand. Ergonomic low-position controls so the user's forearms are supported by the table.
- Monocular head with integrated camera, rotating through 360° with eye piece holder tube inclined at 45°.
- Large field eye piece, 10 X.
- 4 position turret for the following lenses: achromatic lenses 4 X, 10 X, 40 X, 100X (can be immersed).
- Rapid coaxial focusing using a rack.
- Rectangular plate with cover.
- Digital 0.65 opening condenser, with iris diaphragm.
- LED lighting incorporated in the base with on/off switch and intensity adjustment.

7 Camera features

Sensor type	1/3"CMOS
Resolution	1M Pixel
Speed image acquisition	12fps
Capture format	JPEG, bmp, TIFF, PNG, vidéos avi
Login	USB 2.0

8 Maintenance

- Protection from dust: After use, place the cover over the apparatus. If there are dust deposits on the lenses, blow them away with the help of a syringe or dust remover (item no. 574 198) or use a martin hair brush; the lenses are fragile and can be scratched by dust.
- Fingerprints: Do not touch the lenses. If, however, they do have fingerprints on them, remove the fingerprints with a soft cloth or lens paper (item no. 803 092).
- Humidity: Humidity, especially in combination with temperature, is particularly harmful to the assembled optics. It is a good idea to store the microscopes in a dry place, and under especially exposed climate conditions, the microscopes should be stored in their polystyrene boxes with a dessication bag.
- Disassembly: A microscope is a finely-tooled, precision instrument. If it is necessary to take it apart (for maintenance), the instrument should be returned to JEULIN or a qualified repair shop. The base plate that seals the microscope may only be opened by an electrician because the body must remain protected according to electrical safety standards. Do not take the microscope apart unless it is unplugged.

9 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
468, Rue Jacques Monod
CS 21900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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Direct connection for technical support

A team of experts
at your disposal
from Monday to Friday
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request
immediately to provide you
with the right answers regarding
your activity field : Biology, Physics,
Chemistry, Technology.

Free service*

+33 2 32 29 40 50**

* Call cost not included.
** Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux