Observation

Stéréomicroscopes

Observation

Stereomicroscopes

Réf :
188 495
571 267
571 286

Français – p 1

English – p 3

Version : 3107

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LED stereomicroscopes

1. Description

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0 825 563 563 *

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1. Description

Light because of their plastic conception, these stereomicroscope are provided to observe subjects in relief with low magnification.

▲ x25 eyepieces

2. Technical specifications

	Ref 188 495	Ref 571 267	Ref 571 286
Head	Binocular, straight		Binocular, 45° inclined
Eyepieces	WF10X	WF10X - WF25X	
Interpupillar distance	55 - 75 mm		
Dioptric adjustment	No		
Objective	x 2		
Magnification	x 20	x 20 and x 50	
Field of view	15 mm / x 20	15 mm / x 20 - 6 mm / x 50	
Working distance	70 mm		
Macrometric adjustment	40 mm		
Max thickness of subjects	25 mm		
Stage	White disc 50mm - Metal clips with spring		
Illumination	LED incident		
Power supply	2 batteries 1,5V LR06 AA (not included) Switch ON/OFF		
Delivered with	Dust cover, 12 rocks in box		Dust cover
Dimensions	Base: 100 x 135 mm / Height: 225 à 262 mm max		
Weight	Around 600g		

3. Set-up

Remove the stereomicroscope from its packaging.
Insert 2 batteries 1,5 V (not included), considering the polarity.
Turn on the stereomicroscope: it is ready to work.
Place your subject on the white disc and fix it with the clips if necessary.
With the macrometric knob, adjust to get a clear image.

4. Maintenance

- Protection from dust: After use, place the cover over the apparatus. If there are dust deposits on the lenses, blow them away with a syringe or dust remover (item no. 574 198) or use a martin hair brush; the lenses are fragile and can be scratched by dust.
- Fingerprints: Do not touch the lenses. If, however, they do have fingerprints on them, remove the fingerprints with a soft cloth or lens paper (item no. 803 092).
 - Humidity: Humidity, especially in combination with temperature, is particularly harmful for assembled optics. It is suitable to store the stereomicroscopes in a dry place, and under especially exposed climate conditions. The microscopes should be stored in their polystyrene box with a dessiccation bag.
 - Disassembly: A stereomicroscope is a finely-tooled precision instrument. If it is necessary to take it apart (for maintenance), the instrument should be returned to JEULIN or a qualified repair shop.

The shelf life of the LED is of 10,000 hours. In case of non-function of the lighting and after checking that the batteries are charged, send the microscope back to Jeulin.

5. After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
468 Rue Jacques Monod
CS 21900
27019 EVREUX CEDEX France

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries