Observation

Stéréomicroscopes

Observation

Stereomicroscopes

Réf :
571 287

Français – p 1

English – p 3

Version : 3107

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0 825 563 563 *

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1. Description

This monocular stereomicroscope is meant to observe any biological or geological material, which cannot be observed with a microscope due to its thickness and/or its opacity.

2. Technical specifications

Head	Monocular, 45° inclined, 360° rotatable
Eyepiece	Fixed WF10X
Dioptric adjustment	No
Objective	x 2
Magnification	x 20
Field of view	10 mm - x 20
Working distance	70 mm
Height adjustment	70 mm
Max thickness of subject	45 mm
Stage	50 mm diam. Disc, reversed black/white Metal clips with spring to maintain the subject
Illumination	No
Power supply	No
Delivered with	Dust cover

3. Set-up

Remove the stereomicroscope from its packaging.

Turn on the stereomicroscope: it is ready to work.

Place your subject on the white disc and fix it with the clips if necessary.

With the macrometric knob, adjust to get a clear image.

4. Maintenance

Protection from dust: After use, place the cover over the apparatus. If there are dust deposits on the lenses, blow them away with a syringe or dust remover (item no. 574 198) or use a martin hair brush; the lenses are fragile and can be scratched by dust.

- Fingerprints: Do not touch the lenses. If, however, they do have fingerprints on them, remove the fingerprints with a soft cloth or lens paper (item no. 803 092).
- Humidity: Humidity, especially in combination with temperature, is particularly harmful for assembled optics. It is suitable to store the microscopes in a dry place, and under especially exposed climate conditions. The stereomicroscopes should be stored in their polystyrene box with a dessiccation bag.
- Disassembly: A stereomicroscope is a finely-tooled precision instrument. If it is necessary to take it apart (for maintenance), the instrument should be returned to JEULIN or a qualified repair shop.

5. After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
468 Rue Jacques Monod
CS 21900
27019 EVREUX CEDEX France

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries