# Observation

Immunologie

## ***Microscope slides***

*Immunology*

Réf :  
575 578

Français – p 1

English – p 5

Version : 7001

**Groupage sanguin ABO et rhésus**

***Blood tests case***

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NOTES

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The **ABO** blood typing system is based on two criteria:

- one: the absence or presence of A, B and H antigens on the surfaces of the red blood cells,
- two: the absence or presence of anti-A and anti-B antibodies in the serum.

There are four possible cases:

blood group	A	B	AB	O
red blood cell antigens	A	B	A and B	H
serum antibodies	anti-B	anti-A		anti-A and anti-B

Consequently, the blood typing technique consists of two tests that are both required:

- determining the type of red blood cell antigens with test serums: Beth-Vincent test,
- determining the type of serum antibodies with red blood cell markers: Simonin test

Two kinds of reactions may occur:

- if the antigen is in the presence of the corresponding antibody, the red blood cells will agglutinate,
- if the antigen is not in the presence of the corresponding antibody, the red blood cells will not agglutinate.


The possible results are listed in the following table:

	reagents						
	test serums				red blood cells with antigens:		
	anti-A antibody	anti-B antibody	anti-A and anti-B antibodies	anti-H antibody	A	B	O
blood type A	+	-	+	-	-	+	-
blood type B	-	+	+	-	+	-	-
blood type AB	+	+	+	-	-	-	-
blood type O	-	-	-	+	+	+	-


Legend: + agglutination - absence of agglutination

## 2- Exemples :

1<sup>st</sup> line/6<sup>th</sup> column: one adds red blood cells with B-antigens to serum from type A blood: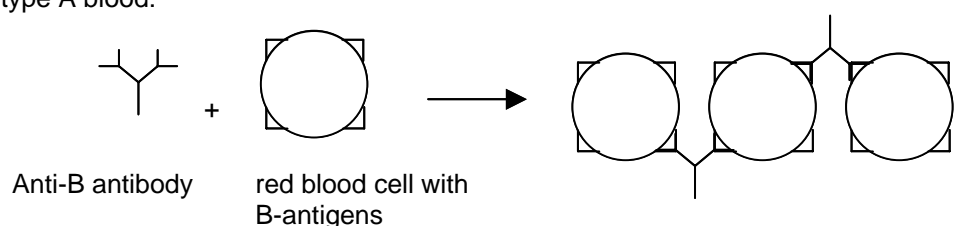

4<sup>th</sup> line/2<sup>nd</sup> column: one adds anti-B antibody test serum to red blood cells from type O blood: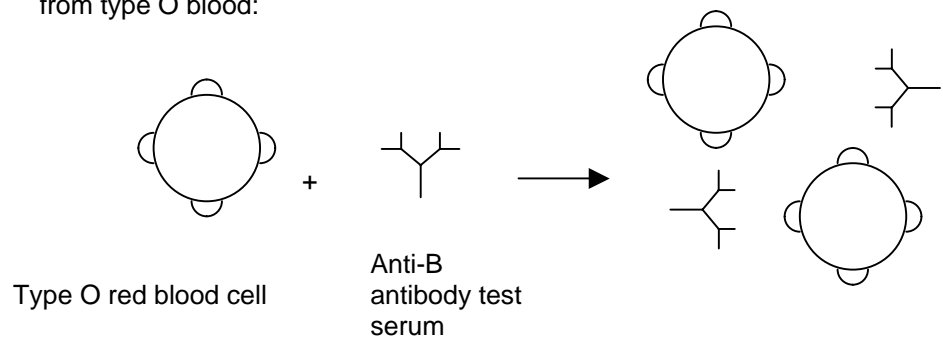

Given the seriousness of transfusion accidents, it is a good idea to classify all of the reactions in this table very precisely in order to conclusively determine the blood type classification: A, B, AB or O. If the slightest anomaly is observed in the seven tests (4 serum tests + 3 red blood cell antigens), the blood being examined is then subjected to more specific, supplementary analyses (type A2, A2B, Bombay: all very rare).



**The Rhesus blood** typing system is based on the absence or presence of the D-antigen on the surface of the red blood cells:

blood type	Rh+	Rh-
red blood cell antigen	D	

(Note: The anti-D antibody is not a normal blood constituent. Only induced anti-D antibodies are present in Rh- blood after immunization.)

Consequently, the blood typing technique (the Diamond method) involves only one test: the detection of the D-antigen on red blood cells by an anti-D antibody test serum.

There are therefore two possible cases:

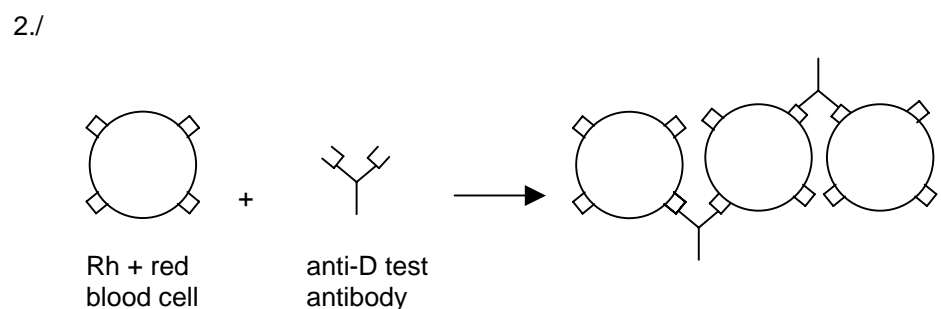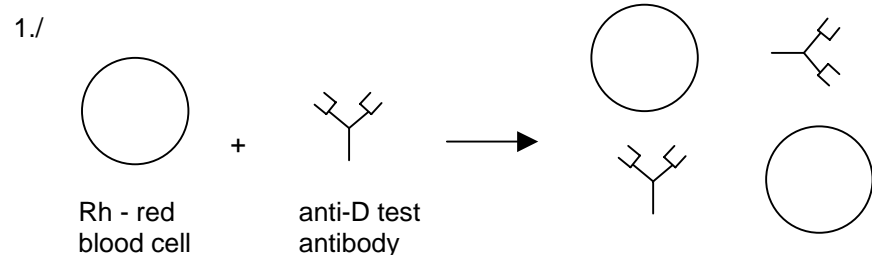The possible results are listed in the following table:

	reagent
	test serum: anti-D antibody
Rh - blood type	-
Rh + blood type	+

Note: If agglutination is not clearly observed (partial or delayed), the blood being examined is then subjected to more specific, supplementary analyses (Du: very rare).

In addition, in order to ensure the validity of the test (absence of auto-agglutination, specificity of the test serum, observation time), it is necessary to verify:

- the absence of agglutination of blood tested by BSA (serum albumin)
- the agglutination of Rh + red blood cells by the anti-D antibody test serum
- the absence of agglutination of Rh - red blood cells by the anti-D antibody test serum.

In the medical analysis laboratory, the observation of agglutinations is macroscopic and extemporaneous. **The NUBLAT MICROSCOPY LABORATORY (CAEN) has developed a technique (dilution, fixation, staining, mounting) for the microscopic and permanent observation of agglutinations.**

### 3- After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

**JEULIN - TECHNICAL SUPPORT**  
**Rue Jacques Monod**  
**BP 1900**  
**27 019 EVREUX CEDEX FRANCE**  
**+33 (0)2 32 29 40 50**

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*Rubrique FAQ*

## Direct connection for technical support

**A team of experts at your disposal from Monday to Friday (opening hours)**

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

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