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Chemistry measurement

**Ref :
701 058**

Français – p 1

English – p 5

Version : 9007

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Electronic scale 2000g/1g

1 Description

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1 Description

This electronic scale model is designed especially for students (high schools, prep schools, etc.). It only has two keys on the front, which makes it especially easy to use. This scale is made of resistant materials and it is also equipped with an internal overload protector. It can be run either on a 9V battery or on the main current with an adapter.

Photo 1
Electronic scale 2000g/1g.

- | | |
|----------------------------|------------------------|
| 1 – PVC weighing platform. | 4 – « OFF » key. |
| 2 – Washable front face. | 5 – « ON / TARE » key. |
| 3 – Digital display. | |

2 Specifications

Measurement range:	2000 g
Precision of the display:	1 g
Off-scale indicator:	« E » displayed on screen
Overload protector:	Sensor protected against excess weights up to 80 kg.
Automatic shutoff: (1)	Automatic shutoff after 90 seconds of idle time
Display:	12 mm LCD.
Dimensions (mm) :	(l×w×h) : 180 x 130 x 45 Ø tray: 105
Power:	Battery: 9V 6F22 Main: 230V/50Hz

(1) This function is automatically deactivated if the device is run on a 230V/50Hz main current with an adapter.

3 Startup and use instructions

a – Use precautions

This scale is not rated for commercial use. It therefore may not be used to weigh anything regulated by law.

This electronic scale is a precision instrument. In order to obtain precise measurements, observe the following recommendations:

- Place the apparatus on a clean, sturdy, level surface.
- When using the electronic scale, avoid vibrations, humidity, intense electromagnetic fields and disturbances, and severe temperature fluctuations.
- Allow the device to warm up for 30 seconds before performing measurements. This time is needed to allow the electronic circuits to stabilize.
- Handle this apparatus carefully. The weighing mechanism is protected against overloads; nevertheless, do not weigh objects with masses in excess of the weighing capacity of the scale. Gently place the objects to be weighed on the platform.

b – Power

This apparatus can either be battery operated or run on the main current.

Battery operation.

Use a 9V 6F22 battery (not included).

To insert the battery, turn the apparatus over and open the battery case cover (Figure 3). Plug the battery into the connector and replace the cover.

The battery needs to be replaced when the symbol « LO » appears on the screen.

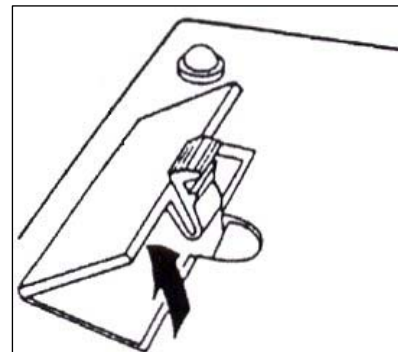
Figure 3

Using the main current.

Use a 230V AC / 9V DC current adapter (Ref. 281 286) with a central negative terminal (adapter not included).

Connect the adapter to the socket located on the back of the scale (Figure 4). The timed shutoff function is automatically deactivated when using the scale on the main current with an adapter.

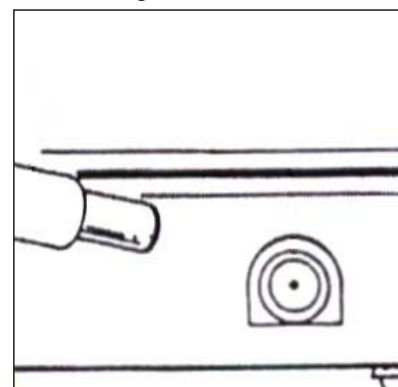
Figure 4

c – Turning the apparatus on and off

Pressing the « I/tare » key turns the scale on. The scale will then display « 8.8.8.8 », which corresponds to zeroing.

Wait until the scale displays « 0 », indicating that it is ready to weigh.

After use, turn the scale off by pressing the « 0 » key.

d – Tare function

Once the scale is running, place the empty container on the weighing platform. Then press the « I/TARE » key. The display will then indicate « 0 ». The « Δ » symbol will then be displayed on the screen, indicating that the tare function is on. Then weigh the desired mass of the substance in the container.

The combined mass of the container and its contents can be measured by pressing the « I/TARE » key again; the « Δ » symbole will then disappear from the screen.

5 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries