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Chemical measurement

**Ref :
703 262**

Français – p 1

English – p 3

Version : 6010

Sonde de température Pt 1000 Inox

Pt 1000 stainless steel

1 Description

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1 Description

The temperature probe Pt 1000 stainless steel is a very robust probe. It is easy to maintain and can replace the expansion thermometer in many applications.

Diagram 1

2 Technical characteristics

Temperature sensor	Pt 1000
Body of the probe	Stainless steel 316L
Cable	Silicon (withstands from -50 to +200 °C). Length: 140 cm.
Dimensions (mm)	Ø Body: 6 mm. Length: 135 mm
Connection	Male jack connector Ø 3.5 mm
Measurement range⁽¹⁾ :	- 50 to + 200 °C
Accuracy⁽¹⁾ :	+/- 0.15 °C + 0.002 T (T = temperature measured).
Stabilisation time:	< 20 s

⁽¹⁾ The measurement range and the accuracy specified here are those of the sensor alone. After connecting the probe to a measurement equipment, the measurement range and the accuracy are defined by the measurement system as a whole.

3 Precautions during use

The temperature probe Pt1000 stainless steel is designed for measuring the temperature of diluted solutions⁽¹⁾, the air, the soil, distillation apparatus...

It is not suitable for:

- Measuring the flame temperature (placing the probe in contact with a flame will result in the destruction of the sensor).
- Measuring the temperature of aggressive chemical products or concentrated solutions. The concentrated mineral acids (hydrochloric, sulphuric, nitric, phosphoric...) and the concentrated oxidants (hypochlorite...) can attack the materials, depending on the concentration, the temperature and the exposure time of the substance.

(1) The solutions of concentration equal to 1 mol.L⁻¹ (approx.) do not involve a risk of corrosion of the probe. It is however advisable to rinse the probe abundantly after each usage.

4 Connection

This probe is equipped with a male jack connector Ø 3.5 mm. It can be connected to any compatible measurement instrument and is interchangeable.

5 Service and maintenance

The probe does not require any particular maintenance. However, for ensuring its long life, it is important to follow these instructions:
Never leave the temperature probe Pt 1000 stainless steel immersed in a solution longer than necessary. Remove it from the solution and rinse it in demineralised water as soon as the measurement is complete.
Rinse the tip of the probe with demineralised water after each usage.
Store the probe in a dry place.

6 After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.
For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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Service gratuit * :
+ 33 (0)2 32 29 40 50

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Rubrique FAQ

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

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