**Ex.A.O.**

Mesure

***Data logging***

Measure

Réf :  
**454 051**

Français – p 1

English – p 3

Version : 5201

**Sonde oxygène air (collège)**

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*\* appel non surtaxé*

## 1. Description

This is a galvanic type sensor with built-in temperature compensation (5°C to 40 °C).

## 2. Installation and use

### 2.1 Connection

Connect the sensor onto the VTT oxymetry adapter via the BNC socket.  
The sensor can also be connected onto a self-contained INITIO oxymeter.

### 2.2 Calibration

#### For measurements in air

With the sensor being placed in air:

- Select the oxymeter position "Measurement in air".
- Set the oxymeter to 21 % (O<sub>2</sub> concentration in ambient air) via the potentiometer located on the adapter housing.

## 3. Technical specifications

- Principle: galvanic type sensor.
- Response time: approx. 15 seconds.
- Usable life: approx. 5 years (stored in air at 20°C).
- Temperature compensation: from 5°C to 40°C.
- Influence of other gases: almost no effect of CO<sub>2</sub> (both gaseous and dissolved), CO, H<sub>2</sub>S, SO<sub>2</sub>.
- The sensor may be affected by high ammonia concentrations, and by ozone.
- Operating pressure: from 800 to 1200 hPa absolute. Note: the output signal variation is proportional to pressure.
- Storage temperature: from - 20°C to 60°C (the electrolyte freezes at -20°C).

## 4. Precautions for use

The electrolyte is isolated from the ambient medium by a thin membrane (behind the protecting grid). In case of membrane damage, the sensor would be definitely unusable.

The tightening or loosening of the sensor head also affects the output signal.  
This may result in the sensor being definitely unusable. Doing so shall render null and void the contractual two-year warranty granted by JEULIN.  
In case of contact with the electrolyte, flush thoroughly with fresh water.

## **5. Maintenance**

Clean with fresh water, avoid strong water jets onto the membrane

## **6. After-sales service**

The device is under a 2-year guarantee, it must be sent back to our workshops.  
For any repairs, adjustments or spare parts please contact:

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