Déclaration de Conformité CE & RoHS
CE & RoHS conformity declaration
Declaración de conformidad CE & RoHS
Declaração de coformidade CE & RoHS

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Regulador de temperatura: 000RG011.

1.- INTRODUCTION

First of all we would like to thank you for your trust in acquiring this **LSCI** equipment. We are sure you will be satisfied with it and its intended use, and you will speak of it highly to other possible customers. From now on, you will have all **LSCI** personal technical and material support for which we give our e-mail:

2.- GUARANTEE

This equipment is warranted for 36 months against any manufacturing defect or defective part, except the parts with a limited hours of life (Lower than 36 months), ever indicated by the parts manufacturer (Motors, bearings, belts, joints, lamps etc.), in this case **LSCI** guarantee just the manufacturer's given period. Please contact your **LSCI** distributor for servicing.

This guarantee is only valable if:

- A) The unit is used for the intended use.
- B) The unit is only manipulated by authorized technical engineers.
- C) The unit is used between the prefixed MWC "Maximum Working Conditions"
- D) The Guarantee form is filled and sent to us by: e-mail, fax, post or **web registry**

LSCI reserves the right to introduce any changes or modifications without previous notice.

Request a Return Goods Authorization (RGA) number before returning any material to **LSCI**. This action prevents delays if acting upon your request any eventual problem with custom clearance.

3 YEARS OF GUARANTEE

When the life of the unit is finished, do not throw it away. Directive 2002/96/EC 27-01-2003 (WEE)

3.- SAFETY

LSCI will be not responsible for any accident or damage produced by the use incorrect or not, or faillure of this product.

GENERAL INSTRUCTIONS TO YOUR OWN PROTECTION

A) Read this operating instructions fully before starting up, and follow these safety instructions.

B) Keep the operating instructions in a place where they can be accessed by everyone.

C) Ensure that only trained people work with this unit.

D) Ever follow the safety instructions, guidelines, occupational health of each country.

E) Ensure that the electric socket is earthed and allows to supply the intensity and voltage required by this unit

F) This unit is not protected against explosions.

G) Keep it separated from a water source conveniently.

H) Keep it separated from an external heat focus conveniently.

I) Before any accessory installation, unplug the unit of the electricity.

J) For a complete disconnection of the electricity, unplug of the power, the main cable.

K) The power supply socket must be placed in a place with an easy and fast access.

L) To unplug the unit of the socket, never pull of the cable, hold the connector and pull it.

M) Use the basic elements for personal protection such as lab. protective glasses, lab coat, masks, protective foots and gloves.

N) Check that the laboratory is provided of the basic safety elements

O) If the experiment generate fumes or dangerous vapours, use a fumes extraction hood.

P) **For educational purposes, this apparatus just must be used by a teacher or under a teacher supervision**

Specific cautions for the apparatus generators of MAGNETICS FIELDS

The effects of the magnetic field are dangerous in few systems e.g. Data carriers, credit cards, cardiac pacemakers, navigation devices, etc

Specific cautions for the MAGNETIC STIRRERS

Splashing dangerous, corrosive, contaminant, biopathogenic and radioactive liquids. To prevent splashing be controlled the speed, the container stability, it must be without horizontal displacement on the working surface.

Choose an adequate stirring bar according to the shape and size container.

Don't introduce the magnetic bar in the container liquid, when the stirring is on, it difficult the magnetic coupling

Do not operate with the stirring without any recipient (No load condition)

The powder generated as consequence of the abrasion process of the magnetic bars when running, can react with the stirred samples.

Specific cautions for the MECANIC MIXERS, SHAKERS, STIRRERS

Do not touch or handle the objects, devices or recipients during the stirring process, it's necessary to stop the motion (Except devices provided with clutch). Be special care with the ties, sleeves and other loose parts of the clothings, do not are in contact with the parts in motion.

Specific cautions for the devices with UV-254 RADIATION

To operate with devices or accessories with UV254nm radiation, it's necessary to use special glasses of complet protection. Be care full with the light reflections as well.

Specific cautions for the devices operating under PRESSION OR VACUUM

Do not operate with presson or vacuum bigger than the resistance of the recipients.

It is necessary to use protection glasses and safety screen.

Specific cautions for HEATING EQUIPMENTS AND DEVICES

Heating Magnetic Stirrers, Hot Plates, Heating elements, Thermostatic Baths, Heating Mantles, Dry Blocks RISK OF BURNS

Exercise caution when touching the housings, hot plates or the heating elements, they can reach high temperature, pay special attention to the residual heat after switching off.

Ensure that the mains power cable isn't in direct contact with any heated element.

Wear your personal protective equipment according to the hazard category of the medium to be processed, otherwise there is a risk of:

Splashing heated liquids.

Projection of projectile solid parts.

Releasing any toxic or combustible gases.

Set the unit in a spacious area on an even, stable, clean, non-slip, dry and the surface must be fireproof.

The feet of the unit must be clean and undamaged.

Check the unit and accessories beforehand for damage each time you use them. Do not use damaged components, parts and accessories.

Increase the speed gradually.

Reduce the speed in the following situations:

If the liquid splashes out of the recipient because the speed is too high.

If the liquid is not running smoothly.

If the recipient moves on the set-up surface.

Handle with care and take caution when process and heat in the following situations:

Processing any product with low flash point, there are danger of ignition.

Processing flammable materials.

Use of glass recipients with incorrect size.

The level of the liquid is too high.

The recipient shape is unsafe.

The recipient position is unsafe.

The processed product can heat up it self when is processed.

The top surface can heat up, (Without using the heating system), due to the action of the internal drive magnets (Foucault effect), when are used at high speed.

Do not use this unit in explosive atmospheres, with hazardous products or under water (ExceptANS-008).

Just process substances that will not react dangerously to the extra energy produced through processing.

This also applicable to any extra energy produced by other ways, e.g by light irradiation.

Do not cover the unit, even partially, and in special the heating plate and the cooling rings and windows, with metallic parts, e.g. Aluminium film, to prevent overheating, **it can destroy the unit**.

Ensure that the working top surface is kept clean.

Do not operate with the unit in a heating way without any recipient on the heating plate it can overheat.

Observe the minimum distances between the devices 100mm, and 100mm between the rear of the unit and the wall; 800mm between the unit top and the roof.

Follow the operating instructions of the used accessories.

Ensure that the external accessories must be securely attached and fixed to the unit, and cannot come off by themselves. Their gravity centre must lie within the set-up surface.

Ensure that the external temperature sensor is inserted in the liquid to a depth at least 20mm.

The safe operation work of the unit, is just guaranted if are used **LSCI** accessories.

Danger when heating products with a low flashing point.

Danger when heating products with a low decomposition temperature point.

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5.-UNPACKING

- . Unpack with care and eliminate the rests of packing material and dust.
- . Ensure that the unit and the accessories are not damaged.
- . Check the packing list in order to know if the equipment is complet.
- . If some part is damaged or incomplete, please contact with your **LSCI** dealer.

6.- DESCRIPTION

It's a heating plate of low twmperature, specially developed to heat biological devices, as slides. microtiter plates, etc

7.-INTENDED USE

- . Place the unit on a laboratory work bench or on a flat surface.
- . Check that the power supply is according with the unit.
- . Check that the main switch placed on the rear panel is OFF (Down position).
- . Connect the main cable, between the unit and the electric shocket.
- . Switch ON the main switch.
- . Press the SET key.
- . Bymeans of the UP & DOWN keys select the working temperature.
- . Press the SET key to confirm.

Attention: Don't keep pressed the SET key more than 5 seconds, it means to open the configuration mode, and can unconfigure the regulator.

Following any interruption of the power supply, the unit will automatically restart in same conditions, previous to the interruption.

8.-MAINTENANCE

- . This unit doesn't need any maintenance, except the periodic cleaning; just use a rag impregnated of cleaning glass liquid not agresive, ecologically correct. Before disconnect the main plug.
- . If a radioactive or biologic contamination is produced, you must contact with an authorized and specialized company.
- . The unit may only be opened by experts electronic technics.
- . Before open disconnect the main plug. **DANGER OF ELECTRO-SHOCK.**
- . Before any maintenace disconnect the main plug.
- . The fuse is placed on the rear part next to the main connector, just replace by another exact (T1A).

9.- SPARE PARTS

To order spare parts, please give the model reference and the serial number.

Cable: 000CN103, Switch: 000CN316, Heating element: 000RC300, Temperature regulator: 000RG011,

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